



CATALOG

GVS Life Sciences 2014





SUMMARY

GVS GROUP	2
1. MICROFILTRATION FILTER DEVICES	4
1.1 - Syringe-Filters	4
1.2 - Vent Filters	16
1.3 - Capsule Filters	18
1.4 - Centrifuge Filters - Centrex	20
1.5 - Bottle-top Filters - ZapCap	21
1.6 - Extractor - Ethidium bromide	21
1.7 - Filter Holders for Membranes	22
1.8 - Mini Vial	25
2. MICROBIOLOGY	27
2.1 - Microbiological monitors & Analytical Funnels	27
2.2 - Savur Analytical Filter Unit	28
2.3 - Nutrient Liquid Media	29
2.4 - SwabCheck	33
2.5 - Dilution Bottles	36
3. FILTRATION MEMBRANES	38
3.1 - Membranes for Filtration	38
3.1.1 - Cellulose Acetate (CA) Membrane	38
3.1.2 - Nitrocellulose (NC) Membrane	39
3.1.3 - Nylon (NY) Membrane	41
3.1.4 - Polyethersulfone (PES) Membrane	43
3.1.5 - Polypropylene (PP) Membrane	44
3.1.6 - Polytetrafluoroethylene (PTFE) Membrane	45
3.1.7 - Polytetrafluoroethylene (PTFE) Hydrophilic Membrane	46
3.1.8 - Regenerated Cellulose (RC) Membrane	46
3.1.9 - Polyvinylidene Fluoride (PVDF) Membrane	46
3.1.10 - Polycarbonate Track Etched (PCTE) Membrane	47
3.1.11 - Polyester Track Etched (PETE) Membrane	51
3.1.12 - Drain Disc	52
3.1.13 - Silver Membrane.....	53
3.2 - Filter Papers	54
3.2.1 - Qualitative Papers	54
3.2.2 - Quantitative Papers	56
3.2.3 - Glass Microfiber	59
3.3 - Glass Fiber Filters with Binder	60
4. TRANSFER MEMBRANES	62
4.1 - Neutral Nylon	62
4.2 - Reprobing Charged Nylon	63
4.3 - Pure Nitrocellulose (MCE)	64
4.4 - Supported Nitrocellulose (MCE)	65
4.5 - Polyvinylidene Fluoride (PVDF)	66
5. FAST - protein array solutions	68
5.1 - FAST Slides	70
5.2 - FAST Pack	71
5.3 - FAST Buffers	71
5.4 - FAST Quant	72
5.5 - FAST Services	73
5.6 - FAST Accessories	74
FILTRATION GUIDE	75
COMPATIBILITY CHART	80
PRODUCT CODE INDEX	82

GVS Life Sciences

With over 30 years of experience GVS Group is one of the world's leading manufacturers of microfiltration devices with applications in the laboratory, medical and pharmaceutical markets.

Medical Filters & Components

The origins of GVS initially focused on medical filters for blood and IV solutions. Today GVS Life Sciences provides a wide range of innovative products including standard and custom devices for laboratory filtration, sample preparation, clarification or fine particulate removal for laboratory applications, protein and nucleic acid binding and analysis and biomarker research.

GVS Life Sciences offers a complete range of

Microfiltration products: Syringe Filters, Vent Filters, Capsule Filters, Centrifuge Filters, Bottle Top, Filter Holders for Membranes, Filter Funnels

Microbiology products: Microbiological Monitors, Analytical Monitors, Nutrient Liquid Media, Swab Kits, Dilution Bottles

Filtration Membranes: Membranes for Filtration (Discs, Sheets, Rolls), Filter Papers

Molecular Transfer Membranes: Nitrocellulose, PVDF, Nylon

Protein Microarray Products: FAST® Slides and Arrays

GVS Life Sciences also offers OEM production of membranes and components.

International expansion

GVS Group's presence in major markets across the world has led to the opening of 16 manufacturing plants located in - Italy (2), UK (5), Brazil (1), USA (4), China (3) and Romania (1), as well as offices in Italy, Germany, UK, Spain, Serbia, USA, México, Brazil, Argentina, China, Japan, Korea, India and Malaysia.

Sophisticated industrial Technology.

GVS's highly innovative technical capabilities include filter material development, hydrophobic and hydrophilic technology, activated carbon filtration, filter surface coating technology, chemical / biological lab for test plastic filter and metal test and polymer development.

Production technologies include: Multi-cavity insert and over-molding, pleating, potting and low compression injection, high-speed automatic assembly, ultrasonic, heat and radio-frequency welding, laser cutting and welding and All in-Mold technology, a revolutionary manufacturing technology combining injection molding and robotic assembly all within the molding tool. All of the products are manufactured in a clean-room environment.

Commitment to Quality.

GVS has obtained ISO 9001:2010 certification, and our Medical Division has qualified for ISO 13485: 2003 certification, plus several of our medical devices have been qualified for CE marking according to E.U. Directive 93/42/EEC. Recently, our Italian headquarters successfully achieved UNI EN ISO 14001:2004 certification for its Environmental Management System (EMS), marking a milestone in GVS Life Sciences's ongoing commitment to reduce its environmental footprint.

Research & Development.

A great part of the know-how incorporated in GVS's products comes from its Research Lab, which ensures that the company's various divisions have access to innovative R&D. With its pioneering tools and facilities and highly sophisticated analytic techniques, this lab also works in close conjunction with a large number of hospitals and academic bodies of international acclaim, in Italy, in the UK and wherever GVS Life Sciences operates. Without it, the group's strongly innovation-oriented policy and commitment to growth would be much less effective.

Maine Manufacturing

Maine Manufacturing LLC joined GVS Group in 2013, offering a full line of laboratory filtration and analysis products. GVS Life Sciences and Maine work together to provide an incomparable experience for customer in the analytical, chemistry and life science laboratories. Maine Manufacturing is a global supplier of membrane-based solutions for life science research. We offer a full line of research tools for filtration, protein and nucleic acid blotting and transfer. Maine Manufacturing also produces custom and OEM solutions for diagnostics, multiplexed immunoassays and sample preparation and analysis.

Your partner for practical laboratory solutions.



MICROFILTRATION



MICROFILTRATION

1. MICROFILTRATION FILTER DEVICES

INTRODUCTION

GVS Life Sciences offers a comprehensive range of disposable syringe filter devices designed to provide fast and efficient filtration of aqueous and organic solutions. They are made with a wide variety of membrane filters. These syringe filters are suitable for numerous applications in pharmaceutical, environmental, biotechnology, food/beverage, and agricultural testing laboratories.



1.1 Syringe-Filters

GVS Life Sciences offer a full range of disposable syringe filter devices designed to provide fast and efficient filtration of aqueous and organic solutions, aerosol separation and venting.

These syringe filters are suitable for numerous applications in pharmaceutical, environmental, biotechnology, food/beverage, and agricultural testing laboratories.

- Polypropylene or Acrylic housing
- Multifunctional connectors - equipped with luer-lock or luer-slip female connections
- Sterile or non-sterile options
- Available in bulk-packages or individual blisters
- Customized product and packaging on request

Filtration Membrane Types

Microporous membranes are commonly used in filtration devices for sample preparation, purification and sterilization. The most common membrane types are described here with physical properties and applications.

Cellulose Acetate (CA)

Hydrophilic membrane. Low protein binding, Ideal for protein, cell culture media and enzymes filtrations, tissue culture media sterilization, biological fluid filtration and other filtration applications where maximum recovery of proteins is critical.

Features and Benefits

- Low protein binding, 3.8 $\mu\text{g}/\text{cm}^2$
- Hydrophilic
- High throughput
- Superior strength and stability
- Uniform pore structure, consistent flow rates
- Burst strength of 130 psi

Applications

- Protein and enzyme filtration
- Biological fluid filtration sterilization
- Tissue culture media sterilization
- Clarification of aqueous and alcohol solutions
- Cell Culture

Detailed Application: Biological sample preparation; Very low protein retention: protein and enzyme filtration with very low protein retention. Filtration of protein-containing solutions with minimal protein loss; Sterilization and clarification; Receptor binding studies; Tissue culture media sterilization; Particle removal of cellular constituents from solution: The 0.22 μm membrane is the filter of choice for sterile filtration of aqueous solutions such as nutrient media, buffers and sera; The 0.45 μm membrane is a very convenient filter type for the reduction of particles and microorganisms in aqueous solutions such as nutrient media, buffers and sera. Enhanced recovery of fastidious gram positive organisms. The 0.8 and 1.2 μm membranes are used for the particle filtration in samples where a low adsorption is required. The 5.0 μm membrane is used for the coarse particle filtration in samples where a low adsorption is required.

1.1 Syringe-Filters

Nylon (NY)

Hydrophilic membrane. Ideal for use in general filtration or medical assays. Superior strength, resistant to a range of organic solvents. Low extractables. High protein binding capacity. Lot-to-lot consistency.

Features and Benefits

- Naturally hydrophilic
- Wide chemical compatibility range
- Strength and dimensional stability
- Low extractables

Detailed Application: Ideally for use in general filtration and HPLC sample preparation of aqueous, solvent-based and organic solutions prior to chromatography or other instrument analyses; Excellent chemical compatibility with esters, bases, and alcohols; Biological sample preparation; High protein binding capacity; Immobilizes antigens, antibodies and other protein.

Applications

- Sterilization, clarification of aqueous and organic solvent solutions
- HPLC sample preparation
- Chromatography
- Hydraulic Fluids and Machined Parts

Polyethersulfone (PES)

Hydrophilic membrane. designed to remove particulates during general filtration, low protein and drug binding characteristics make it ideally suited for use in life science applications. Its strength and durability are advantageous during usage that involves aggressive handling or auto- mated equipment. Low protein and drug binding characteristics maximize recovery of critical drugs used in I.V. therapy, chemotherapy and open-heart surgery.

Features and Benefits

- Very Low Protein Binding
- Fast Flow Rates
- Low Extractables
- Wide Chemical Compatibility Range
- Strength and Dimensional Stability
- Autoclavable

Detailed Application: Ideally for use in life science applications. Preparation of aqueous, biological or protein based solutions prior to chromatography or other instrument analyses. Accurate results for the most sensitive analysis of ionic species. Specifically designed for IC applications. Low drug and protein binding for pharmaceutical filtration.

Applications

- Protein and enzyme filtration sterilization
- Biological fluid filtration sterilization
- Tissue culture media sterilization
- Pharmaceutical sterilizing filtration
- Environmental water studies
- Filtration of Aqueous Solutions
- Analytical Sample Prep, uHPLC
- IC Chromatography
- Sterile Filtration and Clarification
- Cell Culture

Asymmetric Polyethersulfone (PES)

Asymmetric PES enables the fast filtration of aqueous solutions with greater throughput. The PES membrane has low drug and protein binding properties with aqueous solutions. Low ion and metals extractables provide for ideal analysis by ion chromatography and ICP MS. Asymmetric PES is provided both sterile and non-sterile.

Features and Benefits

- Fast flow rate
- High filter capacity
- Low protein binding
- Extended filtration capacity and lifetime
- Higher particle loads and protein concentrations tolerance

Applications

- Prefiltration and Clarification
- Liquid filtration and sterilization
- Ion chromatography
- Gas separation

Nitrocellulose Mixed Esters (MCE)

Hydrophilic membrane. Aqueous clarification and particulate capture. Consistent high flow rate for faster filtration. Uniform pore structure for selectivity. Hydrophilic, inert cellulose nitrate. High binding capacity. Manufactured thickness within 10 microns.

Features and Benefits

- Hydrophilic for aqueous clarification and particulate capture
- Consistent high flow rate for faster filtration
- Uniform pore structure for selectivity
- Hydrophilic, inert cellulose nitrate
- High binding capacity
- Manufactured thickness within 10 microns

Applications

- Filtration of Aqueous and Organic Solutions
- Analytical Sample Prep, uHPLC
- Chromatography
- Clarification

Detailed Application: Used to clarify aqueous samples especially in water studies. Disposable syringe filters for wastewater, food and beverage filtration. Higher protein binding than cellulose acetate for most proteins. HCG; HIV; LH; Chlamydia. Drugs of Abuse Pathogenic Microorganisms, Environmental Contaminants.

MICROFILTRATION

Polyvinylidene Difluoride (PVDF)

Hydrophilic membrane. Ideal for use in Sterilizing and Clarifying filtration of biological solutions. High Flow Rates. Low Extractables. Broad Chemical Compatibility. Very low protein binding.

Applications

- Filtration of Aqueous and Organic Solutions
- Analytical Sample Prep, uHPLC
- Chromatography
- Clarification

Detailed Application: Sterilizing and clarifying filtration of biological solutions. Preparation of protein-containing solutions prior to chromatography or other instrument analyses. Useful for a wide range of applications, including aggressive and non-aggressive solvent-based mobile phase. Offers excellent chemical compatibility, even with aggressive acids and alcohols. Provides high flow rates and throughput, low extractables and broad chemical compatibility. The 0.22 μm pore size assures removal of the contaminant size most likely to plug a UHPLC column and offers significantly better protection of your analytical results.

Polytetrafluoroethylene (PTFE)

Hydrophobic membrane. Ideal for filtration of strong acids and aggressive solutions, venting applications, phase separations, aerosol samplings. Chemically and biologically inert. Superior chemical resistance. Can withstand high temperatures.

Features and Benefits

- Naturally hydrophobic
- Compatible with strong acids and aggressive solutions
- Improved durability and handling

Detailed Application: Preparation of organic solutions prior to chromatography or other instrument analyses. Exceptional chemical and temperature compatibility. Filtration of strong acids and aggressive solutions. Venting applications. Ideal for filtration of gas and/or HPLC organic solvents, aggressive solutions and small venting. Ultimate in chemical compatibility for filtering harsh chemicals that destroy other membrane materials. For solvents (acetone, dimethylformamide, DMSO...) or for aggressive aqueous liquids.

Applications

- Filtration of strong acids and aggressive solutions
- Venting applications
- Phase separations
- Aerosol sampling

Hydrophilic Polytetrafluoroethylene (PTFE 'HP')

PTFE HP membrane is compatible with organic solvents, acids, and basic solutions. Hydrophilic PTFE membrane has low drug and protein binding properties with excellent aqueous and solvent compatibility. High sample recoveries and low ion and UV extractables provide for ideal analysis by uHPLC and LC/MS.

Features and Benefits

- No need to pre-wet the membrane
- No need to flush membrane of pre-wetting chemicals
- No pre-wetting means production time reduction
- Reduce potential interference with biological processes
- Longer shelf life because the filters are stored and shipped dry

Detailed Application: The PTFE HP has extremely low levels of extractables. PTFE HP syringe filters will not contribute extractables that will interfere with the ionization process. Excellent chemical resistance – Use this universal filter for all your uHPLC, Ion Chromatography and LCMS samples. The Hydrophilic PTFE membrane can be used with both organic and aqueous solvents. When used within a polypropylene housing, the membrane offers excellent chemical resistance.

Applications

- Molecular identification
- Structural determination
- Pharmacokinetics
- Drug discovery and development
- Drug testing
- Environmental monitoring
- Food safety monitoring
- Oil composition determination

Glass Fiber (GF)

Hydrophilic material. Used also as a pre-filter to extend membrane life. Eliminate sample contamination. Excellent wet strength for each handling and filter integrity. Ideal for water/air pollution analysis, liquid clarification and cell harvesting.

Features and Benefits

- Acrylic binder
- High dirt holding capacity
- Biologically inert
- Bonding reduces media migration

Applications

- Filtration of Aqueous and Organic Solutions
- Analytical Sample Prep, uHPLC
- Difficult to Filter Solutions
- Fuel Hydraulic Fluids and Machined Parts

Regenerated Cellulose (RC)

Hydrophilic membrane. Resistant to a very wide range of solvents. Suitable for use with either aqueous solutions or organic solvents. Compatible with HPLC solvents. Very low protein binding capacity and hence excellent for protein recovery applications.

Applications

- Filtration of Aqueous and Organic Solutions
- Analytical Sample Prep, uHPLC

Detailed Application: Extremely versatile can be used with almost any laboratory procedure. Hydrophilic, solvent-resistant membrane. Extremely chemically resistant: suitable for use with aqueous solutions and resistant to a very wide range of solvents. Compatible with HPLC solvents: for simple, rapid and reliable ultra cleaning of samples for the HPLC or GC analysis. Chemically resistant with both aqueous and organic samples. Very low binding coefficient.

- Chromatography
- Clarification
- Protein Chemistry

1.1 Syringe-Filters

Polyethylene (PE)

An "Universal" Filter for all analytical filtration requirements. For aqueous and aggressive organic solvent filtration. Highly recommended for filtering HPLC samples: HPLC sample preparation requiring low detection levels. Wide application in sample preparation. Ion chromatography.

Polypropylene (PP)


GVS Life Sciences polypropylene membrane filters are composed of pure polypropylene with absolute pore size ratings. These filters offer broad chemical compatibility allowing its use with aqueous and organic solvents samples. The polypropylene filter has extremely low extractable levels designed to provide accurate, consistent analysis results for sensitive ion chromatography applications while prolonging column life.

Applications

- Aqueous and organic solvent filtration
- HPLC sample preparation requiring low detection levels
- Ion chromatography
- Total digest for heavy metals

MICROFILTRATION

Quick application guide

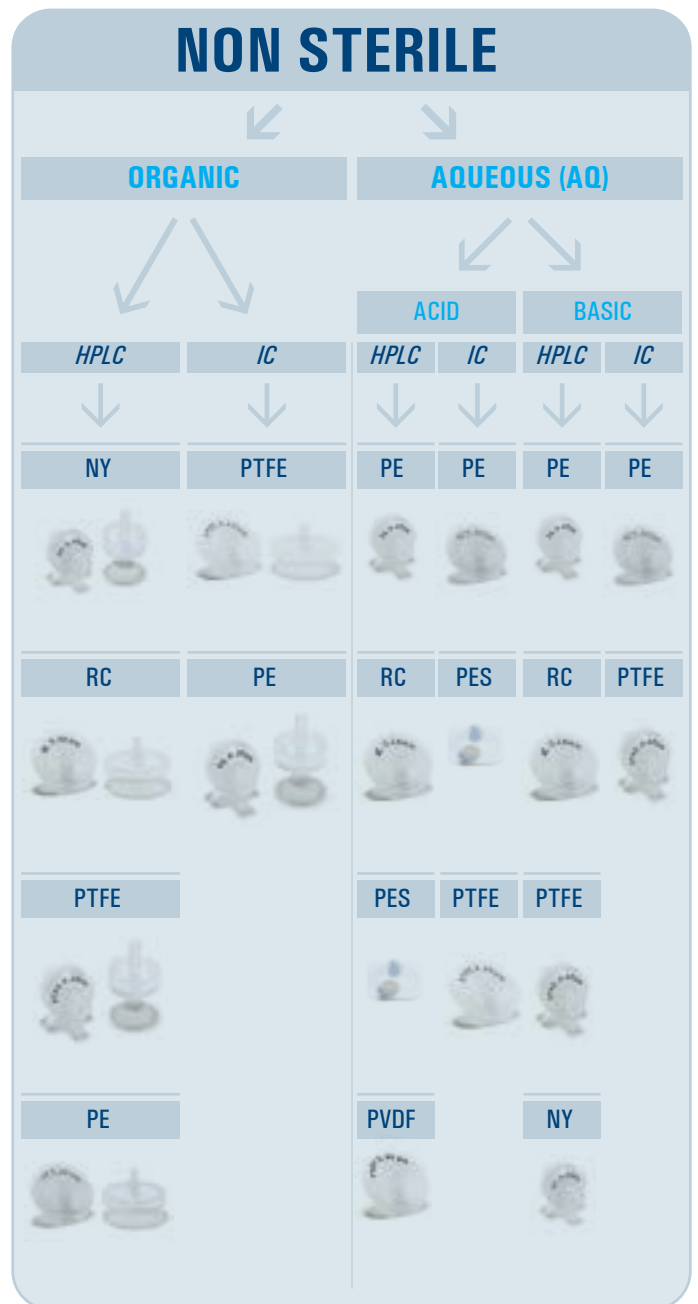

1.1 Syringe-Filters

1.1.1 - 13 mm Syringe Filters ABLUO®



Features and Benefits

Housing Material: Acrylic and Polypropylene

Membrane Material: Cellulose Acetate, Nylon, PES, Nitrocellulose Mixed Esters, Regenerated Cellulose, PVDF, PTFE and PE

Membrane Diameter: 13 mm

Housing Diameter: 18 mm

Effective Filtration Area: 0.6 cm²

Pressure Bar: 5

Sterile: No

Inlet / outlet: FLL/MLL-MLS

Applications

- Filtration of Aqueous, Organic and Alcohol Solutions
- Analytical Sample Prep, uHPLC
- IC Chromatography
- Fuel Hydraulic Fluids and Machined Parts
- Clarification
- Protein Chemistry
- Cell Culture

Specification

- Quick and efficient filtration of samples and all kind of solutions, solvents or gases
- Housing injected in Polypropylene or Acrylic
- Multifunctional Syringe Filters: equipped with luer-lock or luer-slip male connections for different applications
- Sterilized or non-sterile
- Available in bulk-packages or individual blisters
- Customized product and packaging on request

Ordering information

Membrane Material	Pore Size (µm)	Housing Material	Color	Product Code
				Packaging 500/pk
Cellulose Acetate (CA)	0.22	Acrylic	Blue	FJ13ANCCA002DD01
Cellulose Acetate (CA)	0.45	Acrylic	Yellow	FJ13ANCCA004FD01
Cellulose Acetate (CA)	0.80	Acrylic	Green	FJ13ANCCA008ED01
Cellulose Acetate (CA)	1.20	Acrylic	Red	FJ13ANCCA012CD01
Cellulose Acetate (CA)	5.0	Acrylic	Brown	FJ13ANCCA050PD01
Nylon (NY)	0.20	Polypropylene	Transparent	FJ13BNPNY002AD01
Nylon (NY)	0.45	Polypropylene	Transparent	FJ13BNPNY004AD01
Polyethersulfone (PES)	0.22	Polypropylene	Transparent	FJ13BNPPS002AD01
Polyethersulfone (PES)	0.45	Polypropylene	Transparent	FJ13BNPPS004AD01
Nitrocellulose Mixed Esters (MCE)	0.22	Acrylic	Transparent	FJ13BNCNC002AD01
Nitrocellulose Mixed Esters (MCE)	0.45	Acrylic	Transparent	FJ13BNCNC004AD01
Regenerated Cellulose (RC)	0.20	Polypropylene	Transparent	FJ13BNPRC002AD01
Regenerated Cellulose (RC)	0.45	Polypropylene	Transparent	FJ13BNPRC004AD01
Polyvinylidene Fluoride (PVDF)	0.22	Polypropylene	Transparent	FJ13BNPPV002AD01
Polyvinylidene Fluoride (PVDF)	0.45	Polypropylene	Transparent	FJ13BNPPV004AD01
Polytetrafluoroethylene (PTFE)	0.20	Polypropylene	Transparent	FJ13BNPPT002AD01
Polytetrafluoroethylene (PTFE)	0.45	Polypropylene	Transparent	FJ13BNPPT004AD01
Polytetrafluoroethylene Hydrophilic (PTFE HP)	0.22	Polypropylene	Transparent	FJ13BNPPH002AD01
Polyethylene (PE)	0.20	Polypropylene	Transparent	FJ13BNPPE002AD01
Polyethylene (PE)	0.50	Polypropylene	Transparent	FJ13BNPPE005AD01

MICROFILTRATION

13 mm Syringe Filters - Sterile ABLUO®



Features and Benefits

Housing Material: Acrylic
Membrane Material: Cellulose Acetate, PES and PVDF
Membrane Diameter: 13 mm
Housing Diameter: 18 mm
Effective Filtration Area: 0.6 cm²
Pressure Bar: 5
Sterile: Yes
Inlet / outlet: FLL/MLL-MLS

Applications

- Filtration of Aqueous Solutions
- Analytical Sample Prep, uHPLC
- IC Chromatography
- Sterile Filtration and Clarification
- Protein Chemistry
- Cell Culture
- Clarification

Ordering information

Membrane Material	Pore Size (μm)	Color	Product Code
			Packaging 50/pk
Cellulose Acetate (CA)	0.22	Blue	FJ13ASCCA002DL01
Cellulose Acetate (CA)	0.45	Yellow	FJ13ASCCA004FL01
Cellulose Acetate (CA)	0.80	Green	FJ13ASCCA008EL01
Cellulose Acetate (CA)	1.20	Red	FJ13ASCCA012CL01
Cellulose Acetate (CA)	5.0	Brown	FJ13ASCCA050PL01
Polyethersulfone (PES)	0.22	Transparent	FJ13BSCPS002AL01
Polyethersulfone (PES)	0.45	Transparent	FJ13BSCPS004AL01
Polyvinylidene Fluoride (PVDF)	0.22	Transparent	FJ13BSCPV002AL01
Polyvinylidene Fluoride (PVDF)	0.45	Transparent	FJ13BSCPV004AL01

1.1 Syringe-Filters

1.1.2 - 17 mm Syringe Filters CAMEO®



Sample volume: < 12 mL without prefilter; < 18 mL with prefilter
Operating parameters: Maximum Operating Temperature 82°C/180°F, Maximum Operating Pressure 130psi
Inlet / outlet: FLL-MLS

Description and Use

GVS Life Sciences offers a comprehensive range of disposable syringe filter devices designed to provide fast and efficient filtration of aqueous and organic solutions. They are made with a wide variety of membrane filters with a polypropylene housing. These syringe filters are suitable for numerous applications in pharmaceutical, environmental, biotechnology, food/ beverage, and agricultural testing laboratories.

Features and Benefits

Housing: Heat-sealed pure polypropylene without the use of glues or sealants
Media: Nylon, PTFE, PES, PVDF, glass fiber, cellulose acetate and polypropylene
Autoclavable: Yes
Prefilter: 1.0 micron binderless glass-fiber
Filtration area: 1.4 cm²
Housing diameter: 22mm
Membrane diameter: 17mm
Holdup volume: < 15 microliter

Applications

- HPLC sample preparation
- Dissolution testing
- Content uniformity
- Environmental samples
- Composite assays
- Food analysis
- Biofuel analysis

Specification

- Increased throughput and speed of sample preparation
- Lower hold-up volume due to an improved flow channel design and reduced spacing between the supports within the housing for better handling of small sample volumes or costly samples.
- Increases operating pressure up to 130 psi due to the overmold that prevents sample leaking at the seam and keeps the filter unit from bursting in half.
- Strict quality control: syringe filters are integrity-tested to ensure a proper filter fit and weld to eliminate any potential filter by-pass
- Accurate labeling: each filter is labeled with the specific filter material and pore size for easy identification even if the syringe filter is not in its original packaging

Ordering information

Media	Pore Size (µm)	Product Code				
		Packaging 50/pk	Packaging 200/pk	Packaging 500/pk	Packaging 1000/pk	Packaging 5000/pk
Cellulose Acetate (CA)	0.22	1225617	1225618	1225619	1233871	
Cellulose Acetate (CA)	0.45	1225620	1225622	1225623	1233882	
Nylon (NY)	0.22	1224746	1224747	1224748	1229460	1224749
Nylon (NY)	0.45	1224753	1224754	1224755	1229462	1224756
Nylon (NY)	1.2	1224760	1224761			
Nylon (NY)	5.0	1224763	1224764	1224765	1229464	
Polyethersulfone (PES)	0.22	1233547			1233544	3049950
Polyethersulfone (PES)	0.45	1233548			1233545	3019423
Polypropylene (PP)	0.22	1224808	1224809	1224810	1229452	1225602
Polypropylene (PP)	0.45	1224811	1224812	1224813	1229454	1225607
Polytetrafluoroethylene (PTFE)	0.22	1224780	1224781	1224782	1229447	1224783
Polytetrafluoroethylene (PTFE)	0.45	1224787	1224788	1224789	1229449	1224790
Polyvinylidene Fluoride (PVDF)	0.22				3049952	
Polyvinylidene Fluoride (PVDF)	0.45	3023135		3023188	3023187	
Glass Fiber/Nylon (GF/NY)	0.22	1224766	1224767	1224768	1229477	1224769
Glass Fiber/Nylon (GF/NY)	0.45	1224773	1224774	1224775	1229479	1224776
Glass Fiber/Polypropylene (GF/PP)	0.22	1224814	1224815		1229473	
Glass Fiber/Polypropylene (GF/PP)	0.45	1224817	1224818	1229481	1229475	
Glass Fiber/PTFE	0.22	1224794	1224795	1224796	1229469	1224797
Glass Fiber/PTFE	0.45	1224801	1224802	1224803	1229471	1224804

MICROFILTRATION

1.1.3 - 25 mm Syringe Filters ABLUO®



Features and Benefits

Housing materials: Acrylic and Polypropylene

Membrane materials: Cellulose Acetate, Nylon, PTFE, PES, PVDF, Regenerated Cellulose, Nitrocellulose, Polyethylene, Glass Fiber

Membrane Diameter: 25 mm

Pore Size: 0.20 µm, 0.22 µm, 0.45 µm, 0.80 µm, 1.20 µm, 5.0 µm and others.

Housing Diameter: 33 mm

Effective Filtration Area: 4.6 cm²

Pressure Bar: 5

Sterile: No

Inlet / outlet: FLL/MLL-MLS

Applications

- HPLC sample preparation
- Biological fluids
- Buffer solutions
- Sterile filtering of tissue culture media
- Protein aqueous solutions

Specification

- Increased throughput and speed of sample preparation.
- Lower hold-up volume due to an improved flow channel design and reduced spacing between the supports within the housing for better handling of small sample volumes or costly samples.
- Increased operating pressure up to 75 psi due to the over-mold that prevents sample leaking at the seam and keeps the filter unit from bursting in half.
- Strict quality control: Syringe filters are integrity tested to ensure a proper filter fit and weld to eliminate any potential filter by-pass.
- Accurate labeling: Each filter is labeled with the specific filter material and pore size for easy identification even if the syringe filter is not in its original packaging.
- Modified acrylic housing to bidirectionally support the membrane allowing sample injection or aspiration.

Ordering information

Membrane Material	Pore Size (µm)	Housing Material	Color	Product Code
				Packaging 500/pkg
Cellulose Acetate (CA)	0.22	Acrylic	Blue	FJ25ANCCA002DD01
Cellulose Acetate (CA)	0.45	Acrylic	Yellow	FJ25ANCCA004FD01
Cellulose Acetate (CA)	0.80	Acrylic	Green	FJ25ANCCA008ED01
Cellulose Acetate (CA)	1.20	Acrylic	Red	FJ25ANCCA012CD01
Cellulose Acetate (CA)	5.0	Acrylic	Brown	FJ25ANCCA050PD01
Nylon (NY)	0.20	Polypropylene	Transparent	FJ25BNPNY002AD01
Nylon (NY)	0.45	Polypropylene	Transparent	FJ25BNPNY004AD01
Polyethersulfone (PES)	0.22	Polypropylene	Transparent	FJ25BNPPS002AD01
Polyethersulfone (PES)	0.45	Polypropylene	Transparent	FJ25BNPPS004AD01
Polyethersulfone Asymmetric (PES)	0.22	Polypropylene	Transparent	FJ25BNPPX002AD01
Nitrocellulose Mixed Esters (MCE)	0.22	Acrylic	Transparent	FJ25BNCNC002AD01
Nitrocellulose Mixed Esters (MCE)	0.45	Acrylic	Transparent	FJ25BNCNC004AD01
Regenerated Cellulose (RC)	0.20	Polypropylene	Transparent	FJ25BNPRC002AD01
Regenerated Cellulose (RC)	0.45	Polypropylene	Transparent	FJ25BNPRC004AD01
Polyvinylidene Fluoride (PVDF)	0.22	Polypropylene	Transparent	FJ25BNPPV002AD01
Polyvinylidene Fluoride (PVDF)	0.45	Polypropylene	Transparent	FJ25BNPPV004AD01
Polytetrafluoroethylene (PTFE)	0.20	Polypropylene	Transparent	FJ25BNPPT002AD01
Polytetrafluoroethylene (PTFE)	0.45	Polypropylene	Transparent	FJ25BNPPT004AD01
Polytetrafluoroethylene Hydrophilic (PTFE HP)	0.22	Polypropylene	Transparent	FJ25BNPPH002AD01
Polyethylene (PE)	0.20	Polypropylene	Transparent	FJ25BNPPE002AD01
Polyethylene (PE)	0.50	Polypropylene	Transparent	FJ25BNPPE005AD01

1.1 Syringe-Filters

Membrane Material	Pore Size (μm)	Housing Material	Color	Product Code
				Packaging 500/pk
Glass Fiber (GF)	1.00	Polypropylene	Transparent	FJ25BNPGF010AD01
Glass Fiber (GF)	1.20	Polypropylene	Transparent	FJ25BNPGF012AD01
Glass Fiber (GF)	3.10	Polypropylene	Transparent	FJ25BNPGF031AD01
Glass Fiber / Cellulose Acetate (GF/CA)	GF 1.00 + CA 0.45	Polypropylene	Transparent	FJ25BNPGA004AD01
Glass Fiber / Nylon (GF/NY)	GF 1.00 + NY 0.45	Polypropylene	Transparent	FJ25BNPGN004AD01
Glass Fiber / Polytetrafluoroethylene (GF/ PTFE)	GF 1.00 + PTFE 0.45	Polypropylene	Transparent	FJ25BNPGT004AD01
Glass Fiber / Polyethylene (GF / PE)	GF 1.00 + PE 0.45	Polypropylene	Transparent	FJ25BNPGE005AD01
Glass Fiber / Polyvinylidene Fluoride (GF / PVDF)	GF 1.00 + PVDF 0.45	Polypropylene	Transparent	FJ25BNPGV004AD01
Glass Fiber / Regenerated Cellulose (GF / RC)	GF 1.00 + RC 0.45	Polypropylene	Transparent	FJ25BNPGC004AD01

25 mm Syringe Filters - Sterile ABLUO®



Features and Benefits

- Housing Material:** Acrylic
- Membrane Material:** Cellulose Acetate, Polyethersulfone, Polyvinylidene Fluoride
- Membrane Diameter (mm):** 25
- Housing Diameter (mm):** 33
- Effective Filtration Area:** 4.6 cm²
- Sterile:** Yes
- Inlet / outlet:** FLL/MLL-MLS

Applications

- Filtration of Aqueous and Alcohol Solutions
- Sterile Filtration and Clarification
- Cell Culture
- Analytical Sample Prep, uHPLC
- IC Chromatography
- Clarification
- Protein Chemistry
- Filtration of Aqueous and Organic Solutions

Ordering information

Membrane Material	Pore Size (μm)	Color	Product Code
			Packaging 50/pk
Cellulose Acetate (CA)	0.22	Blue	FJ25ASCCA002DL01
Cellulose Acetate (CA)	0.45	Yellow	FJ25ASCCA004FL01
Cellulose Acetate (CA)	0.80	Green	FJ25ASCCA008EL01
Cellulose Acetate (CA)	1.20	Red	FJ25ASCCA012CL01
Cellulose Acetate (CA)	5.0	Brown	FJ25ASCCA050PL01
Polyethersulfone (PES)	0.22	Transparent	FJ25BSCPS002AL01
Polyethersulfone (PES)	0.45	Transparent	FJ25BSCPS004AL01
Polyvinylidene Fluoride (PVDF)	0.22	Transparent	FJ25BSCPV002AL01
Polyvinylidene Fluoride (PVDF)	0.45	Transparent	FJ25BSCPV004AL01

MICROFILTRATION

25mm Syringe Filters- Sterile CAMEO®



Membrane diameter: 25mm
Holdup volume: < 60 microliter
Sample volume: < 100 mL without prefilter; < 150 mL with prefilter
Operating parameters: Maximum Operating Temperature 55°C/131°F, Maximum Operating Pressure 75psi (5.2 bar)
Inlet / outlet: FLL-MLS

Description and Use

GVS Life Sciences offers a comprehensive range of disposable syringe filter devices designed to provide fast and efficient filtration of aqueous and organic solutions. They are made with a wide variety of membrane filters. These syringe filters are suitable for numerous applications in pharmaceutical, environmental, biotechnology, food/beverage, and agricultural testing laboratories.

Features

Housing: Acrylic heat-sealed without the use of glues or sealants

Media: nylon, PES, glass fiber, and cellulose acetate

Filtration area: 4.1 cm²

Autoclavable: No

Prefilter: 1.0 micron binderless fiberglass

Sterile: Gamma irradiated

Housing diameter: 29mm

Benefits

- Increased throughput and speed of sample preparation
- Lower hold-up volume due to an improved flow channel design and reduced spacing between the supports within the housing for better handling of small sample volumes or costly samples.
- Increased operating pressure up to 75 psi due to the over-mold that prevents sample leaking at the seam and keeps the filter unit from bursting in half.
- Strict quality control: syringe filters are integrity-tested to ensure a proper filter fit and weld to eliminate any potential filter by-pass.
- Accurate labeling: Each filter is labeled with the specific filter material and pore size for easy identification even if the syringe filter is not in its original packaging.
- Modified acrylic housing to bidirectionally support the membrane allowing sample injection or aspiration.

Applications

- HPLC
- Biological fluids
- Buffer solutions
- Sterile filtering of tissue culture media
- Proteinaceous solutions

Ordering information

Media	Pore Size (µm)	Product Code			
		Packaging 50/pk	Packaging 200/pk	Packaging 500/pk	Packaging 1000/pk
Cellulose Acetate (CA)	0.22	1213657	1213678	1213667	1224304
Cellulose Acetate (CA)	0.45	1213723	1213739	1213731	1224305
Cellulose Acetate (CA)	0.8	1213758		1213767	1262953
Glass Fiber (GF)	1.00	1212701			
Nylon (NY)	0.1	1213699	1213702	1213701	
Nylon (NY)	0.22	1213708	1213711	1213709	1215268
Nylon (NY)	0.45	1213724	1213726	1213725	1221513
Nylon (NY)	1.2	1213737	1213740	1213738	1239715
Nylon (NY)	5.0	1213744	1213746	1213745	1239369
Nitrocellulose Mixed Ester (MCE)	0.22	1213785	1213806	1213795	1224307
Nitrocellulose Mixed Ester (MCE)	0.45	1213814	1213828	1213822	1224306
Polyethersulfone (PES)	0.22	1214220	1222370	1222369	1215295
Polyethersulfone (PES)	0.45	1215130	1222266	3012605	1223379
Polyethersulfone (PES)	0.8	1222374		1222375	
Glass Fiber / Cellulose Acetate	0.22	1214344	1214349	1214346	1221496
Glass Fiber / Cellulose Acetate	0.45	1214358	1214360	1214359	1215245
Glass Fiber / Cellulose Acetate	0.8	1212630	1213122	1222378	3038584
Glass Fiber / Polyethersulfone	0.22	1222388	1215190	1214904	1215294
Glass Fiber / Polyethersulfone	0.45	1215146			1224085
Glass Fiber / Polyethersulfone	0.8	1222395			

1.1 Syringe-Filters

1.1.4 - 30 mm Syringe Filters - CAMEO®



Description and use

GVS Life Sciences offers a comprehensive range of disposable syringe filter devices designed to provide fast and efficient filtration of aqueous and organic solutions. They are made with a wide variety of membrane filters with a polypropylene housing. These syringe filters are suitable for numerous applications in pharmaceutical, environmental, biotechnology, food/ beverage, and agricultural testing laboratories.

Features

Housing: Pure polypropylene is heat-sealed without the use of glues or sealants

Media: Nylon, PTFE, PES, PVDF, glass fiber, cellulose acetate and polypropylene

Autoclavable: Yes

Prefilter: 1.0 micron binderless glass-fiber

Filtration area: 4.8 cm² (0.74 inch²)

Housing diameter: 33mm

Membrane diameter: 30mm

Holdup volume: < 60 microliter

Sample volume: < 120 mL without prefilter

Inlet / outlet: FLL-MLS

Benefits

- Increased throughput and speed of sample preparation.
- Lower hold-up volume due to an improved flow channel design and reduced spacing between the supports within the housing for better handling of small sample volumes or costly samples.
- Increased operating pressure up to 130 psi due to the over-mold that prevents sample leaking at the seam and keeps the filter unit from bursting in half.
- Strict quality control: Syringe filters are integrity tested to ensure a proper filter fit and weld to eliminate any potential filter by-pass.
- Accurate labeling: Each filter is labeled with the specific filter material and pore size for easy identification even if the syringe filter is not in its original packaging.

Applications

- HPLC sample preparation
- Dissolution testing
- Content uniformity
- Environmental samples
- Composite assays
- Food analysis
- Biofuel analysis

Ordering information

Material w/o Pre-filter	Pore Size (µm)	Product Code				
		Packaging 50/pk	Packaging 200/pk	Packaging 500/pk	Packaging 1000/pk	Packaging 5000/pk
Cellulose Acetate (CA)	0.22	1213641	1213192	1214014	1229443	1270152
Cellulose Acetate (CA)	0.45	1214778	1214932	1214966	1229444	1227275
Cellulose Acetate (CA)	0.8	1226939	1226941	1226940	1229445	
Glass Fiber (GF)	1.00	1227204		1227205	1229451	1226150
Nylon (NY)	0.1	1224100	1224101	1224103	1229459	
Nylon (NY)	0.22	1224104	1224105	1224106	1229461	1224110
Nylon (NY)	0.45	1224112	1224113	1224114	1226917	1224117
Nylon (NY)	1.2	1224119	1224120	1224121	1229463	
Nylon (NY)	5.0	1224124	1224125	1224126	1229465	3019424
Polyethersulfone (PES)	0.22	1233549			1233541	
Polyethersulfone (PES)	0.45	1233550		1233551	1233543	
Polypropylene (PP)	0.22	1224172	1224173	1224174	1229453	1237235
Polypropylene (PP)	0.45	1224310	1224311	1224312	1229458	
Polytetrafluoroethylene (PTFE)	0.22	1224143	1224144	1224145	1229448	1224146
Polytetrafluoroethylene (PTFE)	0.45	1224150	1224151	1237721	1229450	1224153
Polyvinylidene Fluoride (PVDF)	0.22	3038551			3038552	
Polyvinylidene Fluoride (PVDF)	0.45	3020528		3020351	3023084	3049953
Glass Fiber/Cellulose Acetate	0.22	1226942	1226944	1226943	1229466	3050049
Glass Fiber/Cellulose Acetate	0.45	1226945	1226947	1226946	1229467	1270153
Glass Fiber/Cellulose Acetate	0.8		1226951	1226950		
Glass Fiber/Nylon	0.1				1229480	
Glass Fiber/Nylon	0.22	1224127	1224128	1224129	1229478	1224133
Glass Fiber/Nylon	0.45	1224135	1224136	1224137	1226916	1224138
Glass Fiber/Polyethersulfone	0.45	3050121		3050122		
Glass Fiber/Polypropylene	0.22	1224175	1224176	1224177	1229474	
Glass Fiber/Polypropylene	0.45	1224313	1224314	1224315	1229476	
Glass Fiber/PTFE	0.22	1224157	1224158	1224159	1229470	1224160
Glass Fiber/PTFE	0.45	1224164	1224165	1224166	1229472	1224167

MICROFILTRATION

1.2 Vent Filters

40 mm Bacterial and Viral Vent Filter

**Membrane:**

Glass Fiber

Retention:99.999 % Bacterial and 99.9999 % Viral filtration efficiency of 0,027 μm phage**Housing:**

Polypropylene

Connectors:

6 to 8 mm stepped barb

Filter Area:14,6 cm^2 **Air Flow Rate:**

21 L/min at 3 psi

Housing Diameter:

53 mm

Housing Length:

59 mm

Maximum Pressure:

2,7 bar (approx. 40 psi)

Autoclavable up to 10 times

Applications

- Sterile venting of filling vessels and carboys
- Autoclave venting
- Low volume sterile filtration of non-aqueous fluids
- In-line sterilization of and particulate removal from air and gases

Box Dimension: 29 x 38 x 30 cm - **Box Weight:** 1,58 Kg - **Quantity / Box:** 50 Units

Ordering information

Product Code	Description
VF40CBPGF000AL01	Vent filter Ø40mm AS connectors PP transparent membr. GF - 50 pieces
VF40CSPGF000AX01	Vent filter Ø40mm AS connectors PP transparent membr. GF - sterile 10 pieces

50 mm Vent Filter

**Membrane:**

hydrophobic PTFE reinforced with polypropylene

Porosities:0,45 μm or 0,20 μm **Housing:**

Polypropylene

Connectors:

6 to 12 mm stepped barb

Filter Area:19,6 cm^2 **Air Flow Rate:**32 L/min at 1 bar (0,45 μm)27 L/min at 1 bar (0,20 μm)**Housing Diameter:**

63 mm

Housing Length:

53 mm

Maximum Pressure:

3,5 bar (approx. 50 psi)

Sterilization:

Autoclave at 121 C or ETO

Autoclave up to 10 times

Applications

- Sterile venting of filling vessels and carboys
- Autoclave venting
- Low volume sterile filtration of non-aqueous fluids
- In-line sterilization of and particulate removal from air and gases

Box Dimension: 30 x 30 x 35 cm - **Box Weight:** 3,6 Kg - **Quantity / Box:** 100 Units
30 x 8,5 x 14 cm 0,45 Kg 10 Units

Ordering information

Product Code	Description
VF50ANPPT002AC01	Vent filter Ø50mm barb connectors PP transparent membr. PTFE 0.20 μm - 100 pieces
VF50ANPPT004AC01	Vent filter Ø50mm barb connectors PP transparent membr. PTFE 0.45 μm - 100 pieces
VF50ASPPT002AX01	Vent filter Ø50mm barb connectors PP transparent membr. PTFE 0.20 μm - sterile 10 pieces
VF50ASPPT004AX01	Vent filter Ø50mm barb connectors PP transparent membr. PTFE 0.45 μm - sterile 10 pieces

MICROFILTRATION

50 mm PTFE Vent Filter



Membrane:
PTFE
Housing:
Polypropylene
Filter Area:
1735 mm²
Air Flow Rate:
27L and 32L /min@1bar
Outer Diameter:
60mm
Overall Length:
63mm
Operating Pressure:
4 bar

Premier Performance

- PTFE membrane and polypropylene housing have broad chemical resistance
- Large surface area provides greater throughput and high air flow rate
- Reusable and autoclavable
- Disposable
- Hose barb connection inlet /outlet (6-12mm)

Description and Use

GVS Life Sciences 50 mm hydrophobic PTFE membrane filters are ideal for sterilizing gases, venting and degassing sterile containers, and sterilizing or clarifying organic solutions.

Features & Benefits

- Hydrophobic
- Chemically Compatible
- Lot-to-lot Consistency

Applications

- Sterile venting of filling vessels and carboys
- Autoclave venting
- Low volume sterile filtration of non-aqueous fluids
- In-line sterilization of and particulate removal from air and gases

Ordering information

Product Code	Description
95000	0.22um pore size, 50/pk
95001	0.45um pore size, 50/pk

MICROFILTRATION

1.3 Calyx Capsule Filters



Description and Use

GVS Life Sciences capsules are disposable filtration units designed for the removal of particles or bacteria from aqueous or solvent solutions and gas streams. They are ready to use, eliminating the need to disassemble, clean and reassemble filter housings. GVS Life Sciences capsules contain no glue or surfactants and feature serial layer filter design for increased throughput and extended life. Two upstream vents are included to facilitate venting in any position. All capsules containing membrane media are preflushed with purified water to reduce extractables. GVS Life Sciences capsules pass class VI toxicology testing and are integrity tested prior to shipment.

Capsule filters are available in sterile and non-sterile versions.

The capsules are available with the following connections: 3/8 or 1/4 inch hose barb, 1/4 or 1/2 inch NPTM, 1.5 inch sanitary flange.

Table 1: Dimensions

Diameter: 3.5" (9 cm)		
Capsule Size	Effective Filtration Area	Length ¹
Small	0.8 ft ² (748 cm ²)	3.5 - 4.7" (9 - 13 cm)
Medium	3.0 ft ² (2806 cm ²)	7.6 - 8.8" (19 - 23 cm)
Large	5.9 ft ² (5500 cm ²)	11.5 - 12.7" (29 - 33cm)

¹Varies with connection style

Table 2: Operational Limits

Maximum Operational Pressure	80psi (5.5bar) @ 70°F (21°C) in Liquid 50psi (3.8bar) @ 70°F (21°C) in Gas
Maximum Differential Pressure	60psi (4.1bar) @70°F (21°C)
Maximum Operating Temperature	110°F (43°C) @ ≤ 30 psi (2.1bar) Operating Pressure

Table 3: Integrity Testing

Water intrusion is ≤ 1.8cc/min per ft ² of Effective Filtration Area	
Micron Rating	Test Pressure
0.1 μm	40psi (2.8bar)
0.2 μm	30psi (2.1bar)

Ordering information: PES Capsule Filters, Hydrophilic, Polyester housing

Pore Size μm	Filtration area ft ² (cm ²)	Length in (cms)	Adaptors: L= ¼ - ½ in stepped hose barb; M= 3/8 in hose barb; R= ¼ in NPT male; Y= 1.5 in sanitary flange; W= ½ in NPT male				
			LL	MM	RR	WW	YY
0.1	0.8 (748)	3,5 (9)			1222323		
		3,5 (9)			1213608 1214001*		
0.2	0.8 (748)	4,0 (10)	1214225*	1223129			
		4,3 (11)				1225346	
		4,7 (12)					1213956
0.2	3.0 (2808)	8,1 (20)	1214436*				
		8,4 (21)				1215154 1223845*	
		8,8 (22)					1222327
0.2	5.9 (5500)	12,7 (32)					1214083
0.4	0.8 (748)	3,5 (9)			1213610		
		4,0 (10)	1214227*				
0.4	3.0 (2808)	8,1 (20)	1222432*				

*sterile product

Ordering information: PTFE Capsule Filters, Hydrophobic, Polypropylene housing

Pore Size μm	Filtration area ft ² (cm ²)	Length in (cms)	Adaptors: L= ¼ - ½ in stepped hose barb; M= 3/8 in hose barb; R= ¼ in NPT male; Y= 1.5 in sanitary flange; W= ½ in NPT male			
			MM	RR	WW	YY
0.1	0.8 (748)	3,5 (9)		1213160		
		4,0 (10)	1213154			
		4,3 (11)			1212936	
0.1	5.9 (5500)	11,5 (29)		1212982		
0.2	0.8 (748)	3,5 (9)		1213158		
		4,0 (10)	1213155			
		4,3 (11)			1212937	
		4,7 (12)				1212978

MICROFILTRATION

Pore Size μm	Filtration area ft^2 (cm^2)	Length in (cms)	Adaptors: L= $\frac{1}{4}$ - $\frac{1}{2}$ in stepped hose barb; M= $\frac{3}{8}$ in hose barb; R= $\frac{1}{4}$ in NPT male; Y= 1.5 in sanitary flange; W= $\frac{1}{2}$ in NPT male			
			MM	RR	WW	YY
0.2	5.9 (5500)	11,5 (29)		1212987		
0.4	0.8 (748)	3,5 (9)		1213161		
0.4	5.9 (5500)	11,5 (29)		1212992		

Ordering information: PP Capsule Filters, Hydrophobic, Polypropylene housing

Pore Size μm	Filtration area ft^2 (cm^2)	Length in (cms)	Adaptors: L= $\frac{1}{4}$ - $\frac{1}{2}$ in stepped hose barb; M= $\frac{3}{8}$ in hose barb; R= $\frac{1}{4}$ in NPT male; Y= 1.5 in sanitary flange; W= $\frac{1}{2}$ in NPT male				
			LL	MM	RR	WW	YY
0.22	0.8 (748)	3,5 (9)			1213584 1213941		
		4,0 (10)	1212949	1213578			
		4,3 (11)				1212929	
		4,7 (12)					1212971
0.22	3.0 (2808)	8,4 (21)				1213057	
		8,8 (22)					1213058
0.22	5.9 (5500)	11,5 (29)			1213089		
		12,3 (31)				1213090 1213893	
		12,7 (32)					1213091
0.45	0.8 (748)	3,5 (9)			1213596		
		4,0 (10)	1212950	1213591			
		4,3 (11)				1212930	
		4,7 (12)					1212972
0.45	5.9 (5500)	11,5 (29)			1213095		
1.2	0.8 (748)	3,5 (9)			1213611		
		4,0 (10)	1212951	1213605			
		4,3 (11)				1212932	
1.2	5.9 (5500)	11,5 (29)			1213100		
		12,0 (30)		1213098			
		12,3 (31)				1213101	
3.0	0.8 (748)	3,5 (9)			1213157		
		4,3 (11)				1212933	
3.0	3.0 (2808)	8,1 (20)		1213070			
5.0	0.8 (748)	3,5 (9)			1213620		
		4,0 (10)	1212953	1213619			
		4,7 (12)					1212975
5.0	3.0 (2808)	7,6 (20)			1213078		
		8,8 (22)					1213080
5.0	5.9 (5500)	11,5 (29)			1213111		
		12,7 (32)					1213113
10.0	0.8 (748)	3,5 (9)			1213622		
		4,0 (10)	1212954	1213621			
		4,3 (11)				1212935	
10.0	3.0 (2808)	8,1 (20)		1213081			
10.0	5.9 (5500)	11,5 (29)			1213117		
25.0	0.8 (748)	3,5 (9)			1213617		
		4,0 (10)		1213616			
25.0	3.0 (2808)	7,6 (20)			1220684		
25.0	5.9 (5500)	12,3 (31)				1215179	

MICROFILTRATION

Ordering information: NYLON Capsule Filters, Hydrophilic, Polyester housing

Pore Size μm	Filtration area ft^2 (cm^2)	Length in (cms)	Adaptors: L= $\frac{1}{4}$ - $\frac{1}{2}$ in stepped hose barb; M= $\frac{3}{8}$ in hose barb; R= $\frac{1}{4}$ in NPT male; Y= 1.5 in sanitary flange; W= $\frac{1}{2}$ in NPT male				
			LL	MM	RR	WW	YY
0.1	0.8 (748)	3,5 (9)			1213540 3044268*		
		4,0 (10)	1212939	1213529			
0.1	3.0 (2080)	8,8 (22)					1221768
0.1	5.9 (5500)	11,5 (29)			1212899		
0.2	0.8 (748)	3,5 (9)			1213561 3031900*		
		4,0 (10)		1213550 1213757*			
0.2	5.9 (5500)	11,5 (29)			1212905		
0.4	0.8 (748)	3,5 (9)			1213577		
		4,0 (10)		1214457*			
0.4	5.9 (5500)	11,5 (29)			1212910		
		12,0 (30)		1212908			
		12,3 (31)				1212911	

*sterile product

1.4 Centrifuge Filters - Centrex

CENTREX centrifuge filters are supplied with a range of filtration and separation media. 1.5 ml and 5.0 ml sterile and nonsterile versions are available.



Features and Benefits

- Centrifugal filter units with various types of membrane filter
- Rapid and simple preparation of a large number of samples
- More than six samples can be processed at once
- Ideal for automated systems and high-speed batch filtration with robots
- Considerably reduced contamination risk when working with radioactive biologically hazardous material
- Cross contamination avoided

Applications

- 0.45 μm cellulose acetate membrane for the rapid elution of agarose gels
- Nylon and cellulose acetate membranes for the removal of particles and microorganisms from HPLC samples
- Sample preparation for quality control
- Cellulose acetate and nitrocellulose membrane for rapid clearing and filtration of aqueous solutions

Ordering information

Color	Pore size	Membrane	1.5ml ST 50/carton	1.5ml Non-ST 250/carton	5ml ST 50/carton	5ml Non-ST 250/carton
Brown	0.2 μm	Nylon	10467003		10467015	10467010
Tan	0.45 μm	Nylon	10467007	10467002	10467021	10467012
Blue	0.2 μm	Cellulose Acetate	10467004	10467009	10467013	
White	0.45 μm	Cellulose Acetate	10467006	10467011	10467017	
Green	0.8 μm	Cellulose Acetate	10467008			
Pink	0.2 μm	Nitrocellulose	10467001			
Rust	0.45 μm	Nitrocellulose	10467005		10467019	

MICROFILTRATION

1.5 Bottle-top Filters - ZapCap



Description and Use

For filtration of medium volumes, cell culture media and HPLC solutions.

Ordering information

Product Code	Description
10443401	ZapCap-S, cellulose acetate, 0.2 μ m, sterile, 12/pk, housing: Polystyrene
10443411	ZapCap-S, cellulose acetate, 0.45 μ m, sterile, 12/pk, housing: Polystyrene
10443421	ZapCap-CR, nylon, 0.2 μ m, 12/pk, housing: Polypropylene
10443423	ZapCap-CR, nylon, 0.45 μ m, 12/pk, housing: Polypropylene
10443425	ZapCap-CR, PTFE, 0.45 μ m, 12/pk, housing: Polypropylene
10443430	ZapCap-S PLUS, cellulose acetate w/ glass fiber prefilter, 0.2 μ m, sterile, 12/pk, housing: Polystyrene
10443435	ZapCap-S PLUS, cellulose acetate w/ glass fiber prefilter, 0.45 μ m, sterile, 12/pk, housing: Polystyrene

Features & Benefits

- Complete 500 ml units with tubing nozzle for attaching to bottles (bottle-top)
- Connection seals on any standard bottles 33 to 45 mm
- Membrane diameter 76 mm, filter area 39.2 cm²
- ZapCap-S with included borosilicate prefilter for high flow rates
- ZapCap-S Plus with integral borosilicate prefilter for very high flow rates
- ZapCap-CR, the chemical-resistant bottle-top filter
- Can be used up to 50°C

Typical Applications

ZapCap-S – Filtration of cell culture media

1. Cellulose acetate membrane filters (CA) with extremely low protein binding for cell culture media and other aqueous solutions
2. Sterile filtration of solutions which cannot be autoclaved

ZapCap-S Plus – Sterile filtration and clarification of difficult-to-filter aqueous solutions

ZapCap-CR – Filtration of HPLC solutions

1. Polyamide membrane filters (NYL) for the retention of fine particles and microorganisms in HPLC/FPLC solutions when the column packing is 10 μ m
2. PTFE membrane filters for the retention of particles in organic solutions; strong acids or aldehydes

1.6 Extractor - Ethidium bromide (EtBr) waste reduction system



One-step filtration polypropylene funnel device for the rapid removal of ethidium bromide from gel-staining solutions.

This disposable unit contains an activated carbon matrix, which removes > 99% of ethidium bromide from electrophoretic buffer quickly and easily. Each device can decontaminate up to 10 litres of gel-staining solution. After filtration, the decontaminated solution can be safely poured down the laboratory drain.


The extractor funnel device fits most standard laboratory flasks and bottles (neck size 33 to 45 mm), and the unit includes a cap for storage between uses. The polypropylene housing is chemically resistant to organics. Also included in the package are glass fiber prefilters, which remove gel pieces and other debris to avoid premature clogging of the carbon filter.

Ordering information

Product Code	Description
10448030	Ethidium Bromide Extractor Waste System 2/pk, Polypropylene
10448031	Ethidium Bromide Extractor Waste System 6/pk, Polypropylene

MICROFILTRATION

1.7 Filter Holders for Membranes

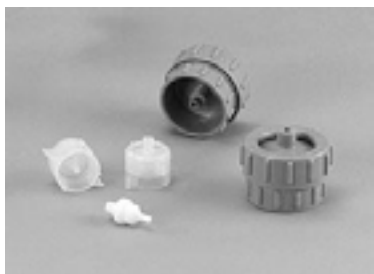

Description and Use

To insure precise filtration, GVS Life Sciences offers a selection of filtration holders and apparatus that are designed to work with GVS Life Sciences membranes and are built to exacting standards. In most applications, the filter holder is just as important as the filter for accurate results every time.


Filter holders are available for a wide variety of applications including air analysis, chemotaxis, tissue culturing and general aqueous and solvent filtration.

We also offer some unique products not found anywhere else like our filter static eliminator and our GVS Life Sciences analytical filter funnel that is so easy to use, you will wonder how you got along without one.

PRODUCT INFORMATION – PRODUCT SELECTION GUIDE

Specifications	13 mm	25 mm	47 mm
Materials	Celcon (acetal copolymer)	Polypropylene - body & support	Polypropylene - body & support
O-rings	PTFE	Silicone	Silicone
Filter Size	13 mm	25 mm	47 mm
Prefilter Cap size	10 mm	22 mm	42 mm
Filtration Area	0.8 cm ²	3.5 cm ²	13.5 cm ²
Diameter	16 mm (0.6 in)	30 mm (1.2 in)	63.5 mm (2.5 in)
Height	35 mm (1.4 in)	30 mm (1.2 in)	50 mm (2.0 in)
Maximum Liquid Temperature	80°C (176°F)	80°C (176°F)	80°C (176°F)
Maximum Operating Pressure	2.8 bar (40 psi)	2.9 bar (42 psi)	1.9 bar (71 psi)
Autoclaving	15 minutes at 121°C (250°F) and 15 psi	20 minutes at 121°C (250°F) and 15 psi	20 minutes at 121°C (250°F) and 15 psi
Connections, Inlet	Female Threaded Luer	Female LuerLok	1/4 inch NPTM, Female Luer Slip
Connections, Outlet	Male Luer Slip	Male Luer Slip	1/4 inch NPTM, Female Luer Slip
Shipping Weight	20 g (0.7 lb)	20 g (0.7 lb)	470 g (1.1 lb)

1.7.1 - 13 mm Filter Holder, Swinney

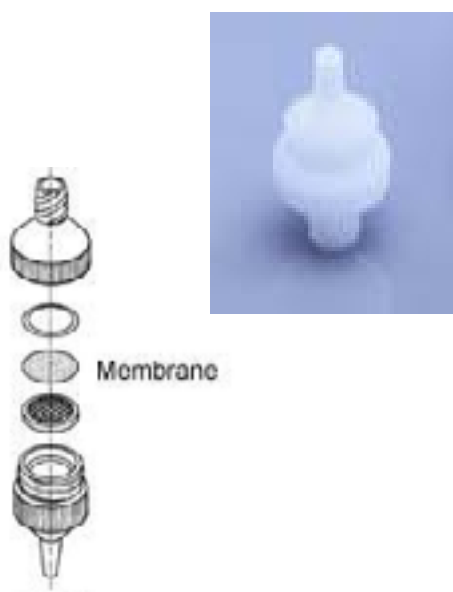

This GVS Life Sciences Swinney 13 mm filter holder is excellent for small volume (1-5 mL) particulate removal from fluids dispensed with a syringe. It is used in filtering biofluids, ophthalmics, gas chromatography samples, and lubricants that must be applied dust free to critical parts such as bearings. The holder is resistant to alcohols, esters, ethers, glycols, aromatic hydrocarbons, halogenated hydrocarbons, ketones, oils, photoresists and many other chemicals. Although suitable for most weak acids and bases, we recommend that you test for compatibility with acids.

Used in filtering

- Biofluids
- Ophthalmics
- Gas chromatography samples
- Lubricants

Ordering information

Product Code	Description	Qty/ pk
1220950	Filter Holder, Swinney, 13mm diameter	5/pk

MICROFILTRATION

1.7.2 - 25 mm Filter Holder, Polypropylene



The GVS Life Sciences polypropylene 25 mm filter holders are very useful for ultra cleaning and sterilizing small volumes of liquids from a syringe. Due to the polypropylene construction they can be used over a wide temperature range with excellent chemical compatibility. In the case of the syringe, the inlet cap locks into the base to prevent twisting damage to the membrane as the cap is tightened. Projection lugs on the base and the cap allow these units to be assembled and sealed quickly and efficiently. Typically, the 25 mm is used to filter up to 50 mL of sample. With the syringe holder type, dual support screens prevent membrane rupture in case back pressure is applied. It also allows for bi-directional sample flow. The polypropylene holder has a broad chemical compatibility range. It can withstand temperatures up to 121°C. and be autoclaved.

Features & Benefits

- Excellent chemical compatibility
- Quick, efficient assembly
- No need for special tools
- Excellent temperature and chemical resistance
- Several syringe holders can be attached together for serial filtration


Applications

- Point of use sampling
- Particulate removal
- Used in filtering chromatography solvents
- General filtration


Ordering information

Product Code	Description	Qty/ pk
1214250	Filter Holder Polypropylene: 25 mm diameter	10/pk
1214526	Filter Holder Polypropylene Support Screen: 25 mm	10/pk

1.7.3 - 47 mm Filter Holder, Polypropylene

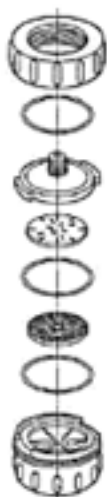

Polypropylene In-Line

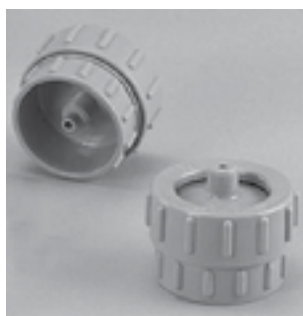

The GVS Life Sciences polypropylene 47 mm filter holder is designed especially for ultra cleaning and sterilizing liquids under positive pressure. In addition this holder can be used for aseptic sampling of liquids or gases at point-of-use or when samples must be collected and processed on-site.

The polypropylene material allows these holders to be used over a wide temperature range with excellent chemical compatibility. Sealing is achieved by simple hand tightening of the locking ring. The 47 mm In-Line holder has dual support screens, which allow for flow in either direction. The inlet cap design and exterior locking ring allow the unit to be assembled quickly and efficiently without tearing the membrane. 3 O-rings help to prevent leaks with all membranes. The 47 mm can filter up to 1 liter depending upon the viscosity of the sample. The polypropylene holder can withstand temperatures up to 121°C and be autoclaved.

Features & Benefits

- Easy to use - unique lock ring design assures proper sealing without damage to the membrane
- Easy to clean
- Conforms with EPA Method 1311 for Toxicity Characteristic Leaching Procedure, 40 CFR, Part 261, 1991 Hazardous Waste Compliance Guide

Applications

- Point of use sampling
- Particulate removal
- Used in filtering chromatography solvents
- General filtration

Ordering information

Product Code	Description	Qty/ pk
1262579	Filter Holder Polypropylene: 47 mm	1/pk
1214260	Filter Holder Polypropylene: 47 mm	10/pk

MICROFILTRATION

1.7.4 - Gravi-Seal Analytical 47 mm Filter Holder



The GVS Life Sciences autoclavable filter holder combines a number of key features and benefits, making it a tremendous value. To begin with, the funnel has only two pieces. There are no clamps or locking devices to manipulate. A unique gravity sealing design allows for one-handed operation with no danger of filter by-pass or sample leakage when using depth filters. And it is stable and very solid with no costly replacement parts. It all adds up to the easiest and most cost-efficient analytical funnel available. GVS Life Sciences analytical funnels are available in polysulfone for aqueous samples. The polysulfone unit is autoclavable and chemically resistant for cell culture and microbiological applications. There are graduations up to 350 mL with 50 mL intervals. The No. 8 stopper mounts in a standard 1-liter filtering flask for individual tests or in three- and six-place stainless steel manifolds for multiple tests to run concurrently.

Features & Benefits

- Durable - break resistant, no extra parts to break or wear out
- Uses a 47 mm depth filter disc
- One-handed operation
- Only two parts
- No clamps, wheel locks, or magnets to wear out
- Solid, stable and easy to use

Applications

- Filtering liquids for sterility
- Particle removal
- General filtration
- Autoclavable

Ordering information

Product Code	Description	Qty/ pk
1213865	Gravi-Seal PES Analytical Filter Holder (complete unit): 47 mm	1/pk
1214124	Gravi-Seal PES Analytical Filter Holder (complete unit): 47 mm	3/pk
1213883	Gravi-Seal PES Analytical Filter Holder, Base Only	1/pk
1213882	Gravi-Seal PES Analytical Filter Holder, Funnel Only	1/pk
1214301	Funnel Extenders, Vaccum, PS	2/pk

MICROFILTRATION



1.8 SEPARA - Mini Vial



Save time and money in sample prep process with GVS Life Sciences filter and vial disposable device.

The single step filtering process is efficient and saves time.

Easy to press, fast and simple to use.

Features & Benefits

- Rapid sample preparation.
- Filtering with a plunger in the vial is a rapid single step, reducing sample loss.
- After filtration, the sample is ready for use in auto-sampling robot
- The pre-slit cap ensures easy and clean transfer of sample.
- Compatible with most standard auto-Samplers

Applications by filtration membrane

PTFE

- Filtration of strong acids and aggressive solutions

Nylon

- Clarification of aqueous and organic solvent solutions
- HPLC sample preparation
- Chromatography

Membrane RC

- Filtration of Aqueous and Organic Solutions
- Analytical Sample Prep, uHPLC
- Chromatography
- Clarification
- Protein Chemistry

Specifications

- Dimensions: 12mm diameter x 33mm height
- Material: Polypropylene, Septa, PTFE and silicone
- Filtering capacity: 0.48mL.
- The force needed to compress the unit is estimated at 8 psi (0.6 bar)
- Maximum operating temperature: 120° F (50° C)



Fig. 1 sample filling



Fig. 2 press down to filter sample



Fig. 3 filtered sample ready for analysis

Ordering information

Pore size	Membrane	Color	100/pk	1000/pk
0.45 μm	Nylon	Blue	MV32ANPNY004UC01	MV32ANPNY004UM01
0.20 μm	Nylon	Light Blue	MV32ANPNY002BC01	MV32ANPNY002BM01
0.45 μm	PTFE	Red	MV32ANPPT004CC01	MV32ANPPT004CM01
0.20 μm	PTFE	Pink	MV32ANPPT002TC01	MV32ANPPT002TM01
0.45 μm	RC	Black	MV32ANPRC004LC01	MV32ANPRC004LM01
0.20 μm	RC	Gray	MV32ANPRC002GC01	MV32ANPRC002GM01
0.45 μm	PVDF	Orange	MV32ANPPV004IC01	MV32ANPPV004IM01
0.20 μm	PVDF	Yellow	MV32ANPPV002FC01	MV32ANPPV002FM01
0.45 μm	PES	Dark Green	MV32ANPPS004WC01	MV32ANPPS004WM01
0.20 μm	PES	Light Green	MV32ANPPS002EC01	MV32ANPPS002EM01



MICROBIOLOGY

2. MICROBIOLOGY

2.1 Microbiological monitors & Analytical Funnels



GVS Life Sciences microbiological monitors provide a complete system solution for liquid sample preparation. Each monitor is a single-use, pre-sterilized filtering unit consisting of a measured filter funnel, base, pad, membrane, removable lid and plug. With no need for flaming or sterilization, GVS Life Sciences monitors cut multiple steps out of the testing process, saving your lab up to 70% in preparation time. These ready-to-use 100 ml units are suited for monitoring contaminants in all types of samples, from raw materials to finished products.

Features and Benefits:

- **Rapid testing:** With no need to sterilize funnels or filter base between samples, testing time can be reduced by up to 70%
- **No flaming required:** Combined filtration unit minimizes the risk of cross-contamination
- **All-in-one system:** Filtration unit easily converts to a Petri dish, which can be labeled and incubated for culturing
- **Reduced contamination:** Single-use materials virtually eliminate cross-contamination between funnel and membrane
- **Reproducible results:** All-in-one filtration unit reduces the chance of external error
- **Easy handling:** Ready-to-use, pre-sterilized monitors are simple to use

Analytical Funnel: The membrane can be removed and transferred to a Petri dish.
Microbiological Monitor: The membrane is fixed to the unit.

Applications:

Microbiological analysis of:

- Potable water
- Beer and wine
- Waste water
- Dairy products
- Soft drinks

Ordering information: Microbiological monitors sterile

Product Code		Description	Qty/ pkg
47 mm	56 mm		
10 497 511	10 497 603	Monitor, Nitrocellulose, 0.2 μ m, white/black grid	50
10 497 500	10 497 600	Monitor, Nitrocellulose, 0.45 μ m, white/black grid	50
10 497 501	n/a	Monitor, Nitrocellulose, 0.45 μ m, white/black grid, packaged ¹	50
10 497 502	10 497 601	Monitor, Nitrocellulose 0.45 μ m, black/white grid	50
10 497 503	10 497 602	Monitor, Nitrocellulose, 0.8 μ m, black/white grid	50

¹ Individually wrapped, sterile

Ordering information: Analytical Funnel

Analytical Funnels 100 ml (removable membrane) – 47 mm, sterile

Product Code	Description	Qty/ pkg
10497504	Funnel, Nitrocellulose, White/Black Grid Sterile 0.45 μ m	50
10497506	Funnel, Nitrocellulose, White/Black Grid Sterile 0.45 μ m, packaged ¹	50
10497507	Funnel, Nitrocellulose, White/Black Grid Sterile 0.2 μ m	50
10497508	Funnel, Nitrocellulose, Black/White Grid Sterile 0.45 μ m	50
10497509	Funnel, Nitrocellulose, Black/White Grid Sterile 0.45 μ m, packaged ¹	50
10497510	Funnel, Nitrocellulose, White/Black Grid Sterile 0.2 μ m, packaged ¹	50

¹ Individually bagged

Ordering information: Manifolds

Product Code	Description	Qty/ pkg
10498761	3-place vacuum manifold	1
10498762	6- place vacuum manifold	1

MICROBIOLOGY

2.2 Savur Analytical Filter Unit



GVS Life Sciences analytical filters are ready to use and available in presterilized individual bags. GVS Life Sciences analytical filter funnels with recoverable membranes are ideal for microbiological analysis and for quick, convenient filtration of aqueous or solvent solutions. Funnel capacity is 115 mL with a 47 mm diameter and 14.4 cm² filtration area.

Complete with low holdup polypropylene support pads, GVS Life Sciences funnels are ideal for quality control, microbiological, bioburden, and water/wastewater testing. Additional applications include patch testing of hydraulic and boiler fluids, sample clarification, particle counting, and gravimetric analysis. Each package of 12 units comes with an extender, which adapts the unit for side-arm flask filtrations. Optional No. 8 rubber stoppers may be ordered for filtering with flask.

The nitrocellulose mixed ester media is suitable for bacteria recovery and count and for water and wastewater testing. Thanks to the chemical compatibility, nylon will be dedicated to solvent clarification.

Polypropylene will be found for total digest of heavy metals.

Ordering information: Savur analytical Filter Unit, Sterile

Media	Pore Size (µm)	Gridded		Plain	
		47mm 12/pkg	47mm 200/pkg	47mm 12/pkg	47mm 200/pkg
Nitrocellulose	0.22		1214278	1214282	
Nylon	0.22			1213751	
Nitrocellulose	0.45	1214286	1214288	1214293	
Nylon	0.45	1213753	1214232	1214319	1229555
Polypropylene	25.0				1214661 1212668*

Product Code	Description	Qty/ pkg
1215262	Savur Funnel Extenders, PS	500/pkg

*non sterile product

2.3 Nutrient Liquid Media

2.3.1 - Liquid Media



2 ml ampouled media

Ready-to-use media considerably reduce the preparation time in quality control laboratories and also effectively reduce the risks of cross contamination. GVS Life Sciences is cooperating closely with quality assurance managers in the industry in the development of its own media and test kits.

This intensive product development has produced a range of products that is being used to monitor production plants and conduct microbiological checks on raw materials through to final product release in laboratories in more than 40 countries.

Features and benefits

- Wide range of products satisfies even special customer requirements
- Optimal media stability, sterility, and reproducibility
- Less time-consuming, higher productivity
- Batch-specific quality certificate in each pack

Applications

Microbiological analysis of:

- Drinking water
- Surface water
- Recreational water
- Purified water

Media variety and flexibility

Do you use ready-to-use media in 2 ml ampoules? Are you looking for nutrient media that are specifically suited to the cultivation of colonies after microfiltration? Would you perhaps like to fill your own agar plates? No matter what your specific needs, our extensive range of ready-to-use or prepared media products is available for just about any application.

Liquid media products come with a host of advantages, too. Aseptic preparation of nutrient media saves you valuable time and keeps costs to a minimum. All media undergo detailed quality control checks in accordance with recognized methods, guaranteeing uniform media preparation at all times. Lastly, our more comprehensive end product tests ensure optimal growth as well as stable and sterile media.

Media Descriptions

Brilliant Green Bile Broth 2%

BGGB contains two inhibitors of both gram-positive and selected gram-negative organisms, namely, oxgall and brilliant green dye. Fermentation is detected by gas production.

Cetrimide Broth

Pseudomonas aeruginosa is characterized by the production of pyocyanin (a blue green, water soluble, non-fluorescent, phenazine pigment) which is stimulated by the inclusion of magnesium chloride and potassium sulfate in the broth. Cetrimide (N-cetyl-NNN-trimethylammonium bromide) is added to inhibit bacteria other than *Pseudomonas aeruginosa*. Its action as a quaternary ammonium cationic detergent causes nitrogen and phosphorous to be released from bacterial cells other than *Pseudomonas aeruginosa*.

Coli Ceck with MDG

Used for presumptive identification of coliforms and the determination of the presence of *E. coli* in water samples by a presence/absence tech.

EC Broth

EC Broth contains casein peptone as a source of nutrients. Lactose provides the carbohydrate fermented by coliform bacteria and *Escherichia coli*. In addition, lactose-positive bacteria metabolize lactose with gas formation. Gram-positive bacteria are inhibited by the mixture of bile salts.

EC Broth with MUG

The presence of fluorescence using a long-wave UV light source confirms the presence of *Escherichia coli* and no further confirmation is required. MUG detects anaerogenic strains, which may not be detected in the conventional procedure. Lactose is a source of energy. Casein peptone provides additional nutrients. The mixture of bile salts is inhibiting for gram-positive bacteria, particularly bacilli and fecal streptococci. The substrate 4-methylumbelliferyl-b-D-glucuronide is hydrolyzed by an enzyme, β -glucuronidase, possessed by most *Escherichia coli* and a few strains of *Salmonella*, *Shigella* and *Yersinia*, to produce a fluorescent end product, 4-methylumbelliferone.



Brilliant Green Bile Broth



EC-Broth: Vial Left: Control; Vial Right: Broth inoculated with *Escherichia coli* ATCC 25922

MICROBIOLOGY

Enterococcus Broth

Enterococcus Broth is a modified version of the improved media described by Slanetz and Bartley with TTC. The membrane filtration method is simple to perform, does not require confirmation and permits a direct count of enterococci in 48 hours.

HPC Broth and HPC Broth with TTC HPC is used to determine total count at incubation temperatures of 35°C. All bacteria develop on HPC with indicator media and produce a red color as a result of the precipitation of formazan following the reduction of 2,3,5- triphenyltetrazolium chloride (TTC) by bacteria.

KF-Streptococcus Broth


KF-Streptococcus Broth is selective for the determination of fecal streptococci in polluted surface waters. Maltose and lactose are fermentable carbohydrates, sodium azide is the selective agent and brom cresol purple is the indicator dye.

Lauryl Sulfate or Lauryl Tryptose Broth

This media was developed for the detection of coliform organisms by the American Public Health Association (APHA). It is now the standard medium of choice in the presumptive phase of the standard coliform MPN test for the microbiological examination of water.

Mannitol Salt Broth

Because of the amount of peptones and beef extract, Mannitol Salt is a nutrient rich medium. Most bacteria (other than staphylococci) are inhibited by the high concentration of sodium chloride. Organisms capable of fermenting mannitol, e.g., *Staphylococcus aureus*, cause a pH change in the media. With phenol red as the pH indicator the colonies appear with a yellow coloration.

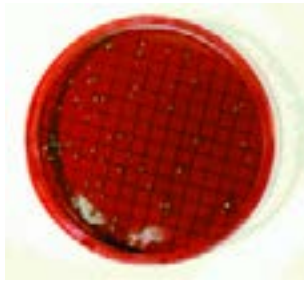

M-Endo Coliform Broth

Membrane Lauryl Sulfate Broth

This media was developed for the detection of coliform organisms and is now the media of choice for the enumeration of total coliforms and *Escherichia coli* in the United Kingdom. This media replaced membrane enriched broth containing 0.4% Teepol 610.

M-Endo Coliform Broth

M-Endo is a red colored media, which needs to be stored in the dark to prevent discoloration. Gram-positive bacteria are inhibited on this media by the desoxycholate and lauryl sulfate. The addition of ethanol increases the antibacterial nature of the formulation. Lactose fermenting organisms form aldehydes, which react with Schiff's reagent (basic fuchsin and sodium sulfite) to give red colored zones around the colonies. Coliform colonies are therefore red with a characteristic metallic sheen.

M-FC Broth

Allows the development of fecal coliforms at elevated temperatures (44.5°C).

M-FC with Rosolic Acid

M-FC with Rosolic Acid acts and functions in the same way as M-FC Broth. Rosolic acid inhibits bacterial growth in general, except for fecal coliforms.

M-Green Yeast and Mold Broth

M-Green Yeast and Mold Broth is an improved modification of the liquid media. The addition of bromocresol green, which diffuses into fungal colonies as an alkaline reaction, allows them to be easily identified. Metabolic by-products from the developing colonies diffuse into the surrounding medium, further reducing the pH which aids in the inhibition of bacterial growth, but also produces an acid reaction that causes residual bromocresol green to change to yellow.

M-Green Select Broth

M-Green Select Broth was developed to improve efficiency of detection and enumeration of fungi in sugar based drinks using the membrane filtration method. This medium has a low pH, which inhibits bacterial growth. The addition of chloramphenicol further inhibits the growth of bacteria to allow for the development and enumeration of yeast and mold.

MI Broth and MI Agar


MI Broth detects the presence of coliform bacteria by the production of β -galactosidase, which cleaves the substrate MU-Gal to produce 4-methylumbelliferone, which fluoresces on exposure to UV light. Non-coliforms do not produce this enzyme and therefore do not fluoresce on the medium. *Escherichia coli* is detected by the compound IBDG. The β -glucuronidase produced by *Escherichia coli* cleaves the substrate to produce a blue indigo color in the colonies. As *Escherichia coli* is also a total coliform, and also produces β -galactosidase, it will also fluoresce. The antibiotic cefsulodin is present to inhibit the growth of gram-positive bacteria and some non-coliform gram-negative bacteria that can cause false positive reactions.

MRS Broth


MRS medium supports luxuriant growth of all lactobacilli, even the slow growing species.

M-TGE Total Count Media

All bacteria develop on TGE media and produce a range of different colored and sized colonies.

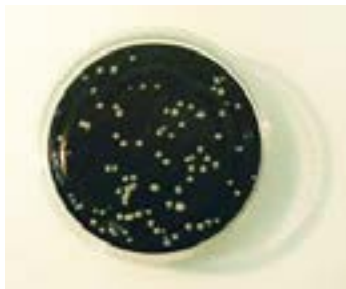

M-Green Yeast and Mold Broth: Typical Growth of *Candida Albicans* ATCC10231 on a Black Membrane

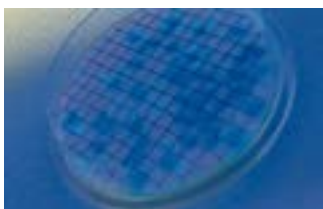

MI-Media: Pure Culture of *Escherichia coli* ATCC 25922 with UV Light

2.3 Nutrient Liquid Media

Orange Serum Media

Organisms known to grow in single strength and concentrated juices are lactic acid and acetic acid bacteria and yeast. Lactobacilli, Leuconostoc and yeast have all been identified as spoilage organisms by numerous authors. Orange serum at pH 5.4 to 5.6 has been reported to yield maximum counts of all types of spoilage organisms in mixed cultures and in single culture comparison tests.

Potato Dextrose Broth and Agar Media

Potato Dextrose Broth is recommended in Standard Methods as the media that gives the most consistent and highest counts for the recoveries of yeast and mold in dairy products. The inclusion of potato extract encourages the growth and development of fungi. Sterile tartaric acid may be added to lower the pH to 3.5 ± 0.2 to further inhibit the growth of conflicting bacteria.

PRY Broth

Preservative Resistant yeast Broth is a low pH selective medium for the detection of spoilage microorganism in beverages and water.

Pseudomonas Broth

Pseudomonas aeruginosa is characterized by the production of pyocyanin (a blue green, water soluble, non-fluorescent, phenazine pigment) which is stimulated by the inclusion of magnesium chloride and potassium sulfate in the broth. Irgasan, an antimicrobial agent, selectively inhibits gram-positive and gram-negative bacteria other than pseudomonads. Glycerol both serves as an energy source and helps in the promotion of pyocyanin.

Sabouraud Dextrose Broth

Peptone in the media is used as a nitrogen source for the development of fungi. Dextrose acts as an energy source for the growth of microorganisms. The low pH is favorable for the development of fungi, especially dermatophytes, but at the same time inhibits the development of contaminating bacteria from clinical specimens.

Standard Methods Agar

All bacteria develop on Standard Methods and produce a range of different colored and sized colonies.

Total Count Media with TTC

All bacteria develop on Total Count Media with indicator and produce a red color as a result of the precipitation of formazan following the reduction of 2,3,5-triphenyltetrazolium chloride (TTC) by bacteria.

Trypticase Soy Broth – Single Strength


General purpose medium used in qualitative procedures for the cultivation of fastidious and non-fastidious microorganisms. Trypticase Soy Broth – Single Strength complies with the demands of the DIN Norm 10167 for the detection of *Escherichia coli* serotype O157:H7 in foods and FDA-BAM for the isolation of enterohemorrhagic *Escherichia coli* (EHEC). In addition the media conforms to the formula of the US Pharmacopoeia.

Trypticase Soy Broth – Double Strength


TSB is a medium that will support the growth of a wide variety of microorganisms including aerobic, facultative, and anaerobic bacteria and fungi.

Wallerstein Nutrient Broth (WL) and WL Differential Broth (WLD)

WL Nutrient Broth is for the cultivation and enumeration of yeast and WL Differential Broth is for determination of bacterial count. Use of the medium at pH 5.5 and incubation at 25°C will give reliable counts for brewer's yeast. Adjustment of the pH to 6.5 and incubation at 30°C allows for the selective growth of baker's and distiller's yeast.

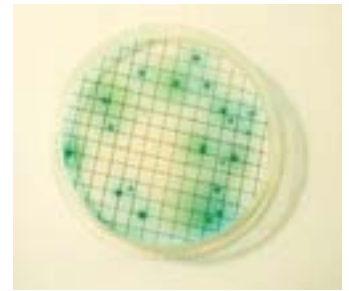

Pseudomonas Media: Typical Growth of *Pseudomonas aeruginosa* ATCC 10145

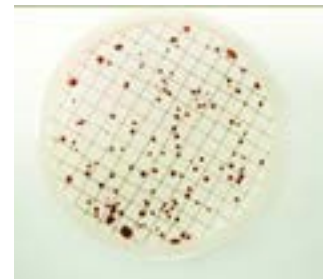

Total Count Media with Indicator. *Escherichia coli* ATCC 25922 and *Staphylococcus aureus* ATCC 25923 can be Easily Detected according to their Red to Pink Colonies



Trypticase Soy Broth Double Strength (not Inoculated)

MICROBIOLOGY

Product Code	Description
2 mL Ampoules	
10496146	Cetrimide Broth, 50/pk
10496120	Enterococcus Broth, 50/pk
10496164	Heterotrophic Plate Count (HPC) Broth, 2 ml , 50/pk
10496151	HPC Broth, 50/pk
10496125	KF-Streptococcus Broth, 50/pk
10496187	Lauryl sulfate PHLS/UK Broth 2ml 50/pk
10496121	Mannitol Salt Broth, 50/pk
10496187	Membrane Lauryl Sulfate PHLS/UK, 50pk
10496103	M-Endo Coliform Broth, 50/pk
10496124	M-FC media, 50/pk
10496114	M-FC Broth with rosolic acid, 50/pk
10496116	M-Green Select Broth, 50/pk
10496101	M-Green Yeast and Mold Broth, 50/pk
10496192	MI-Broth Media, 50/pk
10496112	MRS Broth, 50/pk
10496102	M-TGE Broth, 50/pk
10496104	Orange Serum Broth, 50/pk
10496106	Pry Both 2ml 50/pk
10496119	Pseudomonas Broth, 50/pk
10496113	Total Count Broth with TTC, 50/pk
10496108	Wallerstein Broth, 50/pk
10496109	Wallerstein Differential Broth, 50/pk
9 mL Tubes	
10496710	Brilliant Green Bile Bottled Broth, with Durham tubes, 20/pk
10496714	EC Bottled Broth, with Durham tubes, 20/pk
10496709	EC with MUG, Bottled Broth, 20/pk
10496722	Lauryl Sulfate/ Lauryl Tryplose Broth
Bottled Media	
10496700	M-Endo Coliform Bottled Broth, 50ml 8/pk
10496851	MI Media, Bottled Broth, 50 ml, 1/pk
10496847	MI Media, Bottled Agar, 50 ml, 1/pk
10496705	M-Green Yeast and Mold Bottled Agar, 100 ml, 1/pk
10496713	Orange Serum Bottled Agar, 100ml 1/pk
10496731	Potats Dextros e Bottled Agar, 100ml 1/pk
10496706	Standard Bottled Agar, 100ml 1/pk
10496707	Trypticase Soy Broth (TSB) Single strength, Bottled Broth, 100 ml, 1/pk
10496708	Trypticase Soy Broth (TSB) Double strength, Bottled Broth, 100 ml, 1/pk
Rapid Test Kit	
10496744	ColiCheck with MUG, PIA Test Kit with Sample Bottles, 30/pk

2.4 SwabCheck



The SwabCheck principle

The surface is wiped with a cellulose swab and any bacteria collected are transferred via the swab into a tube containing a special medium with an indicator dye, which is then incubated. A single bacterium is sufficient to cause a color change. This means that SwabCheck is about 1000 times more sensitive than the conventional ATP method. This accuracy is particularly important in the food industry. With this simple method, it is possible to identify microorganisms such as *Listeria monocytogenes*, which must not be present in any concentration in food and beverages.

Features and Benefits:

- The right test for each type of contamination
- Qualitative and semi-quantitative hygiene control
- Sterile packed and ready-for-use
- Easy to handle
- Rapid results
- Long shelf-life

SwabCheck use

Handling is easy.

Open the sterile pack, remove the swab and wipe it over an area of about 10 x 10 cm. Then twist off the cap of the medium tube and insert the swab so that the cap fits tightly. Label the sample tube and incubate at the appropriate temperature.

A change in color indicates the presence of the microorganism in question. The quicker the color change occurs, the higher the bioburden. If no color change has been observed after the maximum incubation period has elapsed, then the corresponding microorganism is not present. GVS Life Sciences offers SwabCheck in packs of 25 pieces. Shelf-life of 12 months.

Neutralizing Buffer Swabs

Neutralizing buffer swabs are used in the monitoring of surfaces for total bacterial count. Neutralizing buffer inactivates the bactericidal and bacteriostatic effects of chlorine and quaternary ammonium detergents. However, neutralizing buffer exhibits no toxic effects on microorganisms. This permits the transfer of swabbed organisms to the laboratory without loss in viability. Neutralizing buffer is not designed to culture and enumerate microorganisms.

Buffer Swabs

Used for the collection of surface contamination from flat or convoluted surfaces prior to transport to a laboratory for culture and enumeration. Buffer swabs contain no bacteriostatic or bactericidal compounds and cannot suppress the action of detergents.

SwabCheck


Used as an indication of hygiene on contact surfaces. SwabCheck changes color from purple to yellow. The color change is based on acid reaction with the indicator. The more rapid the color change, the higher the level of bacteria in the sample. SwabCheck is useful in determining the sanitation levels of preparation surfaces, filling ports, and processing areas in beverage and food processing plants, dairies, restaurants, and healthcare facilities.

Coliform SwabCheck

Escherichia coli and coliforms are used traditionally as indicator organisms for fecal contamination in water and other environmental samples. Detection of these organisms usually points to poor hygiene at some stage in the production process or pollution of water at source. The presence of coliforms is indicated by a color change from brown to yellow. The more rapid the color change the higher the level of coliform bacteria.

Hygiene SwabCheck

Easy to use: The Hygiene SwabCheck shows an obvious color change from red to yellow. The time taken for this change is an indication of the level of contamination. This should be used in conjunction with known specification levels of your process/product. Rapid screening hygiene test is a same day test that will detect gross bacterial and fungal contamination of work surfaces, equipment machinery or other sampling sites.

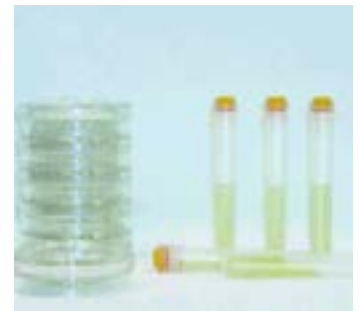

Total Count Swab Kit



Coliform SwabCheck

MICROBIOLOGY

Listeria SwabCheck

Listeria Isolation SwabCheck is designed to be used alongside traditional selective methods to improve the quality system and minimize the risk of Listeria contamination. This simple to use diagnostic test can be applied anywhere in the environment and on foodstuffs where the presence of Listeria species would be critical.

Listeria sp and specifically Listeria monocytogenes are rapidly becoming the most important pathogen in the food industry; regulatory bodies from around the world are insisting that all food products are Listeria free. Listeria Isolation SwabCheck works on an enhanced Esculin media formulation. The hydrolysis of esculin gives a distinctive black/brown precipitate. Inhibitors and antibiotics are present in the media, which will inhibit the growth of non-Listeria species.

SwabCheck Escherichia coli

Used for the detection of Escherichia coli on surfaces. The presence of fluorescence using a longwave UV light source confirms the presence of Escherichia coli and any further confirmation is not required. MUG detects anaerogenic strain that may not be detected in the conventional procedure. Lactose is a source of energy. Casein peptone provides additional nutrients. The mixture of bile salts is inhibiting for gram-positive bacteria, particularly bacilli and fecal streptococci. The substrate 4-methylumbelliferyl-b-D-glucuronide is hydrolyzed by an enzyme, b-glucuronidase, possessed by most Escherichia coli and a few strains of Salmonella, Shigella, and Yersinia, to produce a fluorescent end product, 4-methylumbelliferone. The presence of Escherichia coli is detected by the appearance of fluorescence throughout the tube.

Total Count Swab Kit

Used for the non-selective development and enumeration of all aerobic bacteria on surfaces in accordance with Hazard Analysis and Critical Control Points (HACCP). The kit includes the swabs and culture medium, packaged with a membrane device, providing a quantitative result. All bacteria develop on TGE media and produce a range of different colored and sized colonies. It is not possible using TGE to presumptively identify any bacteria. Identification can only be undertaken using traditional microbiology techniques following initial colony development.

Yeast and Mold Swab Kit

Used for the enumeration of yeast and molds on surfaces in accordance with HACCP. The kit includes the swabs and culture medium, packaged with a membrane device, providing a quantitative result. M-Green yeast and mold is an improved modification of the liquid medium, and was developed to improve efficiency of detection and enumeration of fungi in sugar based drinks using the membrane filtration method. This medium has a low pH, which inhibits bacterial growth. The addition of bromocresol green, which diffuses into fungal colonies as an alkaline reaction, allows them to be easily identified. Metabolic by-products from the developing colonies diffuse into the surrounding medium, further reducing the pH that aids in the inhibition of bacterial growth, but also produces an acid reaction that causes residual bromocresol green to change to yellow. Green opaque colonies against a yellow background are indicative of the growth of yeasts. Mold colonies are green and filamentous.

Polywipe Sponge

Used for the recovery of microorganisms from a surface. Polywipe is a blue sponge that is premoistened with neutralizing buffer to neutralize the effects of surface disinfectants. The sponge material is selected to be free of the preservatives found in commercially available sponges, which can inhibit microorganism growth. Polywipe sponges are biocide free and tested for zero toxicity to microorganisms. Each sponge is individually wrapped in a peel pouch and gamma irradiated to ensure sterility.

Buffer Swabs

Used for the collection of surface contamination from flat or convoluted surfaces prior to transport to a laboratory for culture and enumeration. Buffer swabs contain no bacteriostatic or bactericidal compounds and cannot suppress the action of detergents.


Hygiene SwabCheck



Listeria SwabCheck



Yeast and Mold Swab Kit

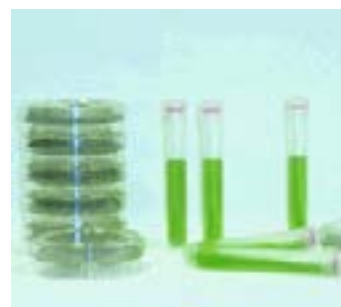

Polywipe Sponge



2.4 SwabCheck

Ordering information

Product Code	Description
10498303	Neutralizing Buffer Swabs, 4 ml, 125/pk
10498304	Neutralizing Buffer Swabs, 4 ml, 500/pk
10498305	Buffer Swabs, 4 ml, 125/pk
10498306	Buffer Swabs, 4 ml, 500/pk
10498404	SwabCheck, 4 ml, 125/pk
10498402	SwabCheck Escherichia coli, 4 ml, 125/pk
10498315	Total Count Swab Kit, 30/pk
10498316	Yeast and Mold Swab Kit, 30/pk
10498406	Coliform SwabCheck, units ready to use, 25/pk
10498407	Hygiene SwabCheck, units ready to use, 25/pk
10498408	Listeria SwabCheck, units ready to use, 25/pk
10498521	Polywipe Sponge, ready to use, single packed,50/pk

MICROBIOLOGY

2.5 Dilution Bottles



Description and Use

Prefilled sterile dilution bottles are designed for sample dilution of water, dairy products, foods, and pharmaceuticals prior to microbiological testing. Final pH for all solutions is 7.2 pH \pm 0.2 pH at 25°C. They come in an easy open, flip-top, plastic container with a tamper-evident seal.

Butterfield's Phosphate Buffer contains monobasic potassium phosphate and is used extensively in the food, dairy, and pharmaceutical industries. Offered in 90 ml and 99 ml volumes for easy 1:10 and 1:100 dilutions. It is recommended as a general diluent in laboratory procedures by the Federal Drug Administrations and in the Bacteriological Analytical Manual. This product is prepared according to Standard Methods for the Examination fo Water and Wastewater for use in water testing.

Phosphate Buffer with magnesium chloride is used asthe diluents for the preparation of dilutions in plate counts in the dairy and food industries. It is recommended y the APHAfor the recovery of injured microorganisms form dairy and food samples. Contains deionized water, monopotassium phosphate, and magnesium chloride.

Ordering information

Product Code	Description
10498503	Dilution Bottle, Butterfield's Buffer, 99 ml, 72/pk
10498504	Dilution Bottle, Butterfield's Buffer, 90 ml, 72/pk
10498505	Dilution Bottle, Phosphate Buffer Magnesium Chloride, 99 ml, 72/pk



FILTRATION MEMBRANES

3. FILTRATION MEMBRANES

3.1 Membranes for Filtration

3.1.1 - Cellulose Acetate (CA) Membrane



Description and Use

GVS Life Sciences Cellulose Acetate (CA) Filtration Membrane is a supported, hydrophilic membrane that is naturally low binding. It is ideal for use in filtration applications where maximal recovery of protein is critical.

Features and Benefits

- **Superior strength:** Can withstand aggressive handling or be used with automated equipment without breaking or tearing.
- **Low extractables:** Ensures tests will be clean with consistent results.
- **Hydrophilic:** Wets out rapidly.
- **Lot-to-lot consistency:** Quality checks ensure consistent flow and diffusion rates for dependable results every time.
- **Nonlysing of cells:** Prevents contamination of critical solutions.

Typical Applications

- Protein and enzyme filtration
- Biological fluid sterilization
- Tissue culture media sterilization
- Cold sterilization

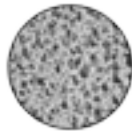


Figure 2: Scanning Electron Micrograph of GVS Life Sciences CA Filtration Membrane

Exceptional Strength for Improved Performance

Produced through a proprietary manufacturing process, GVS Life Sciences CA Membrane meets rigorous quality standards throughout every step of production. This process generates consistent lot-to-lot filtration properties among the membranes and ensures product uniformity.

GVS Life Sciences CA Filtration membranes are composed of pure cellulose acetate that is internally supported by an inert polyester web. This web gives each membrane exceptional strength to prevent cracking, tearing, breaking and distortion when handled or creased. The resulting membrane has dimensional stability that can withstand autoclaving or steam sterilizing leaving the membrane unaffected in temperatures up to 274°F (135°C).

The exceptional dimensional strength and low binding characteristics of GVS Life Sciences CA Filtration Membranes provides higher throughputs than competitive offerings and reduces the amount of filter changes needed during proteinaceous solution filtering. Its uniform pore size and consistent flow rates ensure reliable performance.

Flexibility for OEM Requirements

GVS Life Sciences CA Filtration Membrane is available in rolls from 25mm to 30cm wide, as well as sheets and cut disks that can be customized to meet your application and size requirements. GVS Life Sciences CA Filtration Membranes are manufactured on-site making customization easy and cost effective.

Table 2: Product Characteristics

USP Class VI testing	Passed
Thickness	65 - 110 μm
Maximum Operating Temperature	274°F (135°C)
Sealing Compatibility	Ultrasonics, Heat, Radio Frequency and Insert Molding
Pore Size Range	0.1 to 20 μm

Table 3: Performance Characteristics

Pore Size	0.22 μm	0.45 μm	0.6 μm	0.8 μm	1.2 μm	5.0 μm	10.0 μm	20 μm
Minimum Bubble Point psi (kg/cm ²)	50 (3.50)	30 (2.10)	18 (1.26)	14 (0.98)	11 (0.77)	6 (0.42)	5 (0.35)	3 (0.21)
Typical Flow Rate, mL/min/cm ² @ 10psi (0.7 kg/cm ²)	16.1 (1.13)	54.7 (3.85)	70.9 (4.98)	81.3 (5.72)	180 (12.7)	375 (26.4)	592 (41.6)	1478 (104)

Ordering information: Cellulose Acetate (CA) Membrane

Dimensions Packaging	13mm 100/pk	25mm 100/pk	47mm 100/pk	50mm 100/pk	90mm 25/pk	102mm 25/pk	142mm 25/pk	293mm 25/pk	20x20mm 5/pk	30cmx3m 1/pk
0.22 μm	1212374	1213124	1213804	1221730	1214357		1215074	1215427		1224211
0.45 μm	1215533	1215635	1215676	3052874	1212375	1221546	1212517	1212620		1240382
0.65 μm		1212846	1212942		1213037		1213125		3061196	
0.8 μm	1213305	1213343	1213358				1213516		3034974	3034975
1.2 μm		1213730	1213805				1213958	1214038		3041202
5.0 μm		1214370	1214411		1212648		1214851			3049247

3.1 Membranes for Filtration

3.1.2 - Nitrocellulose (NC) Membrane



Description and Use

GVS Life Sciences Nitrocellulose (NC) Filtration Membrane is an unsupported, hydrophilic membrane. Its rapid flow rate and high throughput make it ideal for use in diagnostic kit manufacturing applications.

Features and Benefits

- **High flow rate:** Provides fast filtration rates
- **Uniform pore structure:** Provides consistent flow and diffusion rates
- **< 4% extractables:** Leads to more consistent results
- **Lot-to-lot consistency:** Produces dependable results every time

Typical Applications

- Aqueous filtration
- Sterility testing
- Gravimetric analysis with ashing technique
- Microbiological and particulate analysis
- Black for food and beverage applications

Consistent Uniformity Improves Control and Performance

Produced through a proprietary manufacturing process, GVS Life Sciences Nitrocellulose meets rigorous quality standards throughout every step of production. This process ensures the membrane has a uniform pore structure and consistent thickness (within 10 μm).

GVS Life Sciences NC Filtration Membranes are composed of a mixture of inert cellulose nitrate and cellulose acetate polymers. The uniform microporous structure of these filters provides the fastest flow rates and highest throughputs available in a membrane filter. Because they are biologically inert, GVS Life Sciences NC Filtration Membranes are ideal for a wide range of clarification, sterilization and analytical applications such as: microbiological analysis, clarification or sterilization of aqueous solutions, industrial hygiene applications, silt density index and particulate-matter analysis. For gravimetric analysis using ashing techniques, GVS Life Sciences NC Membranes yield a residue or less than 0.045% of their initial weight. They are hydrophilic with a noncytotoxic wetting agent and yield extractable levels of less than 4% of their weight. These membranes are autoclavable at 121°C (250°F) for 20 minutes. Sterilized product lifetime is 18 months from sterilization date (1 year warranty).

Flexibility for OEM Requirements

GVS Life Sciences Nitrocellulose is available in rolls from 25 mm (1.0 in) to 40 cm (31 inches) wide. It is also available in sheets and cut disks gridded, plain, white or black, that can be customized to meet your application and size requirements. Slitting tolerance is from +/- 0.039 inches (+/- 1 mm). Because GVS Life Sciences Nitrocellulose Filtration Membrane is manufactured on-site all customization, such as wicking rate and thickness, is easy and cost-effective.

Table 2: Product Characteristics

Sterilization	Gamma Irradiation or Ethylene Oxide (EtO)
USP Class VI testing	Passed
Thickness	Approx. 6 mil (150 μm) +/- 10 μm
Extractables	< 4%
BSA Protein Binding	Approx. 160 $\mu\text{g}/\text{cm}^2$ (depending on pore size)
Maximum Operating Temperature	356°F (180°C)
Sealing Compatibility	Ultrasonic, Heat, Radio Frequency and Insert Molding
Pore Size Range	0.1 to 8.0 μm

Table 3: Performance Characteristics

Pore Size	0.1 μm	0.22 μm	0.45 μm	0.8 μm	1.2 μm	5.0 μm	8.0 μm
Minimum Bubble Point psi (kg/cm ²)	80 (5.62)	52 (3.66)	30 (2.11)	11 (0.77)	9 (0.63)	6 (0.42)	4 (0.28)
Typical Flow Rate, mL/min/cm ² @ 10psi (0.7 kg/cm ²)	6.9 (0.49)	19 (1.34)	51 (3.59)	198 (13.9)	265 (18.6)	650 (45.7)	2316 (163)

Ordering information: Nitrocellulose Gridded Filtration Membrane, White and Black

Pore sizes	Bulk Packaging*				Individual Sterile Package Without Pad*						
	25mm Packaging	25mm	47mm	47mm	47mm	47mm	47mm	47mm	47mm	47mm	50mm
	100/pk	100/pk	100/pk	100/pk	100/pk	100/pk	200/pk	200/pk	1000/pk	1000/pk	1000/pk
Color	white	black	white	black	white	black	white	black	white	black	white
0.22 μm	1214872				1216720		1214861		1214396		
0.45 μm	1215237	1214866	1215249	1213145	1216721	1216719	1215230	1214991	1214923	1213643	1222980
0.7 μm	1215407		1215409		1216722	1216718	1215406	1213331	1215408	1221948	
0.8 μm	1225460				1216724	1216723					1215590

FILTRATION MEMBRANES

		Individual Sterile Package With Pad*			
Dimensions Packaging	47mm 100/pk	47mm 100/pk	47mm 1000/pk	47mm 1000/pk	
Color	white	black	white	black	
0.22 μm	1214872				
0.45 μm	1215237	1214866	1215249	1213145	
0.7 μm	1215407		1215409		
0.8 μm	1225460				

*Larger bulk packaging available, call for details.

Ordering information: Nitrocellulose Filtration Membrane, White and Black

Dimensions Packaging	13mm 100/pk	25mm 100/pk	25mm 100/pk	37mm 100/pk	47mm 100/pk	47mm* 100/pk
Color	white	white	black	white	white	black
0.1 μm		1214527			1214533	
0.22 μm	1214882	1214898			1214909	
0.45 μm	1215257	1215263	1215019	1215272	1215281	
0.65 μm		1215376			1215380	
0.8 μm	1215424	1215425	1215415	1215426	1215428	1215416
1.2 μm	1215438	1215440			1215441	
5.0 μm	1215448	1215450			1215451	
8.0 μm	1214456	1215455			1215456	

*Sterile and bulk packaging available, call for details.

Ordering information: Nitrocellulose Filtration Membrane, White and Black (cont.)

Dimensions Packaging	90mm 25/pk	142mm 25/pk	293mm 25/pk	20x20cm 5/pk	20x20cm 5/pk
Color	white	white	white	white	black
0.1 μm		1214554	1214565		
0.22 μm	1214941	1214950	1214959	3031100	
0.45 μm	1215305	1215316	1215323	1225781	3053082
0.8 μm	1215431	1215432	1215433	3050851	
1.2 μm	1215442	1215443			
5.0 μm	1215452	1215453			
8.0 μm	1215027	1221955	1212631		

3.1 Membranes for Filtration

3.1.3 - Nylon (NY) Membrane



Description and Use

GVS Life Sciences Nylon Filtration Membrane is a supported, naturally hydrophilic membrane designed to wet out evenly and retain its superior strength during use in general filtration or medical assays

Features and Benefits

- **Hydrophilic:** Eliminates the need for wetting agents that can potentially interfere with biological processes
- **Super strength:** Eases handling when used with automated equipment
- **Low extractables:** Ensures tests will be clean and pure leading to more consistent results
- **Lot-to-lot consistency:** Quality checks ensure lot-to-lot consistency, both down and across the polyester web, for dependable results every time

Typical Applications

- Sterilization and clarification of aqueous and organic solvent solutions
- HPLC sample preparation

Versatile Capabilities, Consistent Performance

Produced through a proprietary manufacturing process, GVS Life Sciences Nylon Filtration Membrane meets rigorous quality standards throughout every step of production. This process generates consistent lot-to-lot flow rates among the membranes ensuring product uniformity.

GVS Life Sciences Nylon Filtration Membrane is internally supported with an inert polyester support web giving it added dimensional strength and stability that prevents cracking, tearing, curling and breaking. This added strength and durability is advantageous during usage that involves aggressive handling or automated equipment.

A naturally hydrophilic membrane, GVS Life Sciences Nylon Filtration Membrane does not require wetting agents that can interfere with biological processes. The resulting membrane has a void volume of 70% to 80% for high diffusion and low-flow resistance.

Table 3: Performance Characteristics

Pore Size	0.1 μ m	0.22 μ m	0.45 μ m	0.6 μ m	0.8 μ m	1.2 μ m	5.0 μ m	10.0 μ m	20.0 μ m
Minimum Bubble Point psi (kg/cm ²)	70 (4.92)	50 (3.51)	35 (2.11)	18 (1.27)	13 (0.91)	11 (0.77)	6 (0.42)	5 (0.35)	3 (0.21)
Typical Flow Rate, mL/min/cm ² @ 10psi (0.7 kg/cm ²)	4.0 (0.28)	9.9 (0.70)	26.9 (1.89)	59.3 (4.17)	80.5 (5.66)	180 (12.7)	331 (23.3)	552 (38.8)	1448 (101.9)

Flexibility for OEM Requirements

GVS Life Sciences Nylon Filtration Membrane is available in rolls from 1cm to 33cm wide, as well as sheets, cut disks and pleated packs that can be customized to meet your application and size requirements. Because the GVS Life Sciences Nylon Filtration Membrane is manufactured on-site, all customization is easy and cost-effective.

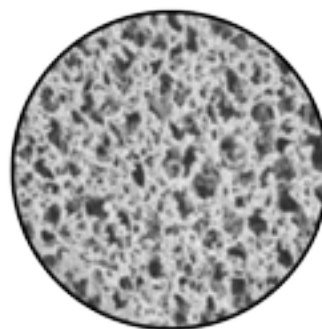


Figure 2: Scanning Electron Micrograph of GVS Life Sciences Nylon Filtration Membrane

Table 2: Product Characteristics

Sterilization	Gamma Irradiation or Ethylene Oxide (EtO)
USP Class VI toxicity	Passed
Thickness	65 - 125 μ m
Extractables	< 0.2% (< 0.0015 mg/cm ²)
BSA Protein Binding	Approx. 120 μ g/cm ²
Maximum Operating Temperature	356°F (180°C)
Sealing Compatibility	Ultrasonics, Heat, Radio Frequency and Insert Molding
Pore Size Range	0.1 to 20 μ m

FILTRATION MEMBRANES

Ordering information: Nylon Membrane

Dimensions Packaging	13mm 100/pk	25mm 100/pk	37mm 100/pk	47mm 100/pk	47mmGRD 100/pk
0.1 μm	1213760	1213761		1213762	
0.22 μm	1213766	1213768		1213769	
0.45 μm	1213774	1213775	1228824	1213776 1220671**	1213825 1213845 ~
0.65 μm		1213782		1213783	
0.8 μm	1213788	1213789	1214881	1213790	3013826
1.2 μm	1213794	1213796	1230356	1213797	1214880
5.0 μm	1213810	1213811	1236904	1213812	3048260
10.0 μm	1213817	1213818		1213819	
20.0 μm	1213801	1213802		1213803	

Dimensions Packaging	90mm 25/pk	142mm 25/pk	293mm 25/pk	200x200mm 5/pk	30cmx3m 1/pk
0.1 μm	1213763	1213764	1213765	1222859	1241477
0.22 μm	1213770	1213771	1213772	1222858	1224690
0.45 μm	1213778	1213779	1213780	1222857	1225982
0.65 μm	1213784	1213786		1222856	3052148
0.8 μm	1213791	1213792	1213793	1222855	
1.2 μm	1213798	1213799	1213800	1222854	1214956
5.0 μm	1213813	1213815	1213816	1222851	1221441
10.0 μm	1213820	1213821	1213823	1222852	
20.0 μm	1213807	1213808	1213809	1222853	

GRD = Gridded

**sterile

3.1 Membranes for Filtration

3.1.4 - Polyethersulfone (PES) Membrane



Description and Use

GVS Life Sciences Polyethersulfone (PES) Filtration Membrane is hydrophilic and cast from pure polyethersulfone polymer. It is designed to remove particulates during general filtration and its low protein and drug binding characteristics make it ideally suited for use in life science applications.

Features and Benefits

- **Hydrophilic:** Eliminates the need for wetting agents that can potentially interfere with analyses
- **Low extractables:** Ensures test results will not be compromised by wetting agents or other extractables
- **Low drug and protein binding:** Maximizes recovery of critical drugs or proteins
- **Wide range of pore sizes:** Pore size range of 0.03 μm to 5.0 μm enables specific pore size selection for given applications
- **Superior burst strength:** Protects the integrity of the membrane under high pressure
- **Lot-to-lot consistency:** Quality checks, both down and across the membrane, ensure dependable results every time

Typical Applications

- Protein and enzyme filtration and sterilization
- Biological fluid filtration and sterilization
- Pharmaceutical sterilization
- Environmental water studies

Table 3: Performance Characteristics

Pore Size	0.03 μm	0.1 μm	0.22 μm	0.45 μm	0.65 μm	0.8 μm	1.2 μm	5.0 μm
Flow Time (secs)	200-500	100-200	35-70	20-40	12-25	80-160	65-130	50-100
Volume/Vacuum (mls/in Hg)	250/20	250/20	250/20	250/20	250/20	500/5	500/5	500/5
Flow Rate (ml/min/cm ² @10psi)	7.95-3.18	15.91-7.95	45.45-22.72	79.53-39.77	132.55-63.63	159-80	196-98	255-127
Bubble Point (psi)	90-110	70-90	50-70	35-50	21-32	13-28	11-22	6-13
Thickness (μm)	110/150	110/150	110/150	110/150	110/150	110/150	110/150	110/150

Ordering information: Polyethersulfone Filtration Membrane

Dimensions Packaging	25mm 100/pk	47mm 100/pk	90mm 25/pk	200x200mm 5/pk	30cmx3m 1/pk
0.03 μm	3032876	3029505	3018505	1235748	3057106
0.1 μm		1214756	1222230	1225881	3026365
0.22 μm	1214193	1214465	1214920	1223871	1226664
0.45 μm	1214532	1214475	1215368	1225882	1226665
0.65 μm		1224487		1225883	1225985
0.8 μm	1214604	1214568	1214669	1225884	3037376
1.2 μm	1222267	1221008	1224492	1223340	1242278
3.0 μm				1232921	
5.0 μm	1224003	1215396	1224496	1236292 1233863*	3030900**

**hydrophobic *30pk

Product Uniformity and High Sensitivity Maximize Performance

GVS Life Sciences PES Filtration Membrane is produced through a proprietary manufacturing process that employs high-quality standards throughout every step of production. This level of quality generates lot-to-lot consistency and product uniformity that results in reliable flow rates, particle retention and fluid purification.

This strong, microporous film membrane is constructed from a high-temperature polyethersulfone polymer that is acid and base resistant. Its strength and durability are advantageous during usage that involves aggressive handling or automated equipment. GVS Life Sciences PES Filtration Membrane is naturally hydrophilic without added wetting agents and has low extractables.

Due to its inherent uniform porosity and controlled pore size, GVS Life Sciences PES Filtration Membrane efficiently removes particulates from solutions during general filtration. Additionally, its low protein and drug binding characteristics maximize recovery of critical drugs used in I.V. therapy, chemotherapy and open-heart surgery.

Flexibility for OEM Requirements


GVS Life Sciences PES is available in rolls 1.3cm - 79cm wide, as well as sheets, cut disks, capsules, pleat packs and cartridges that can be customized to meet your application requirements. Because the GVS Life Sciences PES is manufactured on-site, all customization can be done easily and cost-effectively.

Table 2: Product Characteristics

Endotoxin levels and Class VI toxicity	Passed
Thickness	110 - 150 μm
Extractables	< 2% (< 0.015 mg/cm ²)
Maximum Operating Temperature	266°F (130°C)
Sealing Compatibility	Ultrasonics, Heat, Radio Frequency and Insert Molding
Pore Size Range	0.03 to 5.0 μm

FILTRATION MEMBRANES

3.1.5 - Polypropylene (PP) Membrane

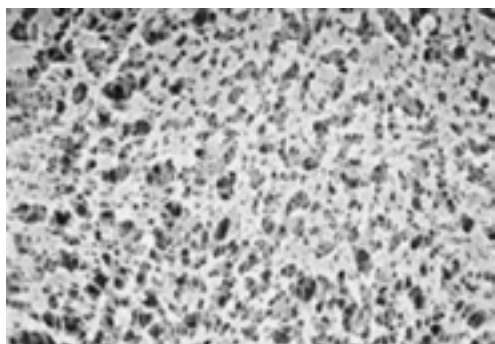

Description and Use

GVS Life Sciences polypropylene membrane filters are composed of pure polypropylene with absolute pore size ratings. These filters offer broad chemical compatibility allowing its use with aqueous and organic solvent samples. The polypropylene filter has extremely low extractable levels designed to provide accurate, consistent analysis results for sensitive ion chromatography applications while prolonging

column life. GVS Life Sciences polypropylene filter is the preferred filter membrane for HPLC applications where the detection levels are below 230 nm. The filters also exhibit negligible protein binding which, is essential for maximum sample recovery of critical, small volume protein samples. GVS Life Sciences filters are available in 0.1, 0.22 and 0.45-micron pore sizes in a wide range of disc diameters.

Features and Benefits

- Broad chemical compatibility
- Hydrophobic
- HPLC applications - detection levels < 230 nm

Typical Applications

- Aqueous and organic solvent filtration
- HPLC sample preparation requiring low detection levels
- Ion chromatography
- Total digest for heavy metals


Ordering information: Polypropylene Membrane

Dimensions Packaging	25mm 100/pk	47mm 100/pk	90mm 25/pk	142 mm 25/pk	200x200 mm 5/pk	254x3000 mm 1/pk
0.1 μm	1222102	1214237	1220824	1222103	1225932	1269206
0.22 μm	1214238	1214239	1214240	1214241	3095433	1224966
0.45 μm	1212379	1212380	1212381	1212383		
0.65 μm		1225490*				
1.0 μm		1225489*				
1.2 μm	1212390	1212391	1212392	1212394		
2.0 μm		3026845*				
5.0 μm		1225487*				
10.0 μm		1225792				

*40/pk

3.1 Membranes for Filtration

3.1.6 - Polytetrafluoroethylene (PTFE) Membrane

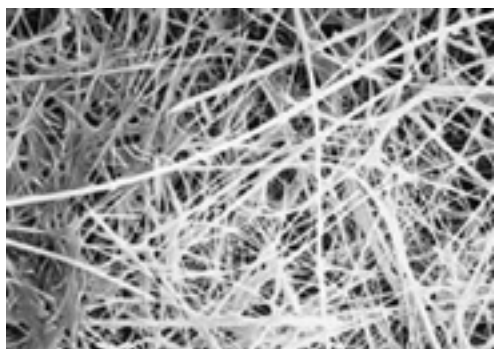

Description and Use

PTFE (fine powder resin) is expanded into a 3-dimensional web-like structure called PTFE which creates billions of microscopic pores. This structure utilizes the inherent hydrophobic (water-resistant) and non-stick nature of PTFE to allow removal of particulate captured on the membrane surface. This allows air to pass

easily through the membrane while collecting particulate as small as 0.1 micron on its surface. PTFE membranes and laminates provide device manufacturers with a consistent, temperature and chemical compatible barrier to microbes and particulate matter. The optimal combination of air flow and water entry pressure adds value to most device designs.

Features and Benefits

- Naturally hydrophobic
- Compatible with strong acids and aggressive solutions
- Improved durability and handling

Typical Applications

- Filtration of strong acids and aggressive solutions
- Venting applications
- Phase separations
- Aerosol samplings

Ordering information: Polytetrafluoroethylene Membrane

Pore sizes	Dimensions	13 mm	25 mm	47 mm
	Packaging	100/pk	100/pk	100/pk
0.22 μm		1215485	1215486	1215487
0.45 μm		1215491	1215492	1215493
1.0 μm		1215502	1215503	1215504 3013362*

Pore sizes	Dimensions	90 mm	142 mm	293 mm	200x200 mm	305x305 mm
	Packaging	25/pk	25/pk	25/pk	5/pk	50/pk
0.22 μm		1215488	1215489	1215490	3026028	1267681
0.45 μm		1215494	1215495	1215496	1237423	3034300
1.0 μm		1215505	1215506	1215507	1214443	1235299

*25/pk

FILTRATION MEMBRANES

3.1.7 - Polytetrafluoroethylene (PTFE) Hydrophilic Membrane



Membrane Material:
Polytetrafluoroethylene
Membrane Diameter (mm):
25 and 47
Sterile:
No
Sterilization:
ETO, autoclave 30 min at 121 C
Max operating temp (°C):
135

General Application:
Filtration of Aqueous and Organic Solutions
Analytical Sample Prep, HPLC
Chromatography
Clarification
Fuel Hydraulic Fluids and Machined Parts

Ordering information: Polytetrafluoroethylene Hydrophilic Membrane

Product Code	Description	Pore Size	Water flow rate (ml/min/cm ²) @10 psi	Water Bubble Point (psi)
ME025NPH002BC01	Membrane Ø25mm PTFE hydrophilic 0.2 µm GVS Life Sciences 100/pc	0.22 µm	16	45
ME047NPH002BC01	Membrane Ø47mm PTFE hydrophilic 0.2 µm GVS Life Sciences 100/pc	0.22 µm	54	27

3.1.8 - Regenerated Cellulose (RC) Membrane



Membrane Material:
Regenerated Cellulose
Membrane Diameter (mm):
25 and 47
Sterile:
No
Sterilization:
Autoclave 121C, or dry heat at 180C for 2 hours, or
Gamma (25 kGy), or ETO
Max operating temp (°C):
134

General Application:
Filtration of Aqueous and Organic Solutions
Analytical Sample Prep, uHPLC
Chromatography
Clarification
Protein Chemistry

Detailed Application: General filtration, particle separation, degassing solvents, filtration requiring a strong supporting membrane. Resistant to many solvents, and to aqueous solutions in the pH range from 3 to 12. Often used for ultra-cleaning and de-gassing solvents and mobile phases for HPLC. Excellent chemical compatibility and resistance to organic solvents. Compatible with almost all solvents. Superior thermal resistance. Features low nonspecific adsorption (bovine serum albumin < 10 micrograms/cm²).

Ordering information: Regenerated Cellulose Membrane

Product Code	Description	Pore Size	Water flow rate (ml/min/cm ²) @10 psi	Water Bubble Point (psi)
ME025NRC002BC01	Membrane, Ø25 mm, RC 0.20 µm	0.20 µm	16	57
ME025NRC004BC01	Membrane, Ø25 mm, RC 0.45 µm	0.45 µm	28	36
ME047NRC002BC01	Membrane, Ø47 mm, RC 0.20 µm	0.20 µm	16	57
ME047NRC004BC01	Membrane, Ø47 mm, RC 0.45 µm	0.45 µm	28	36

3.1.9 - Polyvinylidene Fluoride (PVDF) Hydrophilic Membrane



Membrane Material:
Polyvinylidene Fluoride
Membrane Diameter (mm):
25 and 47
Sterile:
No
Sterilization:
Autoclave (121 °C at 1 bar), Gamma, Beta, ETO
Max operating temp (°C):
85

General Application:
Filtration of Aqueous and Organic Solutions
Analytical Sample Prep, uHPLC
Chromatography
Clarification
Protein Chemistry



Detailed Application: Sterilizing filtration of biological solutions. Excellent chemical compatibility with aggressive solvents, acids and alcohols. HPLC.

Ordering information: Polyvinylidene Fluoride Hydrophilic Membrane

Pore Size	25 mm 100pk	47mm 100pk	90mm 25pk
0.22	3044272	3044270	3044271
0.45	3037802	3037800	3037801

3.1 Membranes for Filtration

3.1.10 - Polycarbonate Track Etched (PCTE) Membrane

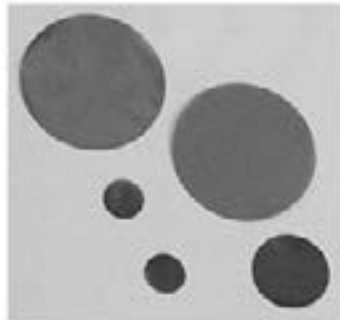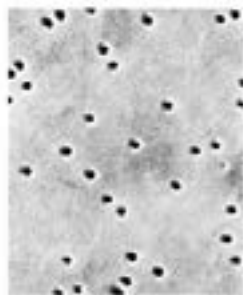

Description and Use

GVS Life Sciences PCTE Membrane is made from a thin, microporous polycarbonate film material. It is ideally suited for use in blood assays and high-purity and general filtration.

Standard membrane is hydrophilic due to PVP treatment. PVP-free hydrophobic membrane also available.

Typical Applications

- General filtration
- Removal of red blood cells from plasma
- Flow control of reagents through assay
- Precise filtration and prefiltration

Table 2: Product Characteristics

Sterilization	Gamma Irradiation or Ethylene Oxide (EtO)
USP Class VI Testing	Passed
Extractables	Very Low
BSA Protein Binding	5 $\mu\text{g}/\text{cm}^2$
Maximum Operating Temperature	284°F (140°C)
Sealing Compatibility	Ultrasonic, Heat, Radio Frequency and Insert Molding
Pore Size Range	0.1 to 20 μm

Features and Benefits

- **Absolute pore size and density:** Provides flow control for liquids moving through the membrane capturing 100 percent of cells larger than pore size
- **Smooth, thin, glass-like surface:** Planar surface makes it ideal for particle identification by microscopy
- **Superior strength:** Tensile strength of 207 bar (> 3000 psi) maintains pore size and density, and will not stretch
- **Low extractables:** Ensures tests will be clean promoting consistent results
- **Low protein binding:** Low binding of < 5 $\mu\text{g}/\text{cm}^2$ minimizes loss of protein analytes
- **Negligible absorption/adsorption:** Maximizes critical solution recovery
- **Available as hydrophilic or hydrophobic:** Allows for a wide range of product applications

Table 3: Nominal Product Characteristics

Thickness	5 - 12 μm
Refractive Indices	Birefringent at 1.584 and 1.625
Water Adsorption (% wt. gain 24-hr immersion)	0.24%
Residual Ash Weight Average	0.92 $\mu\text{g}/\text{cm}^2$
Specific Gravity	0.94-0.97
Autoclavable	Yes
Leachables	Negligible
Wetting Characteristics	Hydrophilic or Hydrophobic
Wetting Agent (hydrophilic)	Polyvinylpyrrolidone (PVP)
Burst Strength Minimum	0.7 bar (10 psi)
Migration of Filter Media	0
Optical Properties	Semi-translucent

Table 4: Performance Characteristics

Pore Size (a) (μm)	Pore Density (b) (pores/ cm^2)	Nominal Thickness (c) (μm)	Min. Bubble Point (d) (psi)	Typical Flow Rates	
				Water (e) ($\text{mL}/\text{min}/\text{cm}^2$)	Air ($\text{L}/\text{min}/\text{cm}^2$)
20	4 x 10 ⁴	3	1	1000	11 (g)
14	5 x 10 ⁴	6	0.2	1400	63.5 (g)
12	1 x 10 ⁵	8	0.4	1250	63.5 (g)
10	1 x 10 ⁵	10	0.5	1150	34.5 (g)
8	1 x 10 ⁵	7	0.7	1000	30 (g)
5	4 x 10 ⁵	10	1.2	700	30 (g)
3	2 x 10 ⁶	9	2	440	37.5 (g)
2	2 x 10 ⁶	10	3	300	16.5 (f)
1	2 x 10 ⁷	11	6	130	20 (f)
0.8	3 x 10 ⁷	9	7	90	18 (f)
0.6	3 x 10 ⁷	9	9	60	7.5 (f)
0.4	1 x 10 ⁸	10	12	33	7.5 (f)
0.2	3 x 10 ⁸	10	20	10	3 (f)
0.1	4 x 10 ⁸	6	30	2.5	1.5 (f)
0.08	4 x 10 ⁸	6	38	0.6	0.75 (f)
0.05	6 x 10 ⁸	6	50	0.4	0.37 (f)
0.03	6 x 10 ⁸	6	NA	0.2	0.075 (f)
0.01	6 x 10 ⁸	6	NA	0.1	0.0075 (f)

(a) Tolerance + 0%, -20%

(b) Tolerance + / - 15%

(c) Tolerance + / - 10%

(d) Measured using Isopropanol (IPA)

(e) Initial flow rates using prefiltered water at 10 psid (0.7 kg/cm²)

(f) Initial flow rates using prefiltered air at 10 psid (0.7 kg/cm²)

(g) Initial flow rates using prefiltered air at 5 psi (0.35 kg/cm²)

FILTRATION MEMBRANES

Precise Pore Size and Density

GVS Life Sciences PCTE Membrane is produced through a two-step, proprietary manufacturing process that employs high quality standards. This unique process allows for increased control over pore size and density to ensure the physical properties of each membrane precisely fit your specifications.

The resulting membrane is a thin, translucent and microporous polycarbonate film with a smooth, flat surface. All particles larger than the pore size are captured on its surface. It is ideal for use when collecting samples for blood assays or for high-purity and general filtration.

Flexibility for OEM Requirements

GVS Life Sciences PCTE Membrane is available in rolls measuring 0.66 - 50 cm (1.0 to 20 inches) wide, or sheets, cut disks, capsules and cartridges, with pore sizes ranging from 0.01 μm to 20 μm . This membrane is also available with a black surface for special use in staining applications, or in its natural hydrophobic state for use in chemotaxis. All varieties of the membrane can be custom manufactured to meet your specific membrane size, pore size and pore densities requirements. Tables 2, 3 and 4 show the product and performance characteristics of the GVS Life Sciences PCTE Membrane.

Ordering information: Polycarbonate Track Etched (PCTE) AOX Membrane

	Dimensions Packaging	25mm 100/pk	47mm 100/pk
Pore sizes	0.4 μm	3026431	1215071

Ordering information: Polycarbonate Track Etched (PCTE) Hydrophilic Black Membrane

	Dimensions Packaging	13mm 100/pk	25mm 100/pk	47mm 100/pk	293 20/pk	203x254mm 100/pk
Pore sizes	0.1 μm	1215311	1215315	1221503		3048982
	0.2 μm	1215185	1215609	1213889	3027176	
	0.4 μm	1215142	1212790	1214567		1227213
	0.6 μm	1222025	1215290	1215198	3048300	3054144*
	0.8 μm	1215236	1215138	1222028	3022140	
	1 μm	1221181	1215161	1222035		
	2 μm		1215297		3033301	
	3 μm		1222452	3032159	3033302	
	5 μm	1221286	1215188	1221230		
	8 μm		1229540			1236363

*100/pk

Ordering information: Polycarbonate Track Etched (PCTE) Hydrophilic Membrane (sheets/rolls)

	Dimensions Packaging	19x42mm 100/pk	25x80mm 50/pk	203x254mm 30/pk	300x3000mm 1/pk
Pore sizes	0.01 μm			1215116	1225184
	0.03 μm			1227264	1239558
	0.05 μm			1215271	3027177
	0.08 μm				
	0.1 μm			1215117	1239556
	0.2 μm			1215118	1239557
	0.4 μm			1215274	
	0.6 μm			1222027	3034261
	0.8 μm			1222030	
	1 μm		1268126	1221429	1267667
	2 μm			1221232	3034567
	3 μm	3019515		1215275	3002536
	5 μm	1221295	1215041	1222080	1264835
	8 μm	1220867	1220686	1222085	3033093
	10 μm			1220823	3033092
	12 μm				1235494
20 μm			1221231		

3.1 Membranes for Filtration

Ordering information: Polycarbonate Track Etched (PCTE) PVDF Hydrophobic Membrane

Dimensions Packaging	13mm 100/pk	25mm 100/pk	47mm 100/pk	90mm 30/pk	203x254mm 100/pk	25x80 50/pk
0.1 μm	1221504	1215059			1232919	
0.2 μm		1222017	1222018		1223036	
0.4 μm		1220835	1215073		1233373	
0.8 μm		1222032				
1.0 μm		1222037	1222038		1224067	
3.0 μm	1215050	1221871	1222077		1228132	1221296
5.0 μm	1215051	1221746	1222081	1222082	1225120	1221331
8.0 μm	1215052	1221293	1215148	1222086	1225783	1215042
10.0 μm	1215053	1222089	1220941		1234298	1215043
12.0 μm	1215055	1221300				1215044
14.0 μm	1221297					

Ordering information: Polycarbonate Track Etched (PCTE) Hydrophilic Membrane (disks)

Dimensions Packaging	13mm 100/pk	19mm 100/pk	25mm 100/pk	37mm 100/pk	47mm 100/pk
0.01 μm	1215046	1227352	1215321		1215068
0.03 μm	1215047		1215057		1215069
0.05 μm	1215048	1221229	1220868		1215070
0.08 μm	1222092	1220668	1215058		1222093
0.1 μm	1215605	1215056	1215606		1215608
0.2 μm	1215610	1220694	1215611		1215612
0.4 μm	1215613	1215147	1215614	1215615	1215617
0.6 μm	1215618		1215619		1215620
0.8 μm	1215621		1215622	1215623	1215624
1 μm	1215625	1227203	1215627	1221302	1215628
2 μm	1215985		1215062		1215629
3 μm	1215049		1215063		1215036
5 μm	1215630		1215631		1215632
8 μm	1215633	3013894	1215634		1215637
10 μm	1221009		1215638		1212661
12 μm	1215054		1215984		3027598
14 μm	1222063	3013893	1222064		1215077
20 μm	1222072		1222073		1215078


FILTRATION MEMBRANES

Ordering information: Polycarbonate Track Etched (PCTE) Hydrophilic Membrane (disks)

Dimensions Packaging	62mm 100/pk	76mm 100/pk	90mm 30/pk	142mm 20/pk	293mm 20/pk
0.01 μm		3032862	1220988		
0.03 μm			1220987		
0.05 μm		1221291	1221227	1221290	1222091
0.08 μm			1222094	1222095	1222096
0.1 μm		1220970	1215150	1215304	1215219
0.2 μm	3038824	1220891	1215151	1215215	1215385
0.4 μm	3023783	1228342	1215303	1215152	1215317
0.6 μm		1224680	1222026	1221485	1220861
0.8 μm		1225894	1215194	1215309	
1 μm		1220860	1215153	1216611	1215145
2 μm			1222070	1222071	1221005
3 μm	3023784	3013824	1222074		1222075
5 μm	3023785	3013825	1221004	1215388	1221329
8 μm		3034848	1215403	1215201	1222084
10 μm		1267014	1222482	1221292	1222088
12 μm		1264834	1239192		
14 μm			1222479		

3.1 Membranes for Filtration

3.1.11 - Polyester Track Etched (PETE) Membrane

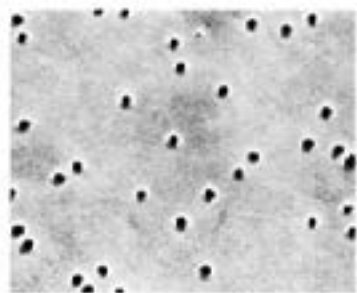

Description and Use

GVS Life Sciences PES is available in rolls up to 31 inches (0.66 - 79cm) wide, as well as sheets, cut disks, capsules, pleat packs and cartridges that can be customized to meet your application requirements. Because the GVS Life Sciences PES is manufactured on-site, all customization can be done easily and cost-effectively.

Features and Benefits

- **Naturally hydrophilic**
- **High range of chemical compatibility:** Resistant to a wide range of chemicals
- **Smooth, thin, glass-like surface:** Captures all particles on the surface larger than the pore size
- **Precise pore size and density:** Provides flow control for liquids moving through the membrane capturing 100 percent of cells larger than pore size
- **Low extractables:** Ensures tests will be clean promoting consistent results
- **Low protein binding:** Low binding prohibits absorption of cells being filtered
- **Negligible adsorption/adsorption:** Maximizes critical solution recovery

Table 2: Product Characteristics

Sterilization	Gamma Irradiation or Ethylene Oxide (EtO)
USP Class VI Testing	Passed
Thickness	10 - 20 μm
Extractables	Low
BSA Protein Binding	< 5 $\mu\text{g}/\text{cm}^2$
Maximum Operating Temperature	284°F (140°C)
Sealing Compatibility	Ultrasonic, Heat, Radio Frequency and Insert Molding
Pore Size Range	0.2 to 10 μm

Typical Applications

- General filtration
- Removal of red blood cells from plasma
- Flow control of reagents through assay
- Precise filtration and prefiltration

Flexibility for OEM Requirements

GVS Life Sciences PETE Membrane is available in rolls measuring 25mm to 50.8cm (1.0 to 20 inches) wide, or sheets, cut disks, capsules and cartridges, with pore sizes ranging from 0,2 μm to 10 μm . The membrane can be custom manufactured to meet your specific membrane size, pore size and pore density requirements.

Product Uniformity and High Sensitivity Maximize Performance

GVS Life Sciences PETE Membrane is produced through a two-step proprietary manufacturing process that employs high quality standards throughout production. This unique process allows for increased control over pore size and pore density to ensure the physical properties of each membrane precisely fit customer specifications.

The resulting membrane is a thin, translucent, microporous polyester film with a smooth, flat surface containing pores of controlled diameter and number. The membrane has better solvent resistance than polycarbonate and captures all particles smaller than the precisely controlled pore size on its surface. GVS Life Sciences PETE is ideal for use when analyzing blood samples or particle analysis by microscopy.

Table 3: Nominal Product Characteristics

Water Adsorption (% wt. gain 24-hr immersion)	0.24%
Residual Ash Weight Average	0.92 $\mu\text{g}/\text{cm}^2$
Specific Gravity	0.94-0.97
Autoclavable	Yes
Leachables	Negligible
Wetting Characteristics	Naturally Hydrophilic
Burst Strength Minimum	0.7 bar (10 psi)
Migration of Filter Media	0
Optical Properties	Semi-translucent

Table 4: Performance Characteristics

Pore Size (a) (μm)	Pore Density (b) (pores/ cm^2)	Nominal Thickness (c) (μm)	Min. Bubble Point (d) (psi)	Typical Flow Rates	
				Water (e) (mL/min/ cm^2)	Air (L/min/ cm^2)
10	1 x 10 ⁵	9	0.5	1150	34.5 (g)
8	1 x 10 ⁵	7	0.7	1000	30 (g)
5	4 x 10 ⁵	10	1.2	700	30 (g)
3	2 x 10 ⁶	9	2	440	37.5 (g)
2	2 x 10 ⁶	10	3	300	16.5 (f)
1	2 x 10 ⁷	11	6	130	20 (f)
0.8	3 x 10 ⁷	9	7	90	18 (f)
0.6	3 x 10 ⁷	9	9	60	7.5 (f)
0.4	1 x 10 ⁸	10	12	33	7.5 (f)
0.2	3 x 10 ⁸	10	20	10	3 (f)

- (a) Tolerance + 0%, -20%
- (b) Tolerance + / - 15%
- (c) Tolerance + / - 10%
- (d) Measured using Isopropanol (IPA)
- (e) Initial flow rates using prefiltered water at 10 psid (0.7 kg/ cm^2)
- (f) Initial flow rates using prefiltered air at 10 psid (0.7 kg/ cm^2)
- (g) Initial flow rates using prefiltered air at 5 psi (0.35 kg/ cm^2)

FILTRATION MEMBRANES

Ordering information: Polyester Track Etched (PETE) Membrane

Dimensions Packaging	13mm 100/pk	25mm 100/pk	47mm 100/pk	90mm 30/pk	142mm* 20/pk	293mm 20/pk	203x254mm 30/pk
0.2 μm	1220969	1221383	1215288	1222240	1221385		1220886
0.4 μm	1221387	1221388	1215373	1220702	1221389		1222242
0.8 μm	1221397	1221398	1215374	1221399		1221401	1222246
1.0 μm	1215379	1215308	1220871	1221402	1222248	1222249	1221334
2.0 μm	1221403	1221404	1221405	1221406			1222251
3.0 μm	1221409	1221410	1215367	1222253	1221411	1221412	1222254
5.0 μm	1215324	1221413	1215183	1221414	1221415	1221416	1222256
8.0 μm	1221417	1221418	1221419	1221420			1222258
10.0 μm	1221423	1220827	1215173	1221424		1221426	1222260

*Bulk packaging available

3.1.12 - Drain Disc

Description and Use

The polyester spun-bonded "drain" type disc prevents "pore blinding" or blockage of the capillary pores in screen membranes resulting in higher flow rates and increased throughputs. The new drain disc also increases flow and capture ability by lifting off of screen supports and exposing all the pores. This ensures efficient performance when placed between two filters in a serial filtration stack too. The spacers prevent air locking of the downstream screen, or function as filters by binding a percentage of pores in the upstream filter .

The spacer may be sized to fit within the diameter of the O-ring in the filter holder. For example , use a 42 mm spacer under a 47 mm filter.

Features and Benefits


- Frequently used with PCTE (Polycarbonatel and PETE (Polyester) membranes to increase flow
- Spacer between stacked membranes

Ordering information: Drain disc

Product Code	Description
1215141	FLT,Drain Disc,25mm,100/pk
1215163	FLT,Drain Disc,47mm 100/pk
1215218	FLT,Drain Disc, 13mm,100/pk
1215500	FLT,Drain Disc,42mm,100/pk
1215522	FLT,Drain Disc, 124mm,25/pk
1215534	FLT,Drain Disc,257mm,25/pk
1221182	FLT,Drain Disc,90mm, 25/pk
3007164	FLT,Drain Disc,293mm, 25/pk
3008749	FLT,Drain Disc,275mm,25/pk
3014223	FLT,Drain Disc,30x30cm,25/pk
3014503	FLT,Drain Disc,50x55mm, 100/pk
1238010	FLT,Drain Disc,37mm,100/pk
Clearing Class Slides	
1215730	Clearing Glass Slides for Track etched membranes 100/pk

3.1 Membranes for Filtration

3.1.13 - Silver Membrane

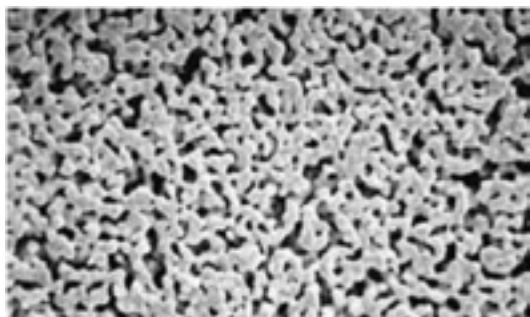

GVS Life Sciences silver membranes are composed entirely of 99.97% pure metallic silver. They combine excellent chemical resistance and high-temperature characteristics with a retention range of 0.2 μ m to 5 μ m. Although silver membranes are used in a large number and variety of applications, their unique chemical and thermal stability is especially valuable for those applications involving aggressive fluids and/or high temperatures. They are ideal collection media for analysis of crystalline silica by X-ray diffraction and for analysis of organic materials by other instrumentation techniques, such as the analysis of Polyaromatic Hydrocarbons (PAH) and Total Organic Carbons (TOC).

Most applications of silver membranes can be divided into two broad categories: clarification/polishing and analytical. In the use of the silver membrane, orientation of the membrane can be important. There is a distinct difference in surface

characteristics of the silver membrane, with one side appearing shinier than the other. For scanning electron microscopy, use the shinier side. For all other filtration applications and analytical work, including X-ray diffraction, the duller of the two sides should be used.

Features and Benefits

- 99.97% pure silver
- High temperature resistance
- Tolerant of chemically or aggressive fluids
- Economical - can be cleaned & reused
- No dissolution or migration of the filter
- Smooth surface for particle capture and easy observation
- Intrinsic bacteriostatic property

Applications

- Airborne asbestos fibers by X-ray diffraction
- Dissolved Organic Carbons (DOC)
- Analysis of airborne silica in foundries, glass plants, quarries, mines, ceramic
- X-ray diffraction applications
- Coke oven emissions analysis
- Carbon and carbon black
- Fly ash - high temperature
- Cleaning liquids, coolants, dopant gases and photoresists

Ordering information: Silver Membrane

	Dimensions Packaging	13mm 100/pk	25mm 50/pk	37mm 25/pk	47mm 25/pk
Pore sizes	0.22 μ m		1145336	1145342	1145348
	0.45 μ m		1145335	1145341	1145347
	0.8 μ m	1145328	1145334		1145346
	1.2 μ m		1145333		1145345
	3.0 μ m		1145332		
	5.0 μ m		1145331		1145337

3.2 Filter Papers

Low Ash / Medium-Slow



Weight (g/m²)

85

Thickness (μm)

180

Retention (μm)

7

Ash content %

< 0,06

Filtration Time for Distilled Water (secs)

60

Diameter (mm)

90, 110, 125, and 150 mm

Wet Bursting Strength (kPa)

> 5

General Application

Qualitative analysis

General filtration

Soil analysis

Air monitoring

Ordering information

Product Code	Description
FP090DMS02QALC01	Qualitative Papers - Low Ash - Medium-Slow, 7.00 μm retention, Ø90 mm, 180 μm thickness
FP110DMS02QALC01	Qualitative Papers - Low Ash - Medium-Slow, 7.00 μm retention, Ø110 mm, 180 μm thickness
FP125DMS02QALC01	Qualitative Papers - Low Ash - Medium-Slow, 7.00 μm retention, Ø125 mm, 180 μm thickness
FP150DMS02QALC01	Qualitative Papers - Low Ash - Medium-Slow, 7.00 μm retention, Ø150 mm, 180 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm - **Box Weight:** Ø90 / Ø110: 110 g - **Quantity / Box:** 100 Units

Ø125: 13,5 x 13,7 x 30 cm

Ø125: 145 g

Ø150: 15,7 x 15,7 x 30 cm

Ø150: 180 g

Detailed Application: Medium-high retention and medium-slow filtration speed. Filtration of fine precipitates; Monitoring of specific contaminants in environment and soil tests.

Low Ash / Medium-Slow/Thick



Weight (g/m²)

200

Thickness (μm)

320

Retention (μm)

5

Ash content %

< 0,06

Filtration Time for Distilled Water (secs)

55

Diameter (mm)

90, 110, 125, and 150 mm

Wet Bursting Strength (kPa)

> 15

General Application

Qualitative analysis

Buchner funnels

High absorbency

Ordering information

Product Code	Description
FP090DMS03QLTC01	Qualitative Papers - Low Ash - Medium-Slow, 5.0 μm retention, Ø90 mm, 320 μm thickness
FP110DMS03QLTC01	Qualitative Papers - Low Ash - Medium-Slow, 5.0 μm retention, Ø110 mm, 320 μm thickness
FP125DMS03QLTC01	Qualitative Papers - Low Ash - Medium-Slow, 5.0 μm retention, Ø125 mm, 320 μm thickness
FP150DMS03QLTC01	Qualitative Papers - Low Ash - Medium-Slow, 5.0 μm retention, Ø150 mm, 320 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm - **Box Weight:** Ø90 / Ø110: 110 g - **Quantity / Box:** 100 Units

Ø125: 13,5 x 13,7 x 30 cm

Ø125: 145 g

Ø150: 15,7 x 15,7 x 30 cm

Ø150: 180 g

Detailed Application: Detailed Application: Medium-high retention and medium-slow filtration speed, with double thickness. Suitable for Buchner funnels. Sample carrier paper for loading greater quantities of solutes due to the high absorbency.

FILTRATION MEMBRANES



Weight (g/m²)

85 - 170

Thickness

170

Retention (μm)

3

Ash content %

< 0,06

Filtration Time for Distilled Water (secs)

300

Diameter (mm)

90, 110, 125, and 150 mm

Wet Bursting Strength (kPa)

> 20

General Application

Qualitative analysis

Clarifying liquids

Water analysis

Soil analysis

Ordering information

Product Code	Description
FP090DXS05QALC01	Qualitative Papers - Low Ash - Very Slow, 3.0 μm retention, Ø90 mm, 170 μm thickness
FP110DXS05QALC01	Qualitative Papers - Low Ash - Very Slow, 3.0 μm retention, Ø110 mm, 170 μm thickness
FP125DXS05QALC01	Qualitative Papers - Low Ash - Very Slow, 3.0 μm retention, Ø125 mm, 170 μm thickness
FP150DXS05QALC01	Qualitative Papers - Low Ash - Very Slow, 3.0 μm retention, Ø150 mm, 170 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm · **Box Weight:** Ø90 / Ø110: 110 g · **Quantity / Box:** 100 Units

Ø125: 13,5 x 13,7 x 30 cm

Ø125: 145 g

Ø150: 15,7 x 15,7 x 30 cm

Ø150: 180 g

Detailed Application: Maximum particle retention. Slow flow rate. High retention of fine particles in chemical analysis. Clarification of cloudy suspensions (wine); Water and soil analysis.

3.2.2 - Quantitative Papers

Ashless / Fast



Weight (g/m²)

80

Thickness (μm)

190

Retention (μm)

20

Ash content %

< 0,007

Filtration Time for Distilled Water (secs)

21

Diameter (mm)

90, 110, 125, and 150 mm

Wet Bursting Strength (kPa)

40

General Application

Quantitative analysis

Air monitorin

Ordering information

Product Code	Description
FP090DFA41QANC01	Quantitative Papers - Ashless - Fast, 20.00 μm retention, Ø90 mm, 190 μm thickness
FP110DFA41QANC01	Quantitative Papers - Ashless - Fast, 20.00 μm retention, Ø110 mm, 190 μm thickness
FP125DFA41QANC01	Quantitative Papers - Ashless - Fast, 20.00 μm retention, Ø125 mm, 190 μm thickness
FP150DFA41QANC01	Quantitative Papers - Ashless - Fast, 20.00 μm retention, Ø150 mm, 190 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm · **Box Weight:** Ø90 / Ø110: 110 g · **Quantity / Box:** 100 Units

Ø125: 13,5 x 13,7 x 30 cm

Ø125: 145 g

Ø150: 15,7 x 15,7 x 30 cm

Ø150: 180 g

Detailed Application: Very fast ashless filter paper. Analytical procedures with large particles or gelatinous precipitates (iron or aluminum hydroxides). Air pollution analysis to determinate gaseous compounds.

3.2 Filter Papers

Ashless / Medium



Weight (g/m²)

80

Thickness (μm)

180

Retention (μm)

14

Ash content %

< 0,007

Filtration Time for Distilled Water (secs)

50

Diameter (mm)

90, 110, 125, and 150 mm

Wet Bursting Strength (kPa)

30

General Application

Quantitative analysis

Food analysis

Soil analysis

Industrial analysis

Ordering information

Product Code	Description
FP090DME43QANC01	Quantitative Papers - Ashless - Medium, 14.00 μm retention, Ø90 mm, 180 μm thickness
FP110DME43QANC01	Quantitative Papers - Ashless - Medium, 14.00 μm retention, Ø110 mm, 180 μm thickness
FP125DME43QANC01	Quantitative Papers - Ashless - Medium, 14.00 μm retention, Ø125 mm, 180 μm thickness
FP150DME43QANC01	Quantitative Papers - Ashless - Medium, 14.00 μm retention, Ø150 mm, 180 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm - **Box Weight:** Ø90 / Ø110: 110 g - **Quantity / Box:** 100 Units

Ø125: 13,5 x 13,7 x 30 cm

Ø125: 145 g

Ø150: 15,7 x 15,7 x 30 cm

Ø150: 180 g

Detailed Application: Medium retention and fast flow. Foodstuff and soil analysis. Air pollution monitoring. Analysis in mining, construction and steel industries.

Ashless / Medium-Slow



Weight (g/m²)

80

Thickness (μm)

14

Retention (μm)

7

Ash content %

< 0,007

Filtration Time for Distilled Water (secs)

50

Diameter (mm)

90, 110, 125, and 150 mm

Wet Bursting Strength (kPa)

36

General Application

Quantitative analysis

Gravimetric analysis

Soil analysis

Air monitoring

Ordering information

Product Code	Description
FP090DMS40QANC01	Quantitative Papers - Ashless - Medium-Slow, 7.00 μm retention Ø90 mm, 170 μm thickness
FP110DMS40QANC01	Quantitative Papers - Ashless - Medium-Slow, 7.00 μm retention, Ø110 mm, 170 μm thickness
FP125DMS40QANC01	Quantitative Papers - Ashless - Medium-Slow, 7.00 μm retention, Ø125 mm, 170 μm thickness
FP150DMS40QANC01	Quantitative Papers - Ashless - Medium-Slow, 7.00 μm retention, Ø150 mm, 170 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm - **Box Weight:** Ø90 / Ø110: 110 g - **Quantity / Box:** 100 Units

Ø125: 13,5 x 13,7 x 30 cm

Ø125: 145 g

Ø150: 15,7 x 15,7 x 30 cm

Ø150: 180 g

Detailed Application: Medium speed and retention. Analysis of components in cements, clays, iron and steel products. Soil analysis. Sediments in milk. Filtration of solutions prior to atomic absorption spectrophotometry; High purity filter in atmosphere analysis.

FILTRATION MEMBRANES



Weight (g/m²)
80
Thickness (μm)
160
Retention (μm)
2
Ash content %
< 0,007
Filtration Time for Distilled Water (secs)
292

Diameter (mm)
90, 110, 125, and 150 mm
Wet Bursting Strength (kPa)
52
General Application
Quantitative analysis

Ordering information

Product Code	Description
FP090DSL44QANC01	Quantitative Papers - Ashless - Slow, 2.00 μm retention Ø90 mm, 160 μm thickness
FP110DSL44QANC01	Quantitative Papers - Ashless - Slow, 2.00 μm retention, Ø110 mm, 160 μm thickness
FP125DSL44QANC01	Quantitative Papers - Ashless - Slow, 2.00 μm retention, Ø125 mm, 160 μm thickness
FP150DSL44QANC01	Quantitative Papers - Ashless - Slow, 2.00 μm retention, Ø150 mm, 160 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm - **Box Weight:** Ø90 / Ø110: 110 g - **Quantity / Box:** 100 Units
 Ø125: 13,5 x 13,7 x 30 cm Ø125: 145 g
 Ø150: 15,7 x 15,7 x 30 cm Ø150: 180 g

Detailed Application: Slow flow rate. Very high retention. Very fine crystalline precipitates.

Ashless / Very Slow



Weight (g/m²)
100
Thickness (μm)
160
Retention (μm)
2
Ash content %
< 0,007
Filtration Time for Distilled Water (secs)
292

Diameter (mm)
90, 110, 125, and 150 mm
Wet Bursting Strength (kPa)
52
General Application
Quantitative analysis
Critical gravimetric analysis.

Ordering information

Product Code	Description
FP090DXS42QANC01	Quantitative Papers - Ashless - Very Slow, 2.00 μm retention Ø90 mm, 160 μm thickness
FP110DXS42QANC01	Quantitative Papers - Ashless - Very Slow, 2.00 μm retention, Ø11, 160 μm thickness
FP125DXS42QANC01	Quantitative Papers - Ashless - Very Slow, 2.00 μm retentionØ125 mm, 160 μm thickness
FP150DXS42QANC01	Quantitative Papers - Ashless - Very Slow, 2.00 μm retentionØ150 mm, 160 μm thickness

Box Dimension: Ø90 / Ø110: 11,7 x 12,0 x 30 cm - **Box Weight:** Ø90 / Ø110: 110 g - **Quantity / Box:** 100 Units
 Ø125: 13,5 x 13,7 x 30 cm Ø125: 145 g
 Ø150: 15,7 x 15,7 x 30 cm Ø150: 180 g

Detailed Application: Highest retention and very slow flow. Extremely difficult filtrations. Analytical precipitates: barium sulphate, metastannic acid and finely precipitated calcium carbonate.

3.2 Filter Papers

3.2.3 Glass Microfiber

Fast



Weight (g/m²)

52

Thickness (µm)

0,26

Retention (µm)

1,6

Filtration Time for Distilled Water (secs)

40

Air Resistance (mbar)

36

Diameter (mm)

47, 55, 90, 110 and 125 mm

Dry Bursting Strength (kPa)

20

Max. Operating Temperature (°C)

500

General Application

General purpose laboratory filtration

Food analysis

Water analysis

Ordering information

Product Code	Description
FP047DFAFAGLFC01	Glass Microfiber - Fast, 1.60 µm retention, Ø47 mm, 0,26 mm thickness
FP055DFAFAGLFC01	Glass Microfiber - Fast, 1.60 µm retention Ø55 mm, 0,26 mm thickness
FP090DFAFAGLFC01	Glass Microfiber - Fast, 1.60 µm retention Ø90 mm, 0,26 mm thickness
FP110DFAFAGLFC01	Glass Microfiber - Fast, 1.60 µm retention Ø110 mm, 0,26 mm thickness
FP125DFAFAGLFC01	Glass Microfiber - Fast, 1.60 µm retention Ø125 mm, 0,26 mm thickness

Box Dimension: Ø47 / Ø55: 13,5 x 13,7 x 30 cm - **Box Weight:** Ø47: 102 g - **Quantity / Box:** 100 Units

Ø90: 11,7 x 12,0 x 30 cm

Ø110 / Ø125: 15,7 x 15,7 x 30 cm

Ø55: 105 g

Ø 90: 108 g

Ø110: 164 g

Ø125: 180 g

Detailed Application: Fine retention with fast flow. Water pollution and air pollution monitoring. Foodstuff analyses. Protein filtration. Filtration of water, algae and bacteria cultures.

Medium



Weight (g/m²)

53

Thickness (µm)

0,26

Retention (µm)

1,2

Filtration Time for Distilled Water (secs)

66

Air Resistance (mbar)

51

Diameter (mm)

47, 55, 90, 110 and 125 mm

Dry Bursting Strength (kPa)

20

Max. Operating Temperature (°C)

500

General Application

Cell culture

Water analysis

Ordering information

Product Code	Description
FP047DMEFCGLFC01	Glass Microfiber - Medium, 1.20 µm retention, Ø47 mm, 0,26 mm thickness
FP055DMEFCGLFC01	Glass Microfiber - Medium, 1.20 µm retention, Ø55 mm, 0,26 mm thickness
FP090DMEFCGLFC01	Glass Microfiber - Medium, 1.20 µm retention, Ø90 mm, 0,26 mm thickness
FP110DMEFCGLFC01	Glass Microfiber - Medium, 1.20 µm retention, Ø110 mm, 0,26 mm thickness
FP125DMEFCGLFC01	Glass Microfiber - Medium, 1.20 µm retention, Ø125 mm, 0,26 mm thickness

Box Dimension: Ø47 / Ø55: 13,5 x 13,7 x 30 cm - **Box Weight:** Ø47: 102 g - **Quantity / Box:** 100 Units

Ø90: 11,7 x 12,0 x 30 cm

Ø110 / Ø125: 15,7 x 15,7 x 30 cm

Ø55: 105 g

Ø 90: 108 g

Ø110: 164 g


Ø125: 180 g

Detailed Application: Fine retention with medium flow. Suspended solids analysis in water and industrial wastes; Cell harvesting. Aqueous clarification and monitoring.

FILTRATION MEMBRANES

3.3 Glass Fiber Filters with Binder

Glass Fiber Filters with Binder

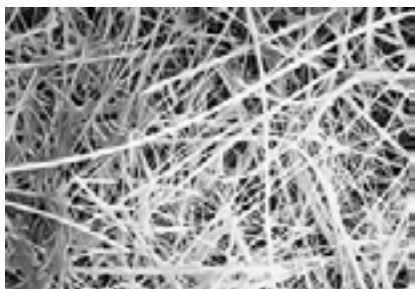

acrylic resin. This bonding produces a filter that reduces media migration and has the strength required for high-volume aqueous filtrations. Glass Fiber membranes with a binder are usually recommended for filtrations of long duration under pressure. Glass Fiber membranes without binders are designed for solvent filtration or gravimetric analysis to avoid binder extractables. Filters without binders are recommended for analytical and gravimetric determinations.

Features and Benefits

- Acrylic binder
- High dirt holding capacity
- Biologically inert
- Bonding reduces media migration

Description and Use

GVS Life Sciences Glass Fiber membranes are biologically inert, autoclavable and highly resistant to oxidizing agents and weak acids. Glass fiber can be used to extend the life of a final filter as a prefilter or they can be used alone for low cost sample clarification. GVS Life Sciences Glass Fiber membranes with binders are composed of borosilicate glass fibers woven into a porous matrix and bonded by an

Ordering information: Glass Fiber Filters

Dimensions Packaging	10mm 100/pk	13mm 100/pk	22 mm 100/pk	25 mm 100/pk	42 mm 100/pk	47mm 100/pk	75 mm 25/pk	90 mm 25/pk
Pore sizes 0.5 μm			1215543	1215544	1215547	1215548	1215549	1215550
1.0 μm (G20)		1215557		1215559	1215561	1215562	1215563	1215564
1.0 μm (G25)	1215570	1215571	1215572	1215573	1215576	1215577	1215578	1215579

Dimensions Packaging	124 mm 25/pk	127 mm 25/pk	142 mm 25/pk	257 mm 25/pk	293 mm 25/pk	24x24 cm 10/pk
Pore sizes 0.5 μm	1215551		1215553	1215554	1215555	1266844
1.0 μm (G20)	1215565	1215566	1215567	1215568	1215569	1268603
1.0 μm (G25)	1215580		1215582	1215583	1215584	

Ordering information: Glass Fiber Filters Binderless

Dimensions Packaging	6 mm 500/pk	7 mm 500/pk	10 mm 500/pk	25 mm 100/pk	37 mm 500/pk	47 mm 100/pk	65mm 100/pk	82 mm 100/pk
Pore sizes 0.7 μm		3029939		1215162		1215540		
1.0 μm	3003072		1214912	1213325*	1215588	1215589*	1221996	1214974

Dimensions Packaging	85 mm 100/pk	90 mm 25/pk	102 mm 100/pk	110 mm 25/pk	142 mm 25/pk	257 mm 100/pk	293 mm 25/pk
Pore sizes 0.7 μm		1215541			1215542		
1.0 μm	3015810	1225509 1212763**	1214671	3034573	3034574	1220678	1220887

*500/pk **100/pk

TRANSFER MEMBRANES

GVS

LIFE SCIENCES



TRANSFER MEMBRANES

4. TRANSFER MEMBRANES

4.1 Neutral Nylon



Description and Use

GVS Life Sciences Neutral Nylon Transfer Membrane is a pure polymer impregnated by an inert polyester web. It is naturally hydrophilic and optimized for protein binding and for high, reproducible binding of nucleic acids.

Features and Benefits

- **Supported:** Has added strength and durability preventing distortion or contamination in multiple reproblings
- **High binding capacity:** With a nucleic acid binding capacity of approximately 350 $\mu\text{g}/\text{cm}^2$, Magna Nylon - Transfer Membrane can bind a wide range of fragment sizes, increasing the efficiency of transfers
- **Hydrophilic:** Eliminates the need for wetting agents that can potentially interfere with biological processes
- **Lot-to-lot consistency:** Quality checks ensure lot-to-lot consistency, both down and across the polyester web, for dependable results every time

Table 2: Product Characteristics

USP Class VI testing	Passed
Thickness	65 - 125 μm
Extractables	< 0.2% (< 0.0015 mg/cm ²)
Nucleic Acid Binding	350 $\mu\text{g}/\text{cm}^2$
Maximum Operating Temperature	356°F (180°C)
Sealing Compatibility	Ultrasonics, Heat, Radio Frequency and Insert Molding
Pore Size Range	0.22 to 0.45 μm

Typical Applications

- Southern transfers
- Northern transfers
- Protein binding
- Microarrays
- Macroarrays
- Dot/Slot blots
- Radiolabeled detection systems
- Non-radiolabeled detection systems
- Colony lifts
- Plaque lifts
- Library screening

Reliable Quality, Increased Efficiencies

Produced through a proprietary manufacturing process, GVS Life Sciences Neutral Nylon Transfer Membrane meets rigorous quality standards throughout every step of production. This process generates consistent lot- to-lot binding among the membranes ensuring product uniformity.

This controlled microporous nylon membrane is cast on an inert, internal support web that gives it added dimensional strength and stability to prevent cracking, tearing, curling and breaking. This added strength and durability is essential in protocols that require aggressive handling, such as colony lifts and plaque lifts. In addition to the dimensional strength and durability of GVS Life Sciences Neutral Nylon Transfer Membrane, its retention of macromolecules can also be enhanced using UV cross-linking. This process can be used to maximize the signal retention of nucleic acids and preserve the integrity of DNA or RNA transfers.

The purity and consistency of GVS Life Sciences Neutral Nylon Transfer Membrane, coupled with its added durability and sensitivity, make it an ideal membrane for use in medical research, scientific studies or test confirmations where precise biological pattern replications, such as DNA and RNA transfers, are integral to the success of the procedure.

Flexibility for OEM Requirements

GVS Life Sciences Neutral Nylon Transfer Membranes are available in rolls up to 13 inches (33cm) wide, as well as sheets, cut disk, and pleat packs that can be customized to meet your application and size requirements. Because the membrane is manufactured on-site, all customization is easy and cost-effective.

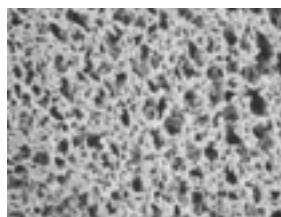


Figure 2: GVS Life Sciences Neutral Nylon Transfer Membrane

Ordering information: Neutral Nylon Transfer Membrane

Dimensions Packaging	82mm 50/pk	85mm 50/pk	132mm* 50/pk	137mm 50/pk	102x133mm 10/pk	115x160mm 10/pk	150x150mm 5/pk	200x200mm 5/pk
Pore sizes								
0.22 μm	1213409	1213410	1213411	1213412	1213422		1213416	1213419
0.45 μm	1213370 1214428*	1213372	1213373 1214509*	1213375 1214245*	1213384	1213391	1213379	1213380 1215310**
	*100/pk		**25/pk					
Dimensions Packaging	220x220mm 5/pk	225x225mm 10/pk	300x300mm 5/pk	300x500mm 5/pk	150x300mm 1/pk	200x300mm 1/pk	300x300mm 1/pk	4.75x11in 10/pk
Pore sizes								
0.22 μm					1213442	1213441	1213405	
0.45 μm	1213382	1224585	1213383	1213395	1213404	1213403	1213364	1214994

TRANSFER MEMBRANES

4.2 Reprobing Charged Nylon



Description and Use

GVS Life Sciences Nylon Reprobing, Charged transfer membrane is an inherently charged nylon membrane, specifically designed to allow for numerous reprobing. GVS Life Sciences Nylon has been tested to provide consistent results through 12 or more reprobing.

The high binding capacity of 450 mg/cm² makes GVS Life Sciences Nylon ideal for all Southern and Northern applications, including alkaline blotting. GVS Life Sciences Nylon is ideally suited for all probes both radioactive and non-radioactive, including chemiluminescent and biotinylated detection systems.

GVS Life Sciences Nylon offers significantly increased binding, maximum “lot-to-lot” consistency, and excellent signal retention. The inherent charge on this nylon membrane along with its hydrophilic nature makes consistent repeatable results a reality for researchers.

After 12 rounds of reprobing, GVS Life Sciences Nylon has a lower background and higher signal.

Features and Benefits

- Supported charged nylon membrane
- Specifically designed for multiple reprobing
- Used for both radiolabelled & non-radiolabelled detection systems
- Can be used for alkaline blotting
- Nucleic acid binding is 450 µg/cm²
-

Typical Applications

- Radiolabelled & non-radiolabelled detection systems
- Northern
- Southern
- Multiple Reprobing
- Alkaline Blotting
- UV Crosslinking

Ordering information: Reprobing Charged Nylon Transfer Membrane

	Dimensions Packaging	82mm 50/pk	82mm 100/pk	85mm 50/pk	132mm 50/pk	132mm 100/pk	137mm 50/pk	137mm 100/pk
Pore sizes	0.45 µm	1226559	1226561	1226560	1226563	1226565	1226562	1226564

	Dimensions Packaging	102x133mm 10/pk	150x150mm 5/pk	150x150mm 25/pk	200x200mm 5/pk	200x200mm 25/pk	220x220mm 5/pk	300x300mm 5/pk
Pore sizes	0.45 µm	1226570	1226566	1226572	1226567	1226573	1226568	1226569

	Dimensions Packaging	300x300mm 25/pk	300x500mm 5/pk	300x500mm 25/pk	150x3000mm 1/pk	200x3000mm 1/pk	300x3000mm 1/pk
Pore sizes	0.45 µm	1226575	1226571	1226574	1226558	1226557	1226556

TRANSFER MEMBRANES

4.3 Pure Nitrocellulose (MCE)



Description and Use

GVS Life Sciences Pure Nitrocellulose Transfer Membrane is the membrane of choice for all protein or immunoblotting applications. The high sensitivity of GVS Life Sciences Pure Nitrocellulose Transfer Membrane ensures excellent results in all transfers, especially in protein blotting.

Features & Benefits

- For procedures that require optimum resolution
- Membrane of choice for protein or immunoblotting applications
- Low background, easily blocked
- BSA binding capacity up to 100 μ g/cm²
- Wets out naturally
- Compatible with all detection systems

Typical Applications

- Westerns
- Protein & immunoblotting
- Northern
- Southern
- Dot/slot blots
- Radiographic, chromogenic and chemiluminescent detection systems

Ordering information: Pure Nitrocellulose Transfer Membrane

Dimensions(mm) Packaging	82 50/pk	82 100/pk	85 50/pk	132 50/pk	137 50/pk	70x84 10/pk	70x100 10/pk	90x120 10/pk	100x100 10/pk	102x133 10/pk	140x160 10/pk	150x150 5/pk
Pore sizes 0.22 μ m	1215459		1215460			1213991	1215394		1213999	1215466		1215463
0.45 μ m	1215472	1214243	1215473	1215474	1215475 1221974*	1213888		1214911	1213314	1215481	1221981	1215476

Dimensions(mm) Packaging	150x150 25/pk	200x200 5/pk	200x200 25/pk	220x220 5/pk	250x250 5/pk	300x300 5/pk	300x500 5/pk	300x500 25/pk	150x3m 1/pk	200x3m 1/pk	300x3m 1/pk
Pore sizes 0.22 μ m		1215464	1215392	1213121		1215465	1215467		1215470	1215469	1215458
0.45 μ m	1215170	1215477	1221976	1215478	1223273	1215480	1215482	1221983	1215484	1215483	1215471

TRANSFER MEMBRANES

4.4 Supported Nitrocellulose (MCE)



Description and Use

GVS Life Sciences Supported Nitrocellulose Transfer Membrane combines the binding characteristics of nitrocellulose membrane with the strength of nylon membrane. In all protocols using nitrocellulose transfer membrane, GVS Life Sciences Supported Nitrocellulose Transfer membrane may be substituted without requiring any changes to the protocol.

Features & Benefits

- Supported for procedures requiring rigorous handling
- Strong - will not curl, bend or crack after baking
- High sensitivities, low backgrounds
- Multiple reprobings
- BSA binding capacity up to 100 μ g/cm²
- Triton Free

Typical Applications

- Northern
- Southern
- Multiple re-hybridizations
- Colony/plaque lifts
- Dot/slot blots
- Radiographic detection systems
- Chemiluminescent detection systems
- Biotinylated detection systems

Ordering information: Supported Nitrocellulose Transfer Membrane

Dimensions(mm) Packaging	82 50/pk	85 50/pk	132 50/pk	137 50/pk	137 100/pk	70x84 10/pk	70x100 10/pk	90x120 10/pk	100x100 10/pk	100x150 10/pk	102x133 10/pk
Pore sizes 0.22 μ m	1212640		1214126					1222287	1214560		1213190
0.45 μ m	1212591 1214247*	1212592	1212594	1215475 1221974**	1214248	1214978	1215319		1213943	1222295	1212600

Dimensions(mm) Packaging	140x160 25/pk	150x150 5/pk	200x200 5 /pk	200x200 25/pk	220x220 5/pk	300x300 5/pk	300x500 5/pk	150x3000 1/pk	200x3000 1/pk	300x3000 1/pk
Pore sizes 0.22 μ m		1212669	1212689			1213195	1213141	1212721	1212690	1212632
0.45 μ m	1214935	1212596	1212597	1214762	1212598	1212599	1212601	1212604	1212602	1212590

* 100/pk ** gridded

TRANSFER MEMBRANES

4.5 Polyvinylidene Fluoride (PVDF)



- **BSA protein binding capacity** : 125 mg/cm²
- **High range of chemical**: Resistant to most commonly used chemicals compatible with chemically aggressive solvents

Typical Applications

- Western blots
- Immunoblotting
- Solid phase assays
- Amino acid or protein analyses

Exceptional Strength and Chemical Compatibility

Exceptional Strength and Chemical Compatibility Produced through a proprietary manufacturing process, GVS Life Sciences PVDF Transfer Membrane meets rigorous quality standards throughout every step of production. This process generates membranes with consistent lot-to-lot protein binding and ensures excellent product uniformity.

GVS Life Sciences PVDF is a naturally hydrophobic, unsupported transfer membrane. It has a high binding capacity, which prevents protein from passing through the membrane, and a low background that provides for an excellent signal-noise ratio. It also has exceptional tensile strength, preventing it from cracking, tearing, breaking or curling.

This membrane also has broad chemical compatibility, which is important when used with common stains such as Amido Black, Colloidal Gold, Coomassie Blue, India Ink and Ponceau-S. GVS Life Sciences PVDF will not degrade, distort or shrink when a high concentration of methanol is used for destaining.

Its exceptional strength, high binding capacity and chemical compatibility make GVS Life Sciences PVDF ideal for use in Western blots, immunoblotting, and solid phase assays and plaque lifts.

Flexibility for OEM Requirements

GVS Life Sciences PVDF Transfer Membrane is available in rolls from 0.6cm - 30cm (12 inches) wide, as well as sheets and cut disks that can be customized to meet your application and size requirements. Because the GVS Life Sciences PVDF Transfer Membrane is manufactured on-site in GVS Life Sciences facilities, all customization is easy and cost-effective.

Description and Use

The GVS Life Sciences PVDF Transfer Membrane is a naturally hydrophobic, unsupported transfer membrane. It has a high binding capacity and low backgrounds and is ideal for use in protein binding applications such as Western blots, solid phase assays and immunoblotting procedures.

Features & Benefits

- **Superior strength**: Can withstand aggressive handling or be used with automated equipment without breaking or tearing
- **Low extractables**: Ensures tests will be clean with consistent results
- **Exceptional sensitivity**: Detects low-level components
- **Hydrophobic**: For high protein binding
- **Lot-to-lot consistency**: Quality checks ensure consistent binding for dependable results every time

Ordering information: PVDF (Polyvinylidene Fluoride) Transfer Membrane

Dimensions(mm) Packaging	70x84 10/pk	90x100 10/pk	90x120 10/pk	100x100 10/pk	100x150 10/pk	100x200 10/pk	150x150 5/pk
Pore sizes 0.22 μm	1214588	1222216	1214423				1215037
0.45 μm	1213992	1214558	1212642	1212644	1212643	1214391	1212636

Dimensions(mm) Packaging	150x150 25/pk	200x200 5/pk	200x200 25/pk	240x240 10/pk	260x500 10/pk	150x3000 1/pk	200x3000 1/pk	300x3000 1/pk
Pore sizes 0.22 μm		1215032				1214842	1214726	1214429
0.45 μm	1222218	1212637	1214802	1214600	1212638	1212781	1212783	1212639

PROTEIN ARRAY



LIFE SCIENCES



PROTEIN ARRAY

5. FAST - protein array solutions



GVS Life Sciences, is proud to reintroduce the entire line of FAST® protein microarray products:

- FAST Slides
- FAST PAK Starter Kits
- FAST Quant® Kits for multiplexed immunoassays
- Full line of Buffers and Accessories
- Scanning, Data Analysis and Development Services

Brought to you by the team who developed and has improved the FAST line of products over the last decade, we welcome your business with products optimized for performance and reliability, using a Quality Management System registered to the ISO 9001 standard.

BACKGROUND

Protein microarrays have begun to fulfill their great potential. They are now recognized as critical tools for proteomics, biomarker research and drug discovery. They allow discovery and analysis of protein expression patterns, posttranslational modifications and protein interactions involved in cell growth and differentiation, environmental and drug responses and disease progression such as in cancer. Cited in more than 150 publications, FAST Slides have proven to be the surface of choice for designing and building protein microarrays. The surface is a proprietary nitrocellulose coating that non-covalently binds proteins maintaining their native structure. Nitrocellulose provides a homogeneous 3-D surface for uniform protein binding and significantly greater binding capacity than 2-D or ultrathin coatings. With sensitivities down to attamoles (10⁻¹⁸ molar) and near quantitative capture over a broad dynamic range of four orders of magnitude, FAST Slides offer unparalleled detection ability, reproducibility and reliability.

FAST PROTEIN ARRAYS ARE IDEAL FOR MANY APPLICATIONS

Protein arrays are now recognized as a key tool for proteomics research. FAST Slides offer unmatched binding capacity, sensitivity and reproducibility ideal for all of your protein array applications. GVS Life Sciences provides a full line of products and services that allow you to apply the FAST Slide technology to any multiplex need.

PROTEIN ARRAYS

The high reproducibility and microporous structure of FAST Slides make them excellent for protein arrays used to diagnose infection and autoimmune diseases and for vaccine development and immunity monitoring. In protein arrays, a purified protein is spotted on the FAST Slide surface and the array is used to detect the presence of antibodies or other binding proteins in clinical or experimental samples.



REVERSE PHASE PROTEIN ARRAYS (RPPAs)

The high binding capacity, sensitivity and reproducibility of FAST Slides make them ideal for Reverse Phase Protein Arrays (RPPAs) used for biomarker discovery and characterization and in clinical trials to monitor drug effectiveness and disease progression.

FAST Slides allow quantitative binding across the broad dynamic range of protein concentration found in complex biological samples, such as cell lysates or tumor aspirates, arrayed directly onto the slides. The expression of specific proteins is detected with antibodies to the biomarkers of interest. GVS Life Sciences's quality-controlled FAST Slide production ensures a consistent surface that allows multiple clinical samples to be interrogated with a variety of antibodies simultaneously.

ANTIBODY ARRAYS or MICRO-SPOT ELISA

FAST Slides are arrayed with multiple antibodies and the presence of specific proteins is detected by applying a complex biological sample to the slide. When used in combination with a standard curve, the unsurpassed binding capacity that retains the native conformation of proteins bound on FAST Slides allows detection and quantitation of multiple proteins in a single sample.

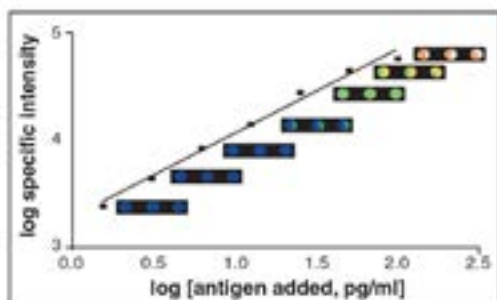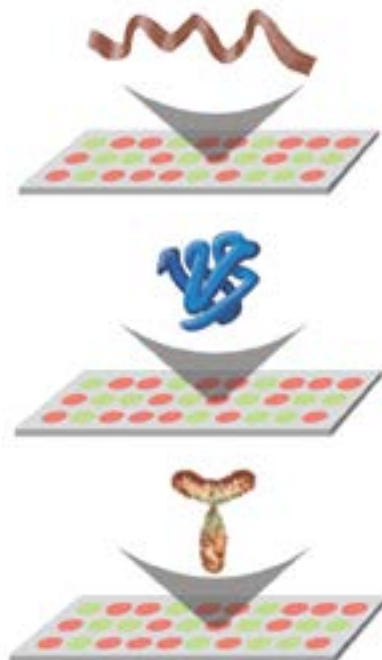

FAST SLIDES PROVIDE AN OPTIMUM IMMUNOASSAY SURFACE

The high binding capacity and broad dynamic range provide the highest sensitivity assuring near-quantitative protein recoveries (94-96%) from complex biological samples such as plasma.


PROTEIN ARRAY

FAST SLIDES OFFER LONG-TERM STABILITY

FAST Slides provide a stabilizing environment such that proteins arrayed with the GVS Life Sciences Protein Arraying Buffer retain their binding characteristics for more than a year. As shown, there is no deterioration in the fluorescent signal on a FAST Quant Slide from 3-12 months in storage. The long-term stability offers the flexibility of being ready whenever you are. For your convenience, GVS Life Sciences offers a number of pre-arrayed FAST Slide kits packaged with the buffers and reagents necessary to accommodate your research needs.



3 Months

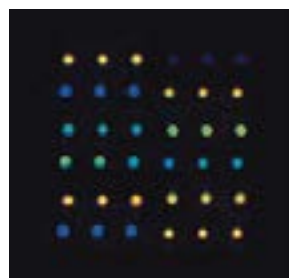

12 Months

FAST SLIDES ARE A BROADLY COMPATIBLE OPEN PLATFORM

FAST Slides are ready for printing (arraying) right out of the box with no activation. They come in five different pad formats compatible with contact and non-contact printing methods and with existing manual arrayers, robots, scanners and all methods of detection including fluorescence, chemiluminescence, colorimetric and isotopic. Their compatibility make FAST Slides easy to set up and use with low costs and minimal start-up time.

Our focus on quality at GVS Life Sciences (registered to the ISO 9001 standard) ensures production of the most consistent surface possible providing the most reproducible results with every slide, every time. The FAST line from GVS Life Sciences provides an unparalleled set of protein analysis products and services for all of your quantitative and non-quantitative multiplexed assays.

FAST Slides

Nitrocellulose-coated glass slides offered in 1-, 2-, 8- and 16-pad formats in 10 or 20 Slide packs.

FAST PAK

Complete FAST Slide kits containing 1-, 2-, 8-, or 16-pad FAST Slides packaged with reaction reagents to create and analyze your own protein microarrays.

FAST Quant® Kits

16-pad FAST Slides pre-arrayed with specific cytokine antibody panels and packaged with reaction buffers and reagents.

1. Human Th1/Th2: panel of antibodies to 9 cytokines associated with the human Th1/Th2 immune response
2. Human Cytokine II: panel of antibodies to 10 commonly requested cytokines
3. Human Angiogenesis: panel of antibodies to 8 cytokines associated with angiogenesis
4. Human Chemokine: antibodies to 9 commonly requested chemokines
5. Human Cytokine 35: panel of antibodies to 35 of the most-requested cytokines



Reaction Buffers

1. Protein Arraying Buffer
2. Protein Array Washing Buffer
3. Protein Array Blocking Buffer

Array Accessories

1. Incubation chambers to fit multi-pad FAST Slides
2. FAST Frame to hold up to four FAST Slides
3. FAST Slide holder (Chip Clip™) for a single FAST Slide
4. MicroCaster™ for printing your own arrays onto FAST Slides
5. Other parts and accessories



Scanning and Data Analysis Service


Slide scanning and data analysis service for FAST Slide or FAST Quant users who do not have access to a fluorescent scanner.

Microarray Assay Development

Contact GVS Life Sciences Customer Care to discuss your protein array development needs.

PROTEIN ARRAY

5.1 FAST Slides - protein array surface

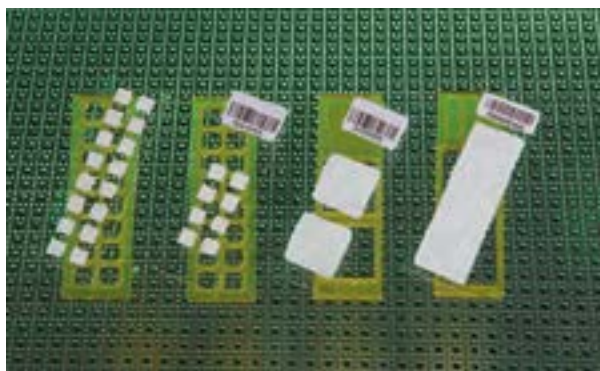

FAST Slides are the premier surface for protein microarray applications. FAST Slides are glass slides coated with a proprietary nitrocellulose polymer. The polymer binds proteins in a noncovalent, irreversible manner and can be probed using the same method as in traditional blots. The three-dimensional surface of a FAST Slide maintains reactivity of proteins and gives excellent reproducible results. It is usable with fluorescent, colorimetric, chemiluminescent or isotopic detection systems and is compatible with microarray scanners and robots. Perhaps the most significant advantage of FAST Slides over modified glass surfaces is that the matrix retains arrayed proteins in near-quantitative fashion. This property translates into antibody arrays with unparalleled sensitivity below 1 pg/ml in antigen concentration. These qualities make FAST Slides the most reliable surface for microarray experiments and provide a high level of confidence.

Array Accessories

- Incubation chambers to fit multi-pad FAST Slides
- Fast Frame to hold up to four FAST Slides
- FAST Slide Holder (Chip Clip™) for a single FAST Slide
- MicroCaster™ for printing your own arrays onto FAST Slides

Features & Benefits

- Superior protein binding capacity
- Highest sensitivity and dynamic range
- Excellent long-term stability of printed proteins
- Compatible with all detection methodologies
- Compatible with commercially available arraying robots

FAST Slides are suitable for many types of protein microarrays including protein arrays, reverse phase protein arrays and micro-spot ELISAs, also known as antibody arrays.

There are tremendous advantages to using FAST Slides for immunoassays over traditional ELISAs including less sample required, better sensitivity, linearity and quantitation.

A major advantage of FAST Slide technology is that hundreds or thousands of antibodies or samples can be screened simultaneously. Compared to other microarray surfaces, FAST Slides provide superior binding properties, allowing quantitative detection of proteins over four orders of magnitude in concentration. Our focus on quality at GVS Life Sciences (registered to the ISO 9001 standard) ensures production of the most consistent surface possible providing the most reproducible results with every slide, every time.

FAST PAK

Complete FAST Slide starter kits containing 1-, 2-, 8- or 16-pad FAST Slides packaged with reaction buffers to create and analyze your own protein microarrays.

Reaction Buffers

- Protein Arraying Buffer
- Protein Array Washing Buffer
- Protein Array Blocking Buffer

Ordering information: FAST Slides

Product Code	Description	Specification
10484182	FAST SLIDE1 20 x 51mm BC, 20 Pk	up to 10,000 spots (150 μ m spot size, 300 μ m pitch)
10486111	FAST SLIDE1 20 x 60mm BC, 20 Pk	up to 10,000 spots (150 μ m spot size, 300 μ m pitch)
10485317	FAST SLIDE2 20X20MM BC, 10Pk	up to 3600 spots per pad (150 μ m spot size, 300 μ m pitch)
10485320	FAST SLIDE8 6X6MM BC, 10Pk	pad spacing: 9 mm (microplate spacing) up to 256 spots (150 μ m spot size, 300 μ m pitch)
10485323	FAST SLIDE16 6x6mm UC, 10Pk	pad spacing: 9 mm (microplate spacing) up to 256 spots (150 μ m spot size, 300 μ m pitch)

PROTEIN ARRAY

5.2 FAST PAK - protein array kit



FAST PAKs (protein array kits) provide the necessary components for researchers to conveniently build and process their own protein microarrays.

FAST PAKs are available in all FAST Slide pad formats (1-, 2-, 8-, and 16-pads). Each kit contains 10 FAST Slides, 10 incubation chambers, 10ml Protein Arraying Buffer (2X), 15ml Protein Array Blocking Buffer, and 125ml Protein Array Washing Buffer (10X). A reusable Slide Holder (sold separately) is also needed, either a FAST Frame for up to four slides or a Chip Clip™ for a single FAST Slide.

Ordering information: FAST Slides

Each kit contains 10 FAST Slides, 10 incubation chambers, 10ml Protein Arraying Buffer (2X), 15ml Protein Array Blocking Buffer, and 125ml Protein Array Washing Buffer (10X).

Product Code	Description
10485262	FAST PAK 1-pad protein array kit 1PK
10485319	FAST PAK 2-pad protein array kit 1PK
10485322	FAST PAK 8-pad protein array kit 1PK
10485325	FAST PAK 16-pad protein array kit 1PK

5.3 FAST BUFFERS- optimized protein array



GVS Life Sciences Protein Array Buffers have been optimized for use on FAST Slides. The buffers include Protein Arraying Buffer, Protein Array Blocking Buffer, and Protein Array Washing Buffer. They are included with several GVS Life Sciences kits, such as the FAST PAKs and FAST Quant, as well as sold separately.

Ordering information: Protein Array Buffers

Product Code	Description
10485331	Protein Arraying Buffer (2X) 10ml 4PK
10485356	Protein Array Blocking Buffer (1X) 100ml 1PK
10485330	Protein Array Washing Buffer (10X) 125ml 4PK

Features & Benefits

- 1, 2, 8, or 16 array pads on each slide
- Suitable for multiplex experiments, side-by-side comparisons, and control experiments all on the same slide
- Multi-well formats allow generation of quantitative arrays
- Reproducible results from slide-to-slide and pad-to-pad
- Increased protein stability and enhanced signal intensity for proteins immobilized with Protein Arraying Buffer
- Proprietary Protein Array Blocking Buffer ensure optimum signal-to-noise ratios and promotes specific binding in immunoassay applications
- Flexibility for detection by fluorescent, chemiluminescent, colorimetric, and isotopic methods

Applications

- ELISA format (sandwich assay) experiments using antibody arrays
- Reverse phase protein arrays using cell or tissue lysates
- Purified protein arrays
- Antigen arrays for antibody screening and autoimmune detection
- Carbohydrate arrays
- Lipids and other materials which can be arrayed on nitrocellulose

PROTEIN ARRAYING BUFFER

Supplied as a 2X concentrate in 10ml plastic bottles.

- Enhances long-term protein stability and molecular recognition activity of arrayed proteins
- Enhances activity of arrayed proteins
- Optimized for use on FAST Slides

PROTEIN ARRAY BLOCKING BUFFER

Supplied neat in 100ml plastic bottles.

- Demonstrates superior blocking of protein microarrays
- Exhibits strong reduction of nonspecific antibody-antibody interactions
- Exhibits minimal effects on specific antibody-antigen interactions
- Results in superior signal-to-noise ratios in protein microarray applications with very low backgrounds
- Compatible with all detection methods

PROTEIN ARRAY WASHING BUFFER

Supplied as a 10X concentrate in 125ml plastic bottles.

- Excellent washing buffer for protein microarrays
- Preserves protein-protein interactions
- Optimized for use on FAST Slides

PROTEIN ARRAY

5.4 FAST Quant™ - human cytokine assays



FAST Quant kits provide high-throughput multiplexed cytokine quantification. Built on FAST technology, FAST Quant combines the high-throughput capabilities of array technology with the quantitative nature of traditional ELISA. With FAST Quant, a researcher can accurately determine the concentration of multiple cytokines in dozens of biological samples simultaneously, using familiar ELISA immunochemistry. Due to the high protein binding capacity of the FAST Slide, FAST Quant immunoassays exhibit greater sensitivity and reproducibility than traditional ELISA

and require significantly smaller sample volumes. Each FAST Quant kit contains 64 arrays, each comprised of 8 to 10 antibodies to human cytokines. Anti-cytokine antibodies are arrayed in triplicate on each array.

Using four 16-pad FAST Slides placed in a FAST Frame (sold separately) FAST Quant offers an 8x12cm footprint - the same as the traditional microplate format. A standard curve can be generated by creating a dilutional series from the recombinant antigen cocktails included in the kit. FAST Quant with the standard curve allows for highly sensitive quantitation of human cytokine concentrations in complex biological samples.

The FAST Quant family of products includes five human cytokine assay kits. Each kit contains four 16-pad FAST Slides pre-arrayed with one of the cytokine antibody panels, four 16-well incubation chambers, recombinant antigen cocktails for generating a standard curve, biotinylated detection antibody cocktails, Protein Array Blocking Buffer, and Protein Array Washing Buffer.

A reusable Slide Holder (sold separately) is also needed: either a FAST Frame for up to four slides or a Chip Clip for a single FAST Slide. Data analysis is seamless using standard data analysis software, or GVS Life Sciences will fluorescently scan the slides and analyze the data for you. Call Customer Care for details.

Ordering information: FAST Quant Kits*

Product Code	Description	Antigens Recognized			
10486031	Human Th1/Th2: antibodies to 9 cytokines associated with the human Th1/Th2 immune response	IL-1 β IL-6 IFN γ	IL-2 IL-10	IL-4 IL-13	IL-5 TNF α
10486060	Human Cytokine II: antibodies to 10 commonly requested human cytokines	IL-1 β IL-8 RANTES	IL-2 IL-10 MCP-1	IL-4 IL-12p70	IL-6 GM-CSF
10486063	Human Angiogenesis: antibodies to 8 cytokines associated with angiogenesis	PDGF-BB MCP-3 MIP-1 α	VEGF MCP-4 MCP-1	FGF β IL-8	Angiogenin Angiopoietin-2
10486064	Human Chemokine: antibodies to 9 commonly requested chemokines	Eotaxin MCP-3 MIP-1 α	RANTES MCP-4	MCP-1 IL-8	MCP-2 IP-10
10486257	Human Cytokine 35: antibodies to 35 of the most requested cytokines	IL-1 α IL-5 IL-12p70 MMP-9 TGF- α ICAM-1 IL-12p40 Fibronectin TNFR-II	IL-1 β IL-6 IL-13 MCP-1 TGF- β 3 IGF-I FGF basic PDGF-BB VEGF	IL-2 IL-8 IFN γ IGF-II Angiogenin IP-10 FAS ligand TGF- β 1 RANTES	IL-4 IL-10 TNF α TNFR-1 Eotaxin GM-CSF MCP-3 TIMP-1

*For Research Use Only

Ordering information: FAST Slide Holders

Required for processing FAST Quant Arrays

Product Code	Description
10486001	FAST Frame Slide Holder for processing up to four FAST Slides INC CHAMBER 2WELL 1/pk
10486081	Chip Clip™ Slide Holder for processing a single FAST Slide 1/pk

PROTEIN ARRAY

5.5 FAST Services - array development and analysis



Scanning and Data Analysis

Using leading-edge automated microarray technologies, GVS Life Sciences provides scanning and data analysis services for FAST Slide users without access to a fluorescent scanner. Whether you are using FAST Quant[®], Serum Biomarker Chip (SBC) kits or have created your own arrays using FAST PAK, GVS Life Sciences will fluorescently scan your FAST Slides and analyze your protein microarray data. You will receive raw fluorescent values and fluorescent values corrected for local background along with some fundamental statistical values such as standard deviation and coefficient of variation. The analysis of FAST Quant results will include generation of antigen specific standard curves and determination of antigen concentrations in experimental samples. SBC analysis will include a comparative ranking of the abundance of all 120 different analytes in your samples. Please call GVS Life Sciences's Customer Care to discuss your specific scanning and data analysis needs.



Microarray Assay Development

Work directly with GVS Life Sciences's protein microarray scientists to create your assay using FAST Slides and either your immunochemicals or immunochemistries provided by GVS Life Sciences. FAST products provide the flexibility, compatibility and reproducible results to meet any assay development requirements. We will work with you to develop and then manufacture a multiplex assay. Project plans will incorporate a gated development approach with go/no go decision points for the most cost effective effort. Typical assay development project results include limits of detection, cross-reactivity and dynamic range for each analyte. GVS Life Sciences's long history and expertise in protein microarray design, development and production make us your ideal partner for all of your microarray assay development needs.

For all FAST Protein Microarray Services, please call GVS Life Sciences's Customer Care.

PROTEIN ARRAY

5.6 FAST Accessories - protein array processing



Increase the ease and convenience of handling and processing FAST Slides and ensure reproducible, consistent results with every slide, every time.

FAST Slide Incubation Chambers

Used in conjunction with the FAST Frame or Chip Clip™ Slide Holder, GVS Life Sciences incubation chambers have a secure gasket design forming a tight, leak-proof seal with the FAST Slides to provide a convenient means to conduct binding reactions on protein microarrays.

Incubation chambers are designed specifically to fit all FAST Slide formats (1-, 2-, 8- and 16-pad). Simply remove the reusable incubation chamber when the reaction is finished.

Features & Benefits

- Superior protein binding capacity
- Highest sensitivity and dynamic range
- Excellent long-term stability of printed proteins
- Compatible with all detection methodologies
- Compatible with commercially available arraying robots

Ordering information: FAST Slide Incubation Chambers

Product Code	Description	Specification
10486137	INC CHMBER SNGLE WELL, 20x51, 10/PK	External dimension: 79 x 25.4 mm Well dimension: 53 x 22 x 4 mm (L x W x D) Volume: 600 to 700 μ l
10486087	INC CHAMBER 2WELL 10/PK	External dimension: 79 x 25.4 mm Well dimension: 21 x 21 x 4 mm (L x W x D) Volume: 300 to 400 μ l
10486046	INC CHMBER 16 WELL 10/PK	External dimension: 79 x 25.4 mm Well dimension: 7 x 7 x 4 mm (L x W x D) Volume: 60 to 100 μ l

Ordering information: FAST Slide Holders

Product Code	Description
10486001	FAST Frame Slide Holder for processing up to four FAST Slides INC CHAMBER 2WELL 10/PK
10486081	Chip Clip™ Slide Holder for processing a single FAST Slide

Ordering information: MicroCaster™ Parts and Accessories

Product Code	Description
10485047	MicroCaster System: 8-pin system hand tool, 8-pin system slide holder, pin conditioner and spare replicator pins, 1/pk
10485061	MicroCaster pin conditioner, 30 ml, 1/pk
10485326	MicroCaster replacement pin, 1/pk
10485370	MicroCaster pad (pin support pad), 1/pk
10486043	Wash and blot station, 1/pk
10486044	96-well microplate indexer, 1/pk

FAST Slide Holders

The Chip Clip securely holds one FAST Slide and incubation chamber for processing multiple arrays simultaneously, ensuring leak-proof barriers around the arrayed pads on the slide. The slide and incubation chamber are easily inserted into and removed from the Chip Clip Slide Holder; side rails hold the chamber firmly against the slide surface. The Chip Clip is compatible with 1 x 3" (25 x 76 mm) slides when used with incubation chambers. The FAST Frame Slide Holder is designed to hold up to four FAST Slides and their corresponding incubation chambers for high-throughput processing of microarrays. The 96-well spacing (9 mm center to center) of the array pads on 16-pad FAST Slides makes the loaded FAST Frame compatible with automated liquid handling systems and 8-channel manual pipettors. Each plate processes up to 64 arrays simultaneously. The rows and columns on each plate are labeled for easy indexing and sample application. The FAST Frame Slide Holder is constructed of autoclavable plastic and is compatible with standard 1 x 3" (25 x 76 mm) glass slides when used with GVS Life Sciences incubation chambers.

MicroCaster™

The MicroCaster is an economical, entry-level manual microarraying system. With the MicroCaster 8-pin hand tool, samples can be loaded from 96-well or 384-well plates.

The MicroCaster Slide Holder accommodates two slides. It has a built-in indexing system that enables precise printing of up to 768 spots in an array of 32 x 24 spots. It is designed for 1-pad FAST Slides with 20 x 51 mm pad size and is compatible with other slide surfaces.

MicroCaster accessories can be used to increase the flexibility of the manual arrayer system by providing accurate source-plate indexing and reliable pin-tool cleaning. The MicroCaster accessories are compatible with standard 96-well microplates and they reduce hassle with pin-tool cleaning.

FILTRATION GUIDE

Membrane information

Filtration through a membrane means that the filter material will stop particles larger than the pore size rating. This enables an absolute pore size rating for the membranes for which they are clearly classified. Bacterial retention claims can be made based on the pore size of the membrane.

Hydrophilic – Hydrophobic Membranes

- **Hydrophilic membranes** have permeability of aqueous solutions and once wetted, they stop gasses. This means that aqueous solutions pass through hydrophilic membranes but gas is stopped when the membrane is wet until the applied pressure exceeds the “bubble point”, at which time the air will evacuate the pore, the liquid is expelled, and the gas will go through. Dry hydrophilic membrane allows gas to pass through. Our HI-FLO PES membranes are hydrophilic membranes.

- **Hydrophobic membranes** have permeability to the gas, but they stop aqueous solutions. In other words, they do the opposite job when compared to hydrophilic membranes. This means that gas will pass through these membranes, but aqueous solutions will be stopped. If air or gas can reach the hydrophobic membrane, it will go through, but if the contact with the hydrophobic membrane is not possible, then the gas will not pass through. The pressure at which aqueous solutions will pass through a hydrophobic membrane is called the water breakthrough (WBT) or water intrusion pressure (WIP). PTFE membranes are hydrophobic membranes. PES membranes are hydrophilic membranes.

Pore size

Pore size is determined by the size of the particle that is expected to be retained with a defined with a high degree of efficiency. Pore size is typically stated in micrometers or microns (μm), and should clearly be designated as either nominal or absolute. Nominal pore size is the ability to retain a majority (60% - 98%) of particles having a specific dimension. Retention efficiency is also depending on such process conditions as concentration, operating pressure etc. Rating parameters can vary among manufacturers. When the pore size, or retention, is “nominal”, it should be stated at a particle size and a percent, i.e., 99.97% retention of 0.3 μm particles. Absolute pore size is the ability to retain the 100% of particles of a specific dimension under defined test conditions (particle size, challenge pressure, concentration, detection method).

Pore Size Challenge Organism

0.1 micron *Acholeplasma laidlawii*


0.2 micron *Brevundimonas diminuta*

0.45 micron *Serratia marcescens*


0.8 micron *Lactobacillus* species

1.2 micron *Candida albicans*

The above table shows proper pore size of hydrophilic membranes to be used to retain the corresponding bacteria. Hydrophobic membranes are about ten times more efficient in retaining bacteria in air than they are in liquids using the same pore size.

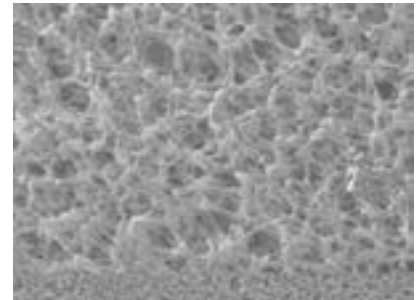

Membrane outside wall

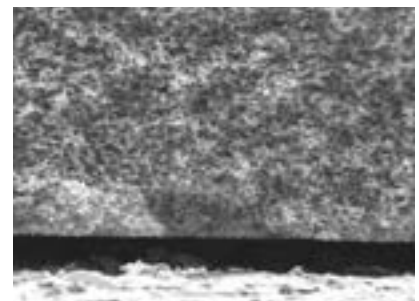

Membrane cross section

Chemical compatibility

This is the ability of the membrane to resist to chemicals without mechanical or chemical damage from chemical exposure. Information about the liquid used with a specific filter material should be outlined before application to determine compatibility, GVS Life Sciences can assist customers in choosing the proper filter (and housing) materials.

Extractables

Extractables are contaminants (typically chemicals) that elute from the filter which might affect quality of the effluent. Wetting agents (surfactants), manufacturing or sterilization residuals are the main cause of undesired extractables. Typical problems caused by extractables are found in the following applications:

- HPLC analysis (strange result)
- Cell culture (cytotoxicity)
- Microbiological analysis (affects the microorganism)
- Environmental analysis (contaminants)

Flushing of the line prior to use can reduce Extractables and their adverse effects.

Binding

This is the property of substances to be filtered having affinity with membranes. This could be a positive effect in some circumstances, but most of the time it can create adverse effects. Particularly it could lead to loss of active components of the liquid to be filtered reducing its beneficial effect. Our PES HI-FLO membrane is low protein binding.

Thermal Stability

This characteristic allows unchanged performance at elevated temperatures. Some membranes can only be sterilized by EtO. Others can be gamma, beta or e-beam sterilized, as well as EtO. Others can be also steam sterilized with no adverse affects. Membrane performance is sometimes reduced at temperature higher than 25°C, and high temperatures can also reduce chemical stability. PTFE membrane is widely stable (any type of sterilization) if the product is designed properly. PES membrane is suggested for EtO and irradiation (no steam sterilization).

Biosafety

These tests are conducted in compliance with ISO-10993 and USP class VI, see specifications Tests that are conducted are: – Cytotoxicity – Sensitization – Irritation intracutaneous reactivity – Systemic toxicity (acute) – Hemocompatibility (Hemolysis)

Pyrogenicity

Pyrogens are chemicals on the filter media and other components that are caused by the waste of dead bacteria. When introduced to a patient, they can elevate the patient's temperature, and can cause complications – even death. Filters that are pyrogenic can make solutions pyrogenic.

They cannot be removed by sterilization, so it is very important that non-pyrogenic filter media and components are used in the production of medical filter devices. The test to determine the pyrogenicity is the LAL test (Limulus Amebocyte Lysate test).

Bubble Point (BP)

Typically this test that is performed on hydrophilic membranes. The BP pressure is the pressure to force air through a wetted hydrophilic membrane. These tests are typically performed with water; however, this test can be conducted on hydrophobic membranes using liquids other than water that will wet the membrane. The BP is an indication of the membrane pore size, as related to actual bacterial retention. This test can also be performed on hydrophobic membranes if the correct solvent (instead of aqueous solution) is used, and is compatible with the entire product.

Water Breakthrough (WBT)

This is the test performed on hydrophobic membranes, and it is also related to the pore size of the membrane. The WBT pressure (sometimes referred to as water intrusion pressure) is the pressure it takes to force an aqueous solution through a hydrophobic membrane.

Water Flow Rate (WFR)

Typically this test is performed on hydrophilic membranes. The BP pressure is the pressure to force air through a wetted hydrophilic membrane, and is an indication of the membrane pore size, as related to actual bacterial retention. This test is



typically performed with water; however, it can be performed with other solutions, as long as the filter media is compatible with the liquid, and it is correlated back to the water bubble point and bacterial retention. The BP test can also be performed on hydrophobic membranes if the correct solvent (instead of aqueous solution) is used, and is compatible with the entire device.

Air Flow (AF)

This is a flow rate typically related to hydrophobic membranes. It is the amount of air that passes through a fixed surface of membrane with a specific applied pressure.

Filter Efficiency (FE)

Quantity of particulate or bacteria retained compared to the total quantity of particulate or bacteria to which the filter is challenged. It is expressed in % and referred to a specific size of particles.

Effective Filtration Area (EFA)

This is the actual filtration area in a device that is subject to filtration. For instance, whereas a 25 mm device may start out with a disc of filter media that is cut to 25 mm, the sealing surfaces should be eliminated from the calculations of the device EFA.

Main membranes used are:

Cellulose Acetate (CA): Hydrophilic membrane. Low protein binding, (reducing filter changes when filtering proteinaceous solutions?). Ideal for protein, cell culture media and enzymes filtrations, tissue culture media sterilization, biological fluid filtration and other filtration applications where maximum recovery of proteins is critical.

Nylon (NY): Hydrophilic membrane. Ideal for use in general filtration or medical assays. Superior strength, resistant to a range of organic solvents. Low extractables. High protein binding capacity. Lot-to-lot consistency.

Polyethersulfone (PES): Hydrophilic membrane. Designed to remove particulates during general filtration, low protein and drug binding characteristics make it ideally suited for use in life science applications. Its strength and durability are advantageous during usage that involves aggressive handling or automated equipment. Low protein and drug binding characteristics maximize recovery of critical drugs used in I.V. therapy, chemotherapy and open-heart surgery.

Nitrocellulose Mixed Esters (MCE): Hydrophilic membrane. Aqueous clarification and particulate capture. Consistent high flow rate for faster filtration. Uniform pore structure for selectivity. Hydrophilic, inert cellulose nitrate. High binding capacity. Manufactured thickness within 10 microns

Polyvinylidene Difluoride (PVDF): Hydrophilic membrane. Ideal for use in Sterilizing and Clarifying filtration of biological solutions. High Flow Rates. Low Extractables. Broad Chemical Compatibility. Very low protein binding.

Polytetrafluoroethylene (PTFE): Hydrophobic membrane. Ideal for filtration of strong acids and aggressive solutions, venting applications, phase separations, aerosol samplings. Chemically and biologically inert. Superior chemical resistance. Can withstand high temperatures.

Glass Fiber (GF): Hydrophilic material. Used also as a pre-filter to extend membrane life. Eliminate sample contamination. Excellent wet strength for each handling and filter integrity. Ideal for water/air pollution analysis, liquid clarification and cell harvesting.

Regenerated Cellulose (RC): Hydrophilic membrane. Resistant to a very wide range of solvents. Suitable for use with either aqueous solutions or organic solvents. Compatible with HPLC solvents. Very low protein binding capacity and hence excellent for protein recovery applications.



Polymers information

Thermoplastics and thermosets are the two basic groups of plastic materials. Thermoplastic resins can be repeatedly melted and solidified by heating and cooling so that any scrap generated in processing can be theoretically reused. No chemical change generally takes place during forming. Usually, thermoplastic polymers are supplied in the form of pellets, which often contain additives to enhance processing or to provide necessary characteristics in the finished product (e.g., color, conductivity, etc.). The temperature service range of thermoplastics is limited by their loss of physical strength and eventual melting at elevated temperatures.

Main Thermoplastic resin used are:

Polypropylene (PP)

It is similar to polyethylene, but each unit of the chain has a methyl group attached. It is translucent, autocavable, and has no known solvent at room temperature. It is slightly more susceptible to strong oxidizing agents than conventional polyethylene because of its many branches (methyl groups, in this case). Polypropylene is noted for its excellent chemical resistance in corrosive environments. This polymer is easily welded and machined.

TYPICAL PROPERTIES OF PP ARE:

- Clean/High Purity
- Good Dimensional Stability
- Good Organoleptic Properties
- High Clarity
- High Flow
- High Stiffness
- Homopolymer
- Low Warpage
- Narrow Molecular Weight Distribution
- Nucleated

Polyethylene (PE) Plastic

Huge family of resins obtained by polymerizing ethylene gas, and it is available in a range of flexibilities. Polyethylene can be formed by a wide variety of thermoplastic processing methods and is particularly useful where moisture resistance is required. Low-density polyethylene (LDPE) has more extensive branching, resulting in a less compact molecular structure. High-density polyethylene (HDPE) has minimal branching, which makes it more rigid and less permeable than LDPE. Linear low-density polyethylene (LLDPE) combines the toughness of low-density polyethylene with the rigidity of high-density polyethylene.

TYPICAL PROPERTIES OF PE ARE:

- Good Processability
- Food Contact Acceptable
- Antioxidant
- High ESCR (Stress Crack Resist.)
- Low Density
- High Impact Resistance

Acrylic-based polymer

Acrylic polymer developed especially for use in the Medical Device Industry. The material is transparent and tough offer gamma and ETO sterilization resistance, and they are easy to process and weld easily to PVC. Typical applications include disposable medical diagnostic devices such as cassettes and cuvettes.

TYPICAL PROPERTIES OF acrylic-based polymer compounds are:

- Excellent chemical resistance to fats and oils
- Excellent bonding and welding capabilities
- Excellent bonding to PVC tubing
- Good impact strength



- Good light transmission
- Good resistance to EtO, gamma and E-beam sterilization
- Superior resistance to lipids and alcohol
- Excellent ductility

Nylon (PA6)

This is a group of linear polymers with repeated amide linkages along the backbone. These are produced by an amidation of diamines with dibasic acids, or polymerisation of amino acids. Nylon is strong and tough. It resists abrasion, fatigue and impact. Nylon offers excellent chemical resistance with negligible permeation rates when used with organic solvents. However, it has poor resistance to strong mineral acids, oxidizing agents and certain salts.

TYPICAL PROPERTIES OF PA 6 ARE:

- Good Chemical Resistance
- Good Colorability
- Good Corrosion Resistance
- Good Processability
- Good Toughness
- Good Wear Resistance
- High Rigidity
- High Strength
- Low Friction



COMPATIBILITY CHART

		Filter Media											Housing			
		Cellulose Acetate	Glass Fiber (binder)	Glass Fiber (no binder)	Polyethersulfone	Nitrocellulose	Nylon	Polycarbonate	Polyester	Polypropylene	PTFE (unlaminated)	PTFE (laminated)	Silver	Modified Acrylic	Polysulfone	Polystyrene
Chemical		ca	gfb	gfn	pes	nc	ny	pc	pet	pp	ptu	ptl	ag	ac	ps	pst
ACIDS	Acetic Acid 5%	R	T	R	R	R	R	R	R	R	R	R	R	N	R	R
	Acetic Acid 10%	N	T	R	R	N	L	R	R	R	R	R	R	N	R	R
	Acetic Acid, Glacial	N	N	R	R	N	N	L	R	R	R	R	R	N	R	R
	Boric Acid	R	T	T	T	R	L	R	R	R	R	R	R	N	R	R
	Hydrochloric, 6N	L	N	R	R	N	N	R	L	R	R	R	R	N	R	R
	Hydrochloric, Conc.	N	N	R	R	N	N	R	N	R	R	R	R	N	R	R
	Hydrofluoric, 10%	N	N	N	T	N	N	R	R	R	R	R	R	T	T	T
	Hydrofluoric, 35%	N	N	N	T	N	N	R	R	T	R	T	R	T	T	T
	Nitric Acid, 6N	L	N	L	N	R	N	R	R	L	R	L	N	N	N	L
	Nitric Acid, Conc.	N	N	L	N	N	N	R	N	N	R	N	N	N	N	N
	Sulfuric Acid, 6N	L	N	R	T	R	N	R	R	L	R	L	N	N	N	N
	Sulfuric Acid, Conc.	N	N	R	N	N	N	N	N	N	R	N	N	N	N	N
ALCOHOLS	Amly Alcohol	R	R	R	N	N	R	R	R	R	R	R	R	N	R	N
	Benzyl Alcohol	L	N	N	N	R	L	L	R	R	R	R	R	R	R	N
	Butyl Alcohol	R	R	R	R	R	R	R	R	R	R	R	R	R	R	T
	Butyl Cellosolve	L	R	R	T	N	R	L	R	R	R	R	R	T	L	T
	Ethyl Alcohol < 80%	R	R	R	R	R	R	R	R	R	R	R	R	L	R	L
	Ethyl Alcohol > 80%	R	R	R	R	L	R	R	R	R	R	R	R	L	R	N
	Ethylene Glycol	R	R	R	R	L	R	R	R	R	R	R	R	T	R	T
	Glycerine (Glycerol)	R	R	R	R	R	R	R	R	R	R	R	R	T	R	T
	Isobutyl alcohol	R	N	N	T	R	R	R	R	R	R	R	R	R	R	R
	Isopropanol	R	R	R	R	L	R	R	R	R	R	R	R	T	R	T
	Methanol	R	R	R	R	N	L	R	R	R	R	R	R	R	R	R
	Methyl Cellosolve	L	R	R	T	L	R	N	R	R	R	R	R	T	R	T
	Propanol	R	R	R	T	R	R	R	R	R	R	R	R	T	R	T
BASES	Ammonium Hydroxide, 6N	N	N	R	R	N	N	N	L	R	R	R	R	R	R	R
	Potassium Hydroxide, 6N	N	N	T	T	N	R	N	N	R	R	R	R	T	R	T
	Sodium Hydroxide, 6N	N	N	T	R	N	N	N	L	R	R	R	R	T	T	T
SOLVENTS	Acetone	N	R	R	N	N	R	L	R	R	R	R	R	N	N	N
	Acetonitrile	N	T	R	R	N	R	L	R	R	R	R	T	N	N	N
	Amyl Acetate	L	N	R	L	N	R	R	R	R	R	R	R	N	N	N
	aniline	N	T	T	R	N	R	N	R	R	R	R	R	T	N	T
	Benezene	L	N	R	R	R	R	L	R	L	R	L	R	N	N	N
	Bromoform	N	R	R	T	R	R	N	R	R	R	R	R	T	N	T
	Butyl Acetate	L	N	R	L	N	R	R	R	R	R	R	R	N	N	N
	Carbon Tetrachloride	L	N	N	R	R	R	L	R	L	R	L	R	N	N	N
	Cellosolve	R	R	R	T	N	R	R	R	R	R	R	R	N	N	T

	Chloroform	N	R	R	N	R	R	N	R	L	R	L	R	N	L	N	
	Cyclohexane	R	R	R	T	R	R	R	R	R	R	R	R	N	R	T	
	Cyclohexanone	N	R	R	N	N	R	L	R	R	R	R	R	N	N	N	
	Diethyl Acetamide	N	R	R	T	N	R	L	R	N	R	N	R	N	N	N	
	Dimethyl Formamide	N	N	R	N	N	R	N	R	R	R	R	R	N	N	N	
	Dimethyl Sulfoxide (DMSO)	N	N	R	N	N	R	N	R	R	R	R	T	N	N	N	
	Dioxane	N	R	R	L	N	R	N	R	R	R	R	R	N	N	N	
	Ethyl Ether	L	T	R	R	L	R	R	R	R	R	R	R	N	L	N	
	Ethylene Dichloride	L	R	R	T	L	R	N	R	R	R	R	R	T	N	T	
	Formaldehyde	L	R	R	R	N	R	R	R	R	R	R	R	N	R	N	
	Freon TF	R	R	R	R	R	R	R	R	R	R	R	R	L	R	N	
	Gasoline	R	R	R	T	R	R	R	R	R	R	R	R	N	R	N	
	Hexane	R	L	R	T	R	R	R	R	R	R	R	R	N	R	N	
	Isopropyl Acetate	N	N	R	T	N	R	R	R	R	R	R	R	N	N	N	
	Kerosene	R	R	R	T	R	R	R	R	R	R	R	R	N	N	N	
	Methyl Acetate	N	N	R	T	N	R	N	R	R	R	R	R	N	N	N	
	Methyl Ethyl Ketone (MEK)	N	R	R	N	N	R	L	R	R	R	R	R	N	N	N	
	Methyl Isobutyl Ketone	N	R	R	T	N	R	L	T	R	R	R	R	N	N	N	
	Methylene Chloride	N	R	R	N	N	L	N	R	R	R	R	R	N	N	N	
	Nitrobenzene	N	N	N	N	N	R	N	R	R	R	R	T	N	N	N	
	Pentane	R	R	R	R	R	R	R	R	L	R	L	R	N	R	N	
	Perchloroethylene	R	N	N	N	R	R	R	T	R	R	R	R	N	L	N	
	Pyridine	N	N	R	N	N	R	N	R	R	R	R	R	N	N	N	
	Tetrahydrofuran	N	T	L	N	N	L	N	R	L	L	L	R	N	N	N	
	Toluene	L	N	R	N	R	R	L	R	L	R	L	R	N	N	N	
	Trichloroethane	L	T	T	R	N	R	N	T	R	R	R	R	N	N	N	
	Trichlorethylene	R	N	N	R	R	R	B	R	L	L	L	R	N	N	N	
	Triethylamine	R	R	R	T	L	R	L	R	R	R	R	R	T	N	T	
	Xylene	R	R	R	L	R	R	R	R	L	R	L	R	N	N	N	
MISCELLANEOUS	Cottonseed Oil	R	L	R	T	R	R	R	T	R	R	R	R	T	R	T	
	Hydrogen Peroxide (30%)	R	R	R	T	R	R	R	R	R	R	R	R	T	R	T	
	Kodak KMER FTFR	N	N	N	T	N	R	R	R	R	R	R	R	N	R	N	
	Peanut Oil	R	R	R	T	R	R	R	R	R	R	R	R	T	R	T	
	Petroleum Oils	T	T	T	L	R	T	R	R	T	R	T	R	T	T	T	
	Sesame Oil	R	R	R	T	R	R	R	R	R	R	R	R	R	T	R	T
	Shibley (AS-111,340,1350)	N	N	N	T	N	R	R	R	R	R	R	R	R	N	R	N
	Silicone Oils	R	R	R	R	R	R	R	R	R	R	R	R	R	T	R	T
	Turpentine	R	R	R	T	R	R	R	R	R	R	R	R	R	T	R	T
	Waycoat 59	N	N	N	T	N	R	R	R	R	R	R	R	R	N	R	N

PRODUCT CODE INDEX

1145328	53	1212701	14	1213155	18	1213700	8	1213791	42	1214227	18	1214726	66	1215069	49
1145331	53	1212721	65	1213157	19	1213701	14	1213792	42	1214232	28	1214756	43	1215070	49
1145332	53	1212763	60	1213158	18	1213702	14	1213793	42	1214237	44	1214762	65	1215071	48
1145333	53	1212781	66	1213160	18	1213705	8	1213794	42	1214238	44	1214778	15	1215073	49
1145334	53	1212783	66	1213161	19	1213706	8	1213795	14	1214239	44	1214802	66	1215074	38
1145335	53	1212790	48	1213190	65	1213708	14	1213796	42	1214240	44	1214842	66	1215077	49
1145336	53	1212846	38	1213192	15	1213709	14	1213797	42	1214241	44	1214851	38	1215078	49
1145337	53	1212899	20	1213195	65	1213710	8	1213798	42	1214243	64	1214861	39	1215116	48
1145341	53	1212905	20	1213305	38	1213711	14	1213799	42	1214245	62	1214866	39	1215117	48
1145342	53	1212908	20	1213314	64	1213721	8	1213800	42	1214247	65	1214872	39	1215118	48
1145343	53	1212910	20	1213325	60	1213722	8	1213801	42	1214248	65	1214880	42	1215130	14
1145345	53	1212911	20	1213331	39	1213723	14	1213802	42	1214250	23	1214881	42	1215138	48
1145346	53	1212929	19	1213343	38	1213724	14	1213803	42	1214260	23	1214882	40	1215141	52
1145347	53	1212930	19	1213358	38	1213725	14	1213804	38	1214278	28	1214898	40	1215142	48
1145348	53	1212932	19	1213370	62	1213726	14	1213805	38	1214282	28	1214904	14	1215145	50
1212374	38	1212933	19	1213372	62	1213730	38	1213806	14	1214286	28	1214909	40	1215146	14
1212375	38	1212935	19	1213373	62	1213731	14	1213807	42	1214288	28	1214911	64	1215147	49
1212379	44	1212936	18	1213375	62	1213737	14	1213808	42	1214293	28	1214912	60	1215148	49
1212380	44	1212937	18	1213379	62	1213738	14	1213809	42	1214301	24	1214920	43	1215150	50
1212381	44	1212939	20	1213380	62	1213739	14	1213810	42	1214319	28	1214923	39	1215151	50
1212383	44	1212942	38	1213384	62	1213740	14	1213811	42	1214344	14	1214932	15	1215152	50
1212390	44	1212949	19	1213391	62	1213742	8	1213812	42	1214346	14	1214935	65	1215153	50
1212391	44	1212950	19	1213409	62	1213743	8	1213813	42	1214349	14	1214941	40	1215154	18
1212392	44	1212951	19	1213410	62	1213744	14	1213814	14	1214357	38	1214946	8	1215161	48
1212394	44	1212953	19	1213411	62	1213745	14	1213815	42	1214358	14	1214950	40	1215162	60
1212517	38	1212954	19	1213412	62	1213746	14	1213816	42	1214359	14	1214956	42	1215163	52
1212590	65	1212971	19	1213416	62	1213751	28	1213817	42	1214360	14	1214959	40	1215164	8
1212591	65	1212972	19	1213419	62	1213753	28	1213818	42	1214370	38	1214966	15	1215170	64
1212592	65	1212975	19	1213422	62	1213757	20	1213819	42	1214391	66	1214974	60	1215173	52
1212594	65	1212978	18	1213516	38	1213758	14	1213820	42	1214396	39	1214978	65	1215179	19
1212596	65	1212982	18	1213529	20	1213760	42	1213821	42	1214411	38	1214991	39	1215183	52
1212597	65	1212987	19	1213540	20	1213761	42	1213822	14	1214423	66	1215019	40	1215185	48
1212598	65	1212992	19	1213550	20	1213762	42	1213823	42	1214428	62	1215027	40	1215188	48
1212599	65	1213037	38	1213561	20	1213763	42	1213825	42	1214429	66	1215032	66	1215190	14
1212600	65	1213057	19	1213577	20	1213764	42	1213828	14	1214436	18	1215036	49	1215194	50
1212601	65	1213058	19	1213578	19	1213765	42	1213845	42	1214443	45	1215037	66	1215198	48
1212602	65	1213070	19	1213584	19	1213766	42	1213865	24	1214456	40	1215041	48	1215201	50
1212604	65	1213078	19	1213591	19	1213767	14	1213882	24	1214457	20	1215042	49	1215215	50
1212620	38	1213080	19	1213596	19	1213768	42	1213883	24	1214465	43	1215043	49	1215218	50
1212630	14	1213081	19	1213605	19	1213769	42	1213888	64	1214475	43	1215044	49	1215219	50
1212631	40	1213089	19	1213608	18	1213770	42	1213889	48	1214509	62	1215046	49	1215230	39
1212632	65	1213090	19	1213610	18	1213771	42	1213943	65	1214526	22	1215047	49	1215236	48
1212636	66	1213091	19	1213611	19	1213772	42	1213956	18	1214527	40	1215048	49	1215237	39
1212637	66	1213095	19	1213616	19	1213774	42	1213958	38	1214532	43	1215049	49	1215245	14
1212638	66	1213098	19	1213617	19	1213775	42	1213991	64	1214533	40	1215050	49	1215249	39
1212639	66	1213100	19	1213619	19	1213776	42	1213992	66	1214554	40	1215051	49	1215257	40
1212640	65	1213101	19	1213620	19	1213778	42	1213999	64	1214558	66	1215052	49	1215262	28
1212642	66	1213111	19	1213621	19	1213779	42	1214001	18	1214560	65	1215053	49	1215263	40
1212643	66	1213113	19	1213622	19	1213780	42	1214014	15	1214565	40	1215054	49	1215268	14
1212644	66	1213117	19	1213629	8	1213782	42	1214038	38	1214567	48	1215055	49	1215271	48
1212648	38	1213121	64	1213641	15	1213783	42	1214083	18	1214568	43	1215056	49	1215272	40
1212661	49	1213122	14	1213643	39	1213784	42	1214124	24	1214588	66	1215057	49	1215274	48
1212668	28	1213124	38	1213648	8	1213785	14	1214126	65	1214600	66	1215058	49	1215275	48
1212669	65	1213125	38	1213657	14	1213786	42	1214144	8	1214604	43	1215059	49	1215281	40
1212689	65	1213141	65	1213667	14	1213788	42	1214193	43	1214661	28	1215062	49	1215288	52
1212690	65	1213145	39	1213678	14	1213789	42	1214220	14	1214669	43	1215063	49	1215290	48
1212692	8	1213154	18	1213699	14	1213790	42	1214225	18	1214671	60	1215068	49	1215294	14

1215295	14	1215467	64	1215570	60	1220702	52	1221414	52	1222096	50	1224113	15	1224760	11
1215297	48	1215469	64	1215571	60	1220823	48	1221415	52	1222102	44	1224114	15	1224761	11
1215303	50	1215470	64	1215572	60	1220824	44	1221416	52	1222103	44	1224117	15	1224763	11
1215304	50	1215471	64	1215573	60	1220827	52	1221417	52	1222216	66	1224119	15	1224764	11
1215305	40	1215472	64	1215576	60	1220835	49	1221418	52	1222218	66	1224120	15	1224765	11
1215308	52	1215473	64	1215577	60	1220860	50	1221419	52	1222230	43	1224121	15	1224766	11
1215309	50	1215474	64	1215578	60	1220861	50	1221420	52	1222240	52	1224124	15	1224767	11
1215310	62	1215475	64	1215579	60	1220867	48	1221423	52	1222242	52	1224125	15	1224768	11
1215311	48	1215476	64	1215580	60	1220868	49	1221424	52	1222246	52	1224126	15	1224769	11
1215315	48	1215477	64	1215582	60	1220871	52	1221426	52	1222248	52	1224127	15	1224773	11
1215316	40	1215478	64	1215583	60	1220886	52	1221429	48	1222249	52	1224128	15	1224774	11
1215317	50	1215480	64	1215584	60	1220887	60	1221441	42	1222251	52	1224129	15	1224775	11
1215319	65	1215481	64	1215588	60	1220891	50	1221485	50	1222253	52	1224133	15	1224776	11
1215321	49	1215482	64	1215589	60	1220941	49	1221496	14	1222254	52	1224135	15	1224780	11
1215323	40	1215483	64	1215605	49	1220950	22	1221503	48	1222256	52	1224136	15	1224781	11
1215324	52	1215484	64	1215606	49	1220969	52	1221504	49	1222258	52	1224137	15	1224782	11
1215367	52	1215485	45	1215608	49	1220970	50	1221513	14	1222260	52	1224138	15	1224783	11
1215368	43	1215486	45	1215609	48	1220987	50	1221546	38	1222266	14	1224143	15	1224787	11
1215373	52	1215487	45	1215610	49	1220988	50	1221730	38	1222267	43	1224144	15	1224788	11
1215374	52	1215488	45	1215611	49	1221004	50	1221746	49	1222287	65	1224145	15	1224789	11
1215376	40	1215489	45	1215612	49	1221005	50	1221768	20	1222295	65	1224146	15	1224790	11
1215379	52	1215490	45	1215613	49	1221008	43	1221871	49	1222323	18	1224150	15	1224794	11
1215380	40	1215491	45	1215614	49	1221009	49	1221948	39	1222327	18	1224151	15	1224795	11
1215385	50	1215492	45	1215615	49	1221181	48	1221955	40	1222369	14	1224153	15	1224796	11
1215388	50	1215493	45	1215617	49	1221182	52	1221974	65	1222370	14	1224157	15	1224797	11
1215392	64	1215494	45	1215618	49	1221227	50	1221976	64	1222374	14	1224158	15	1224801	11
1215394	64	1215495	45	1215619	49	1221229	49	1221981	64	1222375	14	1224159	15	1224802	11
1215396	43	1215496	45	1215620	49	1221230	48	1221983	64	1222378	14	1224160	15	1224803	11
1215403	50	1215500	52	1215621	49	1221231	48	1221996	60	1222388	14	1224164	15	1224804	11
1215406	39	1215502	45	1215622	49	1221232	48	1222017	49	1222395	14	1224165	15	1224808	11
1215407	39	1215503	45	1215623	49	1221286	48	1222018	49	1222432	18	1224166	15	1224809	11
1215408	39	1215504	45	1215624	49	1221290	50	1222025	48	1222452	48	1224167	15	1224810	11
1215409	39	1215505	45	1215625	49	1221291	50	1222026	50	1222479	50	1224172	15	1224811	11
1215415	40	1215506	45	1215627	49	1221292	50	1222027	48	1222482	50	1224173	15	1224812	11
1215416	40	1215507	45	1215628	49	1221293	49	1222028	48	1222851	42	1224174	15	1224813	11
1215424	40	1215522	52	1215629	49	1221295	48	1222030	48	1222852	42	1224175	15	1224814	11
1215425	40	1215533	38	1215630	49	1221296	49	1222032	49	1222853	42	1224176	15	1224815	11
1215426	40	1215534	52	1215631	49	1221297	49	1222035	48	1222854	42	1224177	15	1224817	11
1215427	38	1215540	60	1215632	49	1221300	49	1222037	49	1222855	42	1224211	38	1224818	11
1215428	40	1215541	60	1215633	49	1221302	49	1222038	49	1222856	42	1224304	14	1224966	44
1215431	40	1215542	60	1215634	49	1221329	50	1222063	49	1222857	42	1224305	14	1225120	49
1215432	40	1215543	60	1215635	38	1221331	49	1222064	49	1222858	42	1224306	14	1225184	48
1215433	40	1215544	60	1215637	49	1221334	52	1222070	50	1222859	42	1224307	14	1225346	18
1215438	40	1215547	60	1215638	49	1221383	52	1222071	50	1222980	39	1224310	15	1225460	39
1215440	40	1215548	60	1215676	38	1221385	52	1222072	49	1223036	49	1224311	15	1225487	44
1215441	40	1215549	60	1215730	52	1221387	52	1222073	49	1223129	18	1224312	15	1225489	44
1215442	40	1215550	60	1215984	49	1221388	52	1222074	50	1223273	64	1224313	15	1225490	44
1215443	40	1215551	60	1215985	49	1221389	52	1222075	50	1223340	43	1224314	15	1225509	60
1215448	40	1215553	60	1216611	50	1221397	52	1222077	49	1223379	14	1224315	15	1225602	11
1215450	40	1215554	60	1216718	39	1221398	52	1222080	48	1223845	18	1224487	43	1225607	11
1215451	40	1215555	60	1216719	39	1221399	52	1222081	49	1223871	43	1224492	43	1225617	11
1215452	40	1215557	60	1216720	39	1221401	52	1222082	49	1224003	43	1224496	43	1225618	11
1215453	40	1215559	60	1216721	39	1221402	52	1222084	50	1224067	4	1224680	50	1225619	11
1215455	40	1215561	60	1216722	39	1221403	52	1222085	48	1224085	14	1224690	32	1225620	11
1215456	40	1215562	60	1216723	39	1221404	52	1222086	49	1224100	15	1224746	11	1225622	11
1215458	64	1215563	60	1216724	39	1221405	52	1222088	50	1224101	15	1224747	11	1225623	11
1215459	64	1215564	60	1220668	49	1221406	52	1222089	49	1224103	15	1224748	11	1225781	40
1215460	64	1215565	60	1220671	42	1221409	52	1222091	50	1224104	15	1224749	11	1225783	49
1215463	64	1215566	60	1220678	60	1221410	52	1222092	49	1224105	15	1224753	11	1225792	44
1215464	64	1215567	60	1220684	19	1221411	52	1222093	49	1224106	15	1224754	11	1225881	43
1215465	64	1215568	60	1220686	48	1221412	52	1222094	50	1224110	15	1224755	11	1225882	43
1215466	64	1215569	60	1220694	49	1221413	52	1222095	50	1224112	15	1224756	11	1225883	43

1225884	43	1229458	15	1262579	23	3034574	60	10485317	70	FJ25BNPPE002AD01	12	FP150DSL440ANC01	58
1225894	50	1229459	15	1262953	14	3034848	50	10485319	71	FJ25BNPPE005AD01	12	FP150DXF040ALC01	54
1225932	44	1229460	11	1264834	50	3034974	38	10485320	70	FJ25BNPPH002AD01	12	FP150DXS050ALC01	56
1225982	42	1229461	15	1264835	48	3034975	38	10485322	71	FJ25BNPPS002AD01	12	FP150DXS420ANC01	58
1225985	43	1229462	11	1266844	60	3037376	43	10485323	70	FJ25BNPPS004AD01	12	ME025NPH002BC01	46
1226150	15	1229463	15	1267014	50	3037800	46	10485325	71	FJ25BNPPT002AD01	12	ME025NRC002BC01	46
1226556	63	1229464	11	1267667	48	3037801	46	10485326	74	FJ25BNPPT004AD01	12	ME025NRC004BC01	46
1226557	63	1229465	15	1267681	45	3038551	15	10485330	71	FJ25BNPPV002AD01	12	ME047NPH002BC01	46
1226558	63	1229466	15	1268126	48	3038552	15	10485331	71	FJ25BNPPV004AD01	12	ME047NRC002BC01	46
1226559	63	1229467	15	1268603	60	3038824	50	10485356	71	FJ25BNPPX002AD01	12	ME047NRC004BC01	46
1226560	63	1229469	11	1269206	44	3041202	38	10485370	74	FJ25BNPRC002AD01	12	MV32ANPNY002BC01	25
1226561	63	1229470	15	1270152	15	3044268	20	10486001	72	FJ25BNPRC004AD01	12	MV32ANPNY002BM01	25
1226562	63	1229471	11	1270153	15	3044270	46	10486031	72	FJ25BSCPS002AL01	13	MV32ANPNY004UC01	25
1226563	63	1229472	15	3002536	48	3044271	46	10486043	74	FJ25BSCPS004AL01	13	MV32ANPNY004UM01	25
1226564	63	1229473	11	3003072	60	3048260	42	10486044	74	FJ25BSCP002AL01	13	MV32ANPPS002EC01	25
1226565	63	1229474	15	3007164	52	3048300	48	10486046	74	FJ25BSCP004AL01	13	MV32ANPPS002EM01	25
1226566	63	1229475	11	3008749	52	3048982	48	10486060	72	FP047DFAFAGLFC01	59	MV32ANPPS004WC01	25
1226567	63	1229476	15	3012605	14	3049247	38	10486063	72	FP047DMFCGLFC01	59	MV32ANPPS004WM01	25
1226568	63	1229477	11	3013362	45	3049950	11	10486064	72	FP055DFAFAGLFC01	59	MV32ANPPT002TC01	25
1226569	63	1229478	15	3013824	50	3049952	11	10486081	72	FP055DMFCGLFC01	59	MV32ANPPT002TM01	25
1226570	63	1229479	11	3013825	50	3049953	15	10486087	74	FP090DFA410ANC01	56	MV32ANPPT004CC01	25
1226571	63	1229480	15	3013826	42	3050049	15	10486111	70	FP090DFAFAGLFC01	59	MV32ANPPT004CM01	25
1226572	63	1229481	11	3014223	52	3050121	15	10486137	74	FP090DME010ALC01	54	MV32ANPPV002FC01	25
1226573	63	1229540	48	3014503	52	3050122	15	10486257	72	FP090DME430ANC01	57	MV32ANPPV002FM01	25
1226574	63	1229555	28	3015810	60	3050851	40	10496101	32	FP090DMFCGLFC01	59	MV32ANPPV004IC01	25
1226575	63	1230356	42	3018505	43	3052148	42	10496102	32	FP090DMS020ALC01	55	MV32ANPPV004IM01	25
1226664	43	1232919	49	3019423	11	3052874	38	10496103	32	FP090DMS030LTC01	55	MV32ANPRC002GC01	25
1226665	43	1232921	43	3019424	15	3053082	40	10496104	32	FP090DMS400ANC01	57	MV32ANPRC002GM01	25
1226916	15	1233373	49	3019515	48	3054144	48	10496106	32	FP090DSL440ANC01	58	MV32ANPRC004LC01	25
1226917	15	1233541	15	3020351	15	3057106	43	10496108	32	FP090DXF040ALC01	54	MV32ANPRC004LM01	25
1226939	15	1233543	15	3020528	15	3061196	38	10496109	32	FP090DXS050ALC01	56	VF40CBPFG000AL01	16
1226940	15	1233544	11	3022140	48	3095433	44	10496112	32	FP090DXS420ANC01	58	VF40CSPGF000AX01	16
1226941	15	1233545	11	3023084	15	10443401	21	10496113	32	FP110DFA410ANC01	56	VF50ANPPT002AC01	16
1226942	15	1233547	11	3023135	11	10443411	21	10496114	32	FP110DFAFAGLFC01	59	VF50ANPPT004AC01	16
1226943	15	1233548	11	3023187	11	10443421	21	10496116	32	FP110DME010ALC01	54	VF50ASPPT002AX01	16
1226944	15	1233549	15	3023188	11	10443423	21	10496119	32	FP110DME430ANC01	57	VF50ASPPT004AX01	16
1226945	15	1233550	15	3023783	50	10443425	21	10496120	32	FP110DMFCGLFC01	59		
1226946	15	1233551	15	3023784	50	10443430	21	10496121	32	FP110DMS020ALC01	55		
1226947	15	1233863	43	3023785	50	10443435	21	10496124	32	FP110DMS030LTC01	50		
1226950	15	1233871	11	3026028	45	10448030	21	10496125	32	FP110DMS400ANC01	57		
1226951	15	1233882	11	3026365	43	10448031	21	10496146	32	FP110DSL440ANC01	58		
1227203	49	1234298	49	3026431	48	10467001	20	10496151	32	FP110DXF040ALC01	54		
1227204	15	1235258	8	3026845	44	10467002	20	10496164	32	FP110DXS050ALC01	56		
1227205	15	1235299	45	3027176	48	10467003	20	10496187	32	FP110DXS420ANC01	58		
1227213	48	1235494	48	3027177	48	10467004	20	10496192	32	FP125DFA410ANC01	56		
1227264	48	1235748	43	3027598	49	10467005	20	10496700	32	FP125DFAFAGLFC01	59		
1227275	15	1236292	43	3029505	43	10467006	20	10496705	32	FP125DME010ALC01	54		
1227352	49	1236363	48	3029939	60	10467007	20	10496706	32	FP125DME430ANC01	57		
1228132	49	1236904	42	3030900	43	10467008	20	10496707	32	FP125DMFCGLFC01	59		
1228342	50	1237235	15	3031100	40	10467009	20			FP125DMS020ALC01	55		
1228824	42	1237423	45	3031900	20	10467010	20			FP125DMS030LTC01	55		
1229443	15	1237721	15	3032159	48	10467011	20			FP125DMS400ANC01	57		
1229444	15	1238010	52	3032862	50	10467012	20			FP125DSL440ANC01	58		
1229445	15	1239192	50	3032876	43	10467013	20			FP125DXF040ALC01	54		
1229447	11	1239369	14	3033092	48	10467015	20			FP125DXS050ALC01	56		
1229448	15	1239556	48	3033093	48	10467017	20			FP125DXS420ANC01	58		
1229449	11	1239557	48	3033301	48	10467019	20			FP150DFA410ANC01	56		
1229450	15	1239558	48	3033302	48	10467021	20			FP150DME010ALC01	54		
1229451	15	1239715	14	3034261	48	10484182	70			FP150DME430ANC01	57		
1229452	11	1240382	38	3034300	45	10485047	74			FP150DMS020ALC01	55		
1229453	15	1241477	42	3034567	48	10485061	74			FP150DMS030LTC01	55		
1229454	11	1242278	43	3034573	60	10485262	71			FP150DMS400ANC01	57		

CATALOG GVS Life Sciences 2014
Copyright © 2014 GVS ® S.p.A
All Right Reserved
Printed in Italy

Printing History:
Version: 010314

While every precaution has been taken in the preparation of this catalog, data are subject to change without notice.

Results in specific application of GVS Life Sciences products may vary according to the conditions and applications. GVS Life Sciences assumes no responsibility for damage resulting from incorrect use of our products.





WORLDWIDE

EUROPE

Italy Office
Headquarters
GVS S.p.A.
Via Roma 50
40069 Zola Predosa (BO) - Italy
tel. +39 051 6176311
fax +39 051 6176200
gvs@gvs.com

Spain
Spain Office (Laboratory)
Avila 50, Edificio Miete, 3-6
08005 Barcelona - Spain
tel. +34 93 300 90 57
gvsspain@gvs.com

United Kingdom
GVS Filter Technology UK Ltd.
NFC House
Vickers Industrial Estate
Mellishaw Lane, Morecambe
Lancashire LA3 3EN
tel. +44 (0) 1524 847600
gvsuk@gvs.com

ASIA

China
GVS Technology (Suzhou) Co., Ltd.
Fengqiao Civil-Run Sci-Tech Park,
602 Changjiang Road, S.N.D.
Suzhou, China 215129
tel. +86 512 6661 9880
fax: +86 512 6661 9882
gvschina@gvs.com

Japan
GVS Japan K.K.
KKD Building 4F, 7-10-12 Nishishinjuku
Shinjuku-ku, Tokyo 160-0023 Japan
tel. +81 3 5937 1447
fax +81 3 5937 1448
gvsjapan@gvs.com

India
India Office
V30, 14th Street, Anna Nagar,
Chennai - 600040 India
Tel. +91 98840 58375
gvsindia@gvs.com

AMERICA

U.S.A.
GVS Maine
63 Community Drive
Sanford, ME 04072 - USA
tel. +1 866 7361250
www.mfgmaine.com
info@mfgmaine.com

GVS Filter Technology Inc.
5353 West 79th Street
Indianapolis, IN 46268 - USA
tel. +1 317 471 3700
fax +1 317 471 8370
gvsusa@gvs.com

Brazil
GVS do Brasil Ltda.
Rodovia Conego Cyriaco Scaranello Pires 251
Jd. Progresso, CEP 13190-000
Monte Mor (SP) - Brasil
tel. +55 19 38797200
fax +55 19 38797251
gvs@gvs.com.br

México
México Office
Paseo de Santa Anita 548
Condominio Santa Anita
45645 Tlajomulco de Zuñiga (JAL) - México
tel. +52-33 3110 0844
gvs_mex@gvs.com

Argentina
GVS Argentina S.A.
Parral 246-9° A
1405 Buenos Aires - Argentina
tel. +54 11 49889041
fax +54 11 49889042
gvsarg@gvs.com